

FIRMWARE COMPATIBILITY

MP150 Firmware—AcqKnowledge for Windows Software	Compatible with AcqKnowledge
v 1.1.22† † Minor fix for averaging rejection functionality	AcqKnowledge 4 for Windows Released 04.15.2008
v 1.1.14 – v 1.1.21† † v1.1.21 fixed Mac compatibility for older stimulation functions.	AcqKnowledge 3.9.1 for Windows Released 6.29.2007 AcqKnowledge 3.9.0 for Windows Released 11.17.2006
v 1.1.14 – v 1.1.21† † v1.1.21 fixed Mac compatibility for older stimulation functions.	AcqKnowledge 3.8.2 for Windows Released 02.21.2006
v 1.1.12 (v 1.1.10, v 1.1.11) For AcqKnowledge 3.8.1 installed with MP150 loaded with firmware v1.1.12, install AcqKnowledge 3.8.2 and execute the firmware update from installed AcqKnowledge 3.8.2. <ul style="list-style-type: none"> <li>• <b>DO NOT</b> just download the latest MP150Tools.exe to execute a firmware update because the installer is not compatible with AcqKnowledge 3.8.1.</li> </ul> With better DHCP client?	AcqKnowledge 3.8.1 for Windows Released 08.20.2004
v 1.1.9 For AcqKnowledge 3.7.3 installed with MP150 loaded with firmware v1.1.9, user needs to purchase a software upgrade to AcqKnowledge 3.8.2 and execute the firmware update from installed AcqKnowledge 3.8.2. <ul style="list-style-type: none"> <li>• <b>DO NOT</b> just download the latest MP150Tools.exe to execute a firmware update because the installer is not compatible with AcqKnowledge 3.8.1.</li> </ul>	AcqKnowledge 3.7.3 for Windows Released 12.25.2002
v 1.1.8 If you have AcqKnowledge 3.7.2 or previous installed (DLC and USB support ONLY) and want to upgrade MP150 firmware from DLC-support ONLY to UDP-support, you must follow a very specific procedure and you will need to have the original 3.7.2 CD or use the MPCCommType utility to switch modes. <ul style="list-style-type: none"> <li>• See Support &gt; Updates at www.biopac.com for detailed instructions.</li> </ul>	AcqKnowledge 3.7.2 for Windows Released 2001
v 1.1.8 ONLY <ul style="list-style-type: none"> <li>• First release with UDP</li> </ul>	AcqKnowledge 3.7.1 for Windows Released 2000
v 1.1.7 and below	Please contact BIOPAC to review firmware and software compatibility and upgrade options.
BioHarness™ Firmware	v2.0.0.14 Upgrade released 2010