

MP150 Firmware Upgrade—AcqKnowledge 3.9


Before performing a Firmware Upgrade, make certain that no other users are connected to or using the MP150 you plan to upgrade. Any attempts to communicate with the MP150 via any user's *AcqKnowledge* software may affect the viability of the upgrade. Fluctuations in the power source during an upgrade may also affect the outcome.

To use an MP150 with UDP communication on a network with a non-Windows DHCP server, you must use firmware 1.1.12 or greater for the MP unit to properly be assigned an IP address. This is also true for any DHCP system with non-Windows operating systems, such as Unix, Linux, Mac OS X, and other DHCP-aware devices.

- *AcqKnowledge* 3.8.2 use MP150TOOLS.EXE to update; see separate instructions.

MP150 Firmware Update Procedure

1. Launch *AcqKnowledge* 3.9.0 or higher.
2. Select MP150 menu > Update Firmware...
3. Click "OK" at the confirmation prompt.
4. Navigate to the MP150.mpb file and click "Open."
 - The default location is C:\Program Files\BIOPAC Systems Inc\AcqKnowledge 3.9.0.
5. Click "Start" to upload the selected firmware file.
6. Wait for the confirmation prompt (usually less than 30 seconds) and then click "OK."

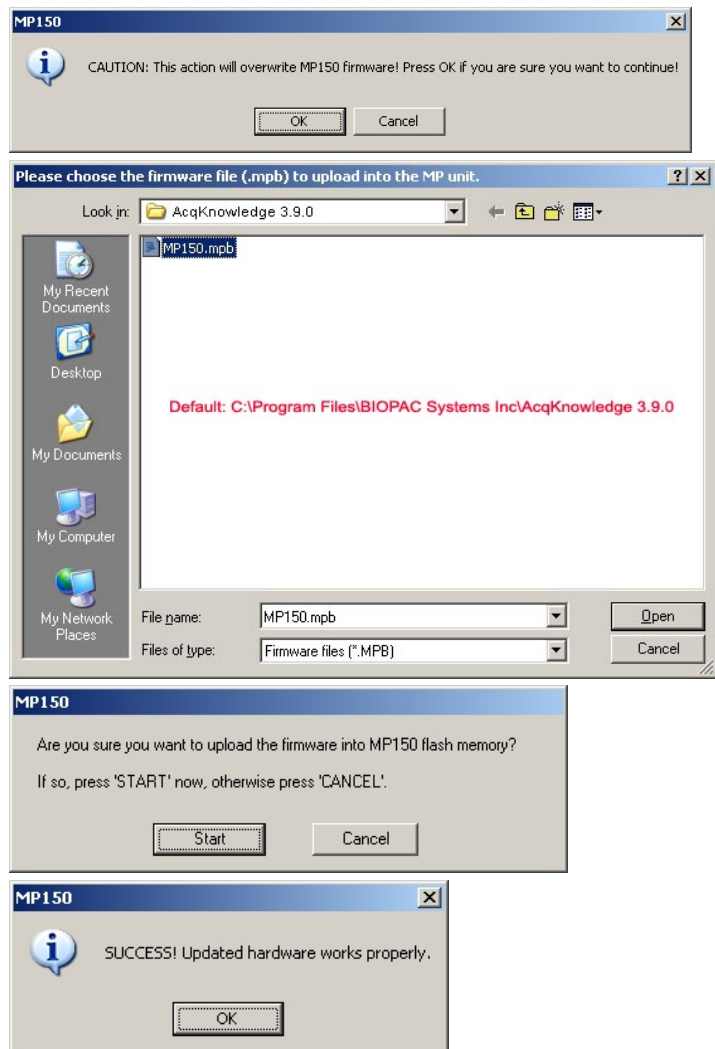

The MP150 is delivered with a factory-programmed firmware version that is not affected by the upgrade procedure—you can always restore the original firmware with the Firmware Rollback Switch.