

## REQUEST FOR QUOTE

Hi,

We've ordered from you before, but some of our items/quantities don't fit your plant so I wanted to check if you could be competitive for this run:

BSL Catalog (see our current catalog here if it helps to quote: [BSL Catalog RFQ.pdf](#))

- Page size 11x8.5
- Pages 52 (cover +48)
- Paper 80# text, gloss text
- Cover 100# text, gloss text
- 4-color both sides
- Qty 1,000 or 1,500
- Turn-around 7 business days
- Proof - Electronic only
- Aqueous coating front and back
- saddle stitched
- no fold
- shipping to Goleta, CA 93117

Competitive pricing is currently approx. \$4,725 for 1,000 or \$5,185 for 1,500.

Please advise if you would like to submit a quote.

Thanks!

**Jocelyn Mariah Kremer**

jocelyn@biopac.com



805.685.0066 x128



biopac

# student lab

an integrated system of hardware, software and a lab manual that **revolutionizes** life science education

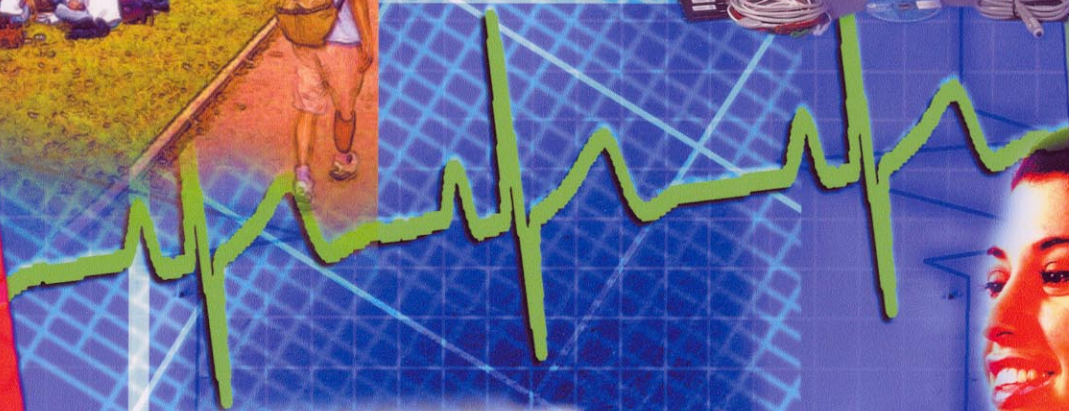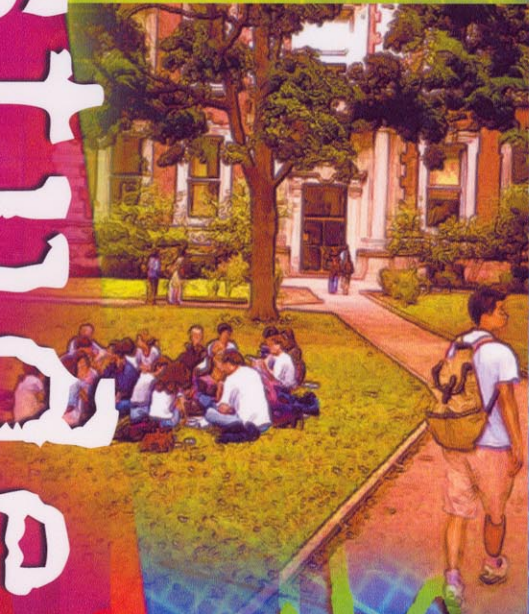

**BIOPAC**  
Systems, Inc.

Registered to ISO 9001:2000




# biopac student lab

Research-Quality Tools for Education



*Dynamic lessons engage students*

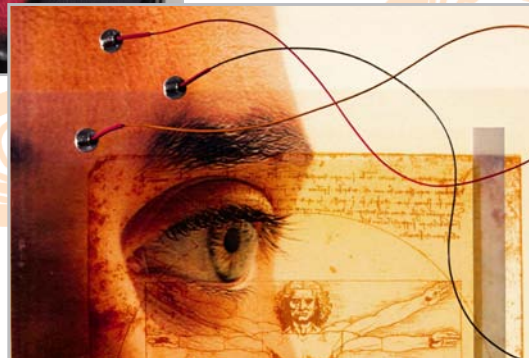

The #1 Choice for Your Lab  
Just Keeps Getting Better

- Curriculum
- Laboratory Disciplines
- Active Learning
- System Level Solutions
- State-of-the-Art Technology

# Table of Contents

Contact BIOPAC today and join thousands of satisfied users, like this one:

“Of the five major players in the market whose products we selected to examine, either by purchasing and using one complete system, or by demonstration, only one measured up. The Biopac Student Lab.”

—Comment from the HAPS List Server



New budget-beating packages

## All New BSL System

- **MP35 Hardware**
- **New Software**
- **New Lessons**

	Page
<b>S BSL System Overview</b>	<b>2</b>
<b>A Applications</b>	<b>8</b>
<b>Lab Manuals</b>	<b>8</b>
<b>Cross-Discipline</b>	
Basic System	8
Advanced System	9
Ultimate System	9
<b>Core Discipline</b>	
Human Physiology	10
Animal Physiology	12
Biology	14
Exercise Phys. & Biomechanics	16
Psychophysiology & Neurophysiology	18
Biomedical Engineering (BME)	20
Pharmacology & Toxicology	22
<b>H BSL Hardware</b>	<b>23</b>
Data Acquisition Unit	24
Stimulators	25
Transducers (“SS Series”)	26
Stimulus Presentation	32
Noninvasive Blood Pressure	32
Tissue Bath Stations	33
Airflow & Gas Analysis	33
Electrodes & Accessories	36
Recording Options	36
Stimulating Options	38
Interface Cables	39
Transducer Interface Options	39
Package Overview	40
Core Package Components	41
Transducer Accessory Packs	42
<b>L Lessons Summary</b>	<b>43</b>
Lesson Hardware Guide	46
<b>I Index</b>	<b>47</b>

 **BIOPAC** Systems, Inc.



QUALITY SYSTEM  
REGISTERED TO  
ISO 9001:2000

42 Aero Camino, Goleta, CA 93117 USA  
Tel (805) 685-0066 Fax (805) 685-0067  
www.biopac.com info@biopac.com

# the #1 solution for life science education just keeps getting better!



The Biopac Student Lab System is an integrated solution that includes hardware and software that students use to record data from their own bodies, animals or tissue preparations. Also includes 60+ life science lesson experiments and easy instructions to modify them or integrate your existing syllabus. The BSL is #1 in CLASS!



## Curriculum

- Total of 60+ lesson experiments
- Comprehensive support from published Lab Manuals—  
Marieb: PEARSON Benjamin Cummings, Fox: WCB/McGraw-Hill,  
Wood: PEARSON Benjamin Cummings, Pflanzler: Kendall/Hunt
- Video demonstration clips
- On-line lessons
- Distance learning
- Suitable for 2- & 4-year programs, medical schools and nursing programs
- Lesson development support from BIOPAC

## Laboratory Disciplines

- Physiology (animal and human)
- Exercise Physiology and Biomechanics
- Psychophysiology and Neurophysiology
- Pharmacology and Toxicology
- Human Anatomy
- Biology
- Bioengineering
- And more!



*New Gas Analysis lessons*

## Active Learning

- Students as subjects
- Inquiry-based, hypothesis-driven lessons
- Easy lesson customization for student experiment design
- Application Notes for advanced experiment design

**total support**

## BIOPAC STUDENT LAB — *Total Solutions...with Total Support!*

**On-line Help** – Quick Guide, Tutorials, searchable manuals (pdf), *PRO* Lessons

**Video Help** – movie clips of procedures and sample data

**Data Examples** – files for male and female subjects

**Web Link** – use your active browser to link to the full spectrum of support at [www.biopac.com](http://www.biopac.com)

**Desktop Streaming** – like having an on-site rep...this option allows remote viewing and mouse control so we can see your data and review your setup for targeted Tech Support!

**Phone Support** – qualified BSL Specialists provide telephone support

**Lab Manuals** – use the BSL Laboratory Manual or choose from major published A&P Lab Manuals





## System Level Solutions

- Hardware—Data acquisition unit, transducers & electrodes
- Software—Building-block structure for guided lessons to open-ended discovery
- Advanced Programming Options—Application Programming Interfaces available to control MP35 hardware or access BIOPAC data files
- Curriculum Library & Extensive support materials
- Grant-friendly solutions—Recognized by the NSF, NIH, The Whitaker Foundation, and most major grant foundations
- Worldwide customer base—Thousands of users
- Multi-lingual support: English, French, Spanish, Japanese...and more on the way!



New Gas Analysis Module



New Tissue Bath



New Stimulator

## State-of-the-Art Technology

- Students learn real-world skills using clinical recording techniques
- Industry standard transducers—Over 60 currently available
- Tissue bath stations—Modular system with integrated heater
- Gas analysis module for CO<sub>2</sub> and O<sub>2</sub> measurements
- Stimulus generation functionality
- Digital control channels—8 Input/8 Output
- Visual presentation options—SuperLab, E-Prime, MediaLab, DirectRT, Inquisit
- Support for the latest operating systems
- New MP35 acquisition unit: 24-bit resolution, USB, high speed
- Interface with all major amplifier and transducer manufacturers—Use our ready-made connectors or the custom kit
- Continuous product development

## dynamic lessons

The student-friendly BSL System reduces setup time by up to 90%, so students collect excellent data and focus on scientific principles. The multi-level learning features of the BSL software let you control the material and method of each experiment—you can even use the BSL for graduate programs and advanced research.

- Use the 18 guided BSL Lessons for introductory concepts
- Choose from 40+ PRO Lessons for introductory & advanced concepts
- Add analysis tools to existing Lessons
- Let students develop experiment protocols
- Modify a Lesson or incorporate existing lab procedures
- Easily create new experiments

BSL lesson experiments are included in a number of the leading published Lab Manuals and have been successfully used to study:

- |       |                               |                      |                  |         |
|-------|-------------------------------|----------------------|------------------|---------|
| • ECG | • EDA (GSR)                   | • Heart Sounds       | • Reaction Time  | • Force |
| • EEG | • Biofeedback                 | • Nerve Conduction   | • Respiration    |         |
| • EMG | • Bioimpedance/Cardiac Output | • Pulmonary Function | • Temperature    |         |
| • EOG | • Blood Pressure              | • Gas Analysis       | • Metabolic Rate |         |

- Engage students with dynamic lessons as they record data from their own bodies
- New Budget-Beating packages
- Customize lessons from BIOPAC's extensive Curriculum Library
- Incorporate clinical techniques in the lab—Students learn real-world skills
- Combine physiological data with stimulus presentation (Analog, Digital, and Calculation channels)
- No complicated knobs or dials
- No messy ink and paper

# the gold standard for

## BIOPAC STUDENT LAB The Total Solution



Engage your students!

## over 1.6 million successful lab hours

Since its introduction in 1995, the Biopac Student Lab has evolved based on input from many thousands of existing users and physiology instructors. To see for yourself why the BSL offers the most meaningful, flexible lab experience for students and teachers, visit [www.biopac.com](http://www.biopac.com) and check out demo video clips and sample lessons.

Give your students a proven advantage...give them the Biopac Student Lab!

## New software tools

### new hardware features


- Increased performance (24-bit A/D converter)
- Expanded amplifier settings
- New Stimulator and Output Controls
- Improved "Electrode Check" for electrode impedance
- New digital input & output capabilities
- Increased triggering options

- New signal Presets to simplify lesson creation—now more than 100 available
- Expanded Help menu includes more student prep and support tools
- More Powerful Markers—distinguish appended segments from events, pre-label markers and assign to function keys, layer printed markers to prevent print overlap, etc.
- Rewind button—delete last or all recorded segments
- "Snap-to" cursor for one data point in X/Y mode
- Output Control enhancements (for Stimulator, Pulses, Digital, Voltage, Analog CH), including toolbar controls
- Save multiple stimulus paradigms within a Template or data file
- Improved real-time zoom functionality for fast-response data
- New overlap appended segments display mode
- New standard curve tools for colorimetry data
- Lesson journal text auto-sized for monitor resolution
- Lesson calibration held for consecutive subjects
- Save options include network servers so students can access files outside of the lab
- Access privileges installed to all users
- Digital registration for support and update info

## Instructor support materials

### Help your students get good data

- The Biopac Student Lab System includes videos of lesson procedure and recorded data, tutorials, sample data files, onscreen prompts and a detailed lab manual to help students acquire high-quality data
- Extensive curriculum library for human, animal, organ, and tissue lesson experiments
- Easy customization to suit your curriculum—create your own lessons with no programming required
- Language options: English, French, Spanish, Japanese
- Biopac Student Lab Specialists available for Technical Support
- **Lab Manual Options**
  - Use the Biopac Student Lab Manual included with each BSL System for 17 scientific lessons
  - Use digital text of the lessons and create a custom lab manual with your own curriculum
  - Select from major published Lab Manuals using BIOPAC lesson experiments (see page 8).
- **Data comparison tools for increased analysis**
  - Normative values provided for FEV<sub>1,2,3</sub>, MVV, BP and ECG
  - Arrhythmia lesson, including sample data files from the NIH PhysioNet database
  - Sample data files for male and female subjects
  - Easily export for statistical analysis or archive student data

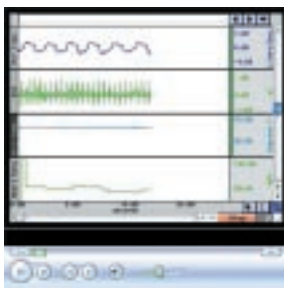

Streaming data videos help students identify signal characteristics for each lesson



Videos to help students prepare for lab time are included in the Help menu and online

when you get the bsl...they'll get science!

# life science labs



simplify your labs & empower your students

## The BSL System helps students focus on scientific principles with...


- Unmatched guidance, power and flexibility
- Familiar point-and-click interface
- Extensive support materials to prepare for lab time, including videos & sample data files
- Onscreen prompts—additional instructions help students record the highest quality data
- On-line instructions—guide the student through the lesson
- Simple sensors—validate transducer and signal connections
- Multiple data displays—view data in industry standard formats: X/Y, overlap, clinical grids, chart recorder, oscilloscope, standard curve, dot display
- Adaptable lessons—follow the lesson tasks or assign new tasks to suit your course requirements
- Adaptable protocols—students can easily design experiments to develop and test their hypothesis with the BSL
- Data analysis and measurement tools—extract baseline and experimental response values and compare them to normative values
- Professional data reports—copy screen shots into lab reports and measurement values into statistical programs, or create high quality printouts with clinical grids, markers and full color
- Analysis options—distribute the analysis software so students can analyze data outside the lab—at home, library, etc.
- Reporting and analysis options—automatically paste measurements and graphs to the journal or other programs.
- Automated calibration—optimizes the display settings for the subject's own signal
- Autosave feature for lessons—includes student's name and lesson number and allows instructors to set file storage location

*Intuitive interface with buttons and pop-up menus. Show or hide data channels and control size, color, label, scale, etc. Add markers and measurements.*


*On-screen prompts and expanded Help menu help students get good data with instructions and videos for setup, calibration, recording and file management.*

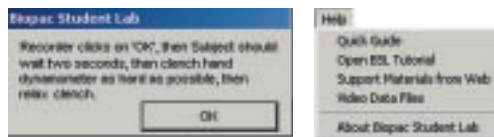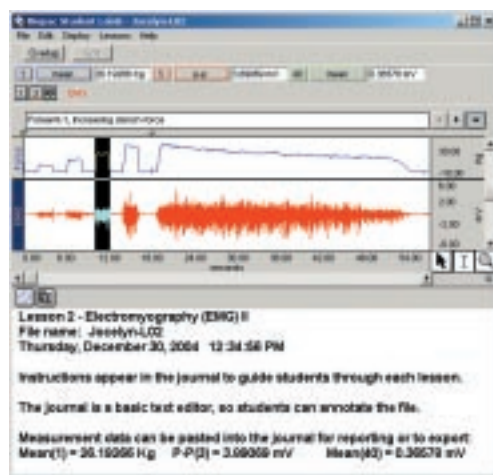

*Sophisticated measurement tools help students extract—and easily report—meaningful data*

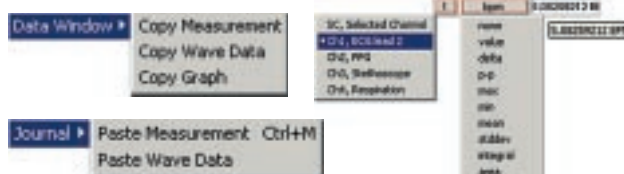

*Data viewing options allow students to optimize display for signal type and include measurements, markers, grids, etc. for analysis or reporting*

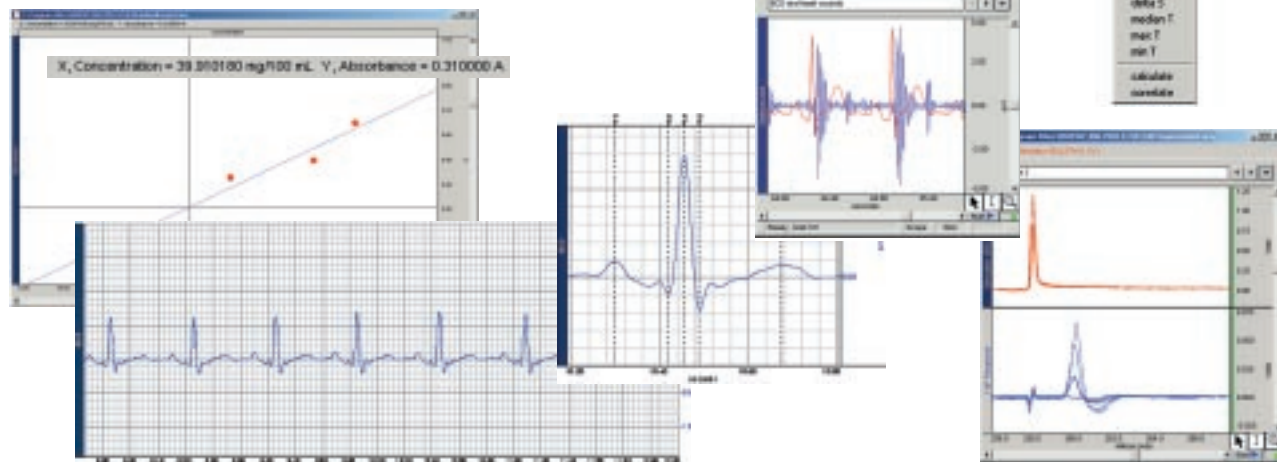



# facilitate learning & enhance



A BSL system was used in a study aboard the "Vomit Comet" in the Zero-G program at the NASA Johnson Space Center by Dr. Rick Puzdrowski and students from the University of Houston—Clear Lake School of Natural and Applied Sciences. Students Carolyn Jessop (left) and Sonya Morgan with JSC techs (in blue).

Source: NASA Johnson Space Center. <http://zerog.jsc.nasa.gov>

The BSL System helps students get on track right away with data acquisition and analysis. The hardware has no knobs or dials and the lesson software uses automatic calibration and onscreen prompts. These features eliminate confusion and frustration and save a lot of time in the lab. More students can record data, which makes the whole laboratory experience much more productive.

For open-ended learning, use the BSL System for student projects and research, or use it to quickly create your own lessons.

- Adapt student task protocols to develop critical thinking skills
- Modify or expand BSL Lesson segments for hypothesis-driven studies

## What will you do with the Biopac Student Lab?

Quick-click curriculum options let you use the lessons any way you want...

- Use existing BSL Lessons to study the fundamentals of physiological systems like cardiovascular and pulmonary function
- Use 40+ BSL PRO Lessons developed by instructors to explore principles like Habituation, Impedance Cardiography, and Nerve Conduction—downloads include lesson setup, procedure and data template
- Customize BSL lessons to incorporate your protocols, add advanced analysis tools, and/or use existing equipment
- Develop lessons with new or existing curriculum and BSL software tools like FFT, Histogram, Trigger, Averaging, and Rate Detection. Add setup & onscreen instructions and then save as a template
- Allow students to design their own research projects. Use BSL Lessons as a starting point for inquiry-based experiments, then modify or expand recording parameters and protocols to test hypotheses

## Lesson development made easy—for you and your students

Create your own lesson in just a few simple steps...

- Pick from more than 100 signal and calculation channel Presets
- Establish the recording parameters (sample rate, duration of experiment/lesson)
- Add instructions to the journal—students will follow your lesson plan
- Save the file as a template—students open the file with your settings in place




Presets simplify setup for custom projects

- Presets include all amplifier and filter settings for the signal to optimize data collection and display
- Presets can be used as is, or customized and saved under a new preset name for your protocol
- Organize and/or limit the list of Presets displayed to further simplify setup



Signal Presets

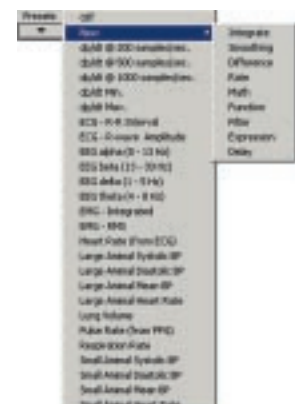


Calculation Presets

# the lab experience

the biopac student lab—student-friendly & research-compatible


The multi-level learning features of the Biopac Student Lab, along with its proven success in thousands of labs worldwide, make the BSL the ideal teaching solution. BSL Lessons cover scientific fundamentals and simplify recording and analysis so your students can focus on lab objectives. As students advance, BSL recording and analysis options can be expanded to allow for sophisticated measurements, student-directed experiments, alternate curriculum and research-level studies.

Show or hide menus and analysis features based on the lesson requirements

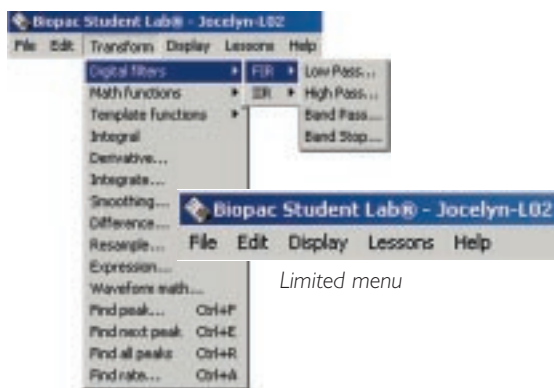


Limited menu

Expanded menu with increased analysis tools

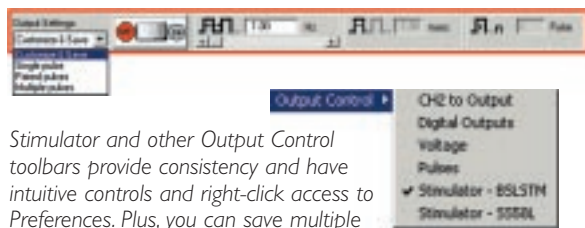

Stimulator and other Output Control toolbars provide consistency and have intuitive controls and right-click access to Preferences. Plus, you can save multiple stimulus paradigms within a file.

Markers layer for printing to prevent overlap. Automatic stimulus markers log pulse width and rate.

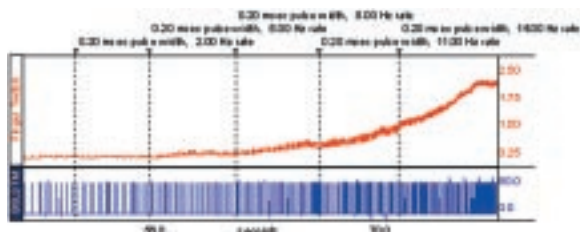

To enhance any lesson, consider these options...

## Control:

- Listen to the sound of EMG and other signals to emphasize a concept
- Utilize multiple display modes to optimize visual feedback for the signal type, in real time or offline. Select chart, scope, or X/Y plot, overlap segments or channels, plot standard curve, histogram or FFT, display as vertical or horizontal bar graphs or numeric values, show filtered EEG data or calculate heart rate.
- Emulate standard lab equipment displays—chart recorder, oscilloscope, X/Y plotter, and metabolic cart
- Set up digital I/O channels to receive trigger information from products like SuperLab or E-prime, or control pumps, lights, etc.
- Select automatic file saving—save to a network server so students can access and analyze data outside the lab
- Control output functions such as sound, digital pulses and stimulation (low and high voltage)
  - Use stimulation features for human or animal subjects (visual, auditory, mechanical, or electric)
  - Set the stimulator to automatically start with the recording or for manual start and stop within the experiment
  - Set automatic stimulator event markers to note the stimulus frequency and pulse width
- Use filtering options to allow students to develop and display the filter response—online and offline
- Choose External or Analog (from any channel) triggering options

## Analysis:

- Perform online measurements of heart rate, systolic, diastolic, mean blood pressure, dP/dt max and min with the rate detection features
- Identify frequency components of a signal—i.e., EEG alpha, beta, delta—with FFT functions
- Use the Equation Generator to calculate values such as RER and stroke volume
- Extract visual and auditory responses with signal averaging tools
- Use advanced integration features for pulmonary function studies
- Make metabolic studies easy with BSL software tools and the CO<sub>2</sub> & O<sub>2</sub> Analysis Module (GASSYS2, page 34)
- Calculate stroke volume and cardiac output with data from the Cardiac Output Sensor (SS31L, page 29)
- Select tools for spike sorting, spike counting and automatic measurements

## Customization:

- Customize menu and toolbar options to show or hide analysis and editing controls based on the lesson requirements
- Add “BSL PRO” tools to any lesson for advanced analysis
- Customize the grid display—easily create clinical ECG grids
- Select the measurements that students must use to analyze the data
- Add text to the Journal so students have on-line access to your instructions
- Establish marker options to predefine recording segments or allow students to add comments to files
  - Predefine segment marker text—the marker describes each new recording segment (sit up, lay down)
  - Predefine event markers—set function keys to mark and describe events/segments in the experiment
- Set up acquisition and recording automation—repeat, trigger, autosave



# scientific applications & core packages

The Biopac Student Lab System is the total solution for a wide variety of scientific applications, including human, animal, organ and tissue studies. BSL Systems are available for general science—Basic, Advanced and Ultimate Systems—and core disciplines. Each Biopac Student Lab System includes a high performance data acquisition unit plus electrodes, transducers and consumable items required to run a specified number of lessons for each particular discipline. Any package can be used as a starting point to develop a custom system. Electrodes and transducers can be purchased separately in order to perform additional lessons not covered with the core package, and hardware upgrade kits are available.

- The BASIC, ADVANCED and ULTIMATE Systems have been offered since the Biopac Student Lab was launched in 1995, and are now in use in thousands of labs worldwide in 2-year and 4-year programs, medical schools, nursing programs, veterinary programs, chiropractic colleges, etc.
- The new CORE packages are offered to meet the diverging requirements of specific departments:

Human Physiology, page 10  
 Animal Physiology, page 12  
 Biology, page 14  
 Exercise Physiology & Biomechanics, page 16  
 Psychophysiology & Neurophysiology, page 18  
 Biomedical Engineering, page 20  
 Pharmacology & Toxicology, page 22

Suggested applications are outlined on the following pages, along with a CORE package with hardware targeted for each discipline. A BSL Lab Manual is included with each system, or you can choose from major published lab manuals that incorporate BSL experiments. Plus, *PRO* Lessons are available online, and you and your students can develop new applications.



BIOPAC MANBSL3S  
 One manual is included with each system. Covers BSL Lessons 1-17 (page 40).




ISBN: 0-7575-0386-1



Main 7/e ISBN: 0-8053-5514-6  
 Cat 8/e ISBN: 0-8053-5516-2  
 Pig 8/e ISBN: 0-8053-5518-9



Main ISBN: 0-8053-7368-3  
 Cat ISBN: 0-8053-7369-1  
 Pig ISBN: 0-8053-7370-5

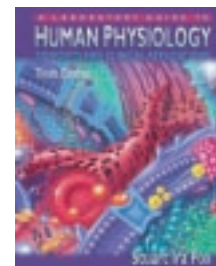

ISBN: 0-07-244087-2  
 Also ISBN: 0-07-027236-0 (9th)



Auteur : Sous la direction de Michel DAUZAT  
 Editeur : SAURAMPS  
 ISBN : 2840234041

A

8

## basic system

These hardware suggestions will enable you to perform a variety of lessons for human subjects. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the Basic System or select items à la carte. If you start with the Basic System, hardware upgrades are available to create Advanced or Ultimate Systems. See BSL Hardware (page 23) for all available transducers, electrodes and accessories.

### Basic System BSLBSC-W (Win) or BSLBSC-M (Mac)

Data Acquisition Unit	MP35, p. 24	Electrode Lead Sets (2x)	SS2L, p. 25	USB Cable	CBLUSB, p. 40
BSL Software	3.7 for Lessons & <i>PRO</i>	Disposable Electrodes (100/pk)	EL503, p. 37	Power Transformer	AC300A, p. 40
BSL Lab Manual	MANBSL3S, p. 40	Abrasive Pads (10/pk)	ELPAD, p. 36		

### Perform 19 or more lessons with the BASIC System:

#### Muscular

<b>BSL1</b>	Standard & Integrated EMG
<b>H07</b>	EMG Contractions—Active Learning
<b>H27</b>	Facial EMG
<b>H34</b>	Electrogastrogram (from human)

#### Cardiovascular

<b>BSL5</b>	Components of the ECG (Lead II)
<b>BSL6</b>	Leads I, II, III & Einthoven's Law
<b>H08</b>	Dive Reflex—Active Learning
<b>H23</b>	Signal Averaged ECG
<b>H32</b>	Heart Rate Variability

#### Bioengineering

<b>H20</b>	Filtering
<b>H33</b>	FFT Fast Fourier Transform

#### Neurophysiology

<b>BSL3</b>	EEG Relaxation & Brain Rhythms
<b>BSL4</b>	Alpha Rhythms in the Occipital Lobe
<b>BSL10</b>	Eye Movement, Saccades & Fixation
<b>H10</b>	EEG & Hemispheric Asymmetry
<b>H12</b>	EOG Saccades & Displacement
<b>H13</b>	EOG Visual Tracking
<b>H14</b>	Ocular Fixation while reading
<b>H15</b>	Ocular Fixation while viewing an image

# advanced system

These hardware suggestions will enable you to perform a variety of lessons for human subjects. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the Advanced System or select items à la carte. A hardware upgrade kit is available to upgrade to the Ultimate System. See BSL Hardware (page 23) for all available transducers, electrodes and accessories.

<b>Advanced System BSLADV-W (Win) or BSLADV-M (Mac)</b>	Airflow Nose Clips (10/pk)	AFT3, p. 34	Surgical Tape	TAPE1, p. 36
Basic BSL System	Calib. Syringe (600 ml)	AFT6, p. 34	Respiratory Effort Trans.	SS5LB, p. 26
Airflow Transducer	EDA (GSR) Lead	SS57L, p. 26	Temp Trans.	SS6L, p. 26
Airflow Filters (10/pk)	EDA Electrodes (100/pk)	EL507, p. 37	Hand Switch	SS10L, p. 26
Airflow Mouthpieces (10/pk)	Colored Paper	PAPER1	Headphones	OUT1, p. 26
			Pulse Plethysmograph	SS4LA, p. 26

## Perform 33 or more lessons with the Advanced System:

### Muscular

BSL1	Standard & Integrated EMG
H07	EMG Contractions—Active Learning
H27	Facial EMG
H34	Electrogastrogram (from human)
H36	Muscular Biofeedback

### Cardiovascular

BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
H05	WAnT Wingate Test
H08	Dive Reflex—Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

### Pulmonary Function

BSL8	Respiratory Cycle
BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV
BSL15	Aerobic Exercise Physiology

### Neurophysiology

BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL9	GSR and Polygraph
BSL10	EOG Eye Movement, Saccades & Fixation
BSL11	Reaction Time & Learning
BSL14	Biofeedback: Relaxation & Arousal

### H10

H10	EEG & Hemispheric Asymmetry
H11	Mirror Test—EDA Sensory motor learning
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image
H16	Reaction Time
H24	Habituation

### Bioengineering

H02	Compartmental Modeling
H20	Filtering
H33	FFT Fast Fourier Transform

See page 43-45 for a description of all available lessons.

Contact BIOPAC to discuss your specific needs. [www.biopac.com](http://www.biopac.com) • [info@biopac.com](mailto:info@biopac.com)

# ultimate system

These hardware suggestions will enable you to perform a variety of lessons for human and animal subjects. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the Ultimate System or select items à la carte. See BSL Hardware (page 23) for all available transducers, electrodes and accessories.

<b>Ultimate System BSLULT-W (Win) or BSLULT-M (Mac)</b>	EDA Electrodes (100/pk)	EL507, p. 37	Pressure Transducer	SS13L, p. 27
Basic BSL System	Colored Paper	PAPER1	BP Cuff Transducer	SS19L, p. 27
Airflow Transducer	Surgical Tape	TAPE1, p. 36	Hand Dynamometer	SS25LA, p. 27
Airflow Filters (10/pk)	Respiratory Effort Trans.	SS5LB, p. 26	Force Transducer	SS12LA, p. 27
Airflow Mouthpieces (10/pk)	Temp Trans.	SS6L, p. 26	Electronic Stethoscope	SS30L, p. 28
Airflow Nose Clips (10/pk)	Hand Switch	SS10L, p. 26	Multi-lead ECG Cable	SS29L, p. 27
Calib. Syringe (600 ml)	Headphones	OUT1, p. 26	BSL Stimulator	BSLSTM, p. 25
EDA (GSR) Lead	Pulse Plethysmograph	SS4LA, p. 26	Stim. Electrode for humans	HSTM01, p. 28
	Electrode Gel	GEL1, p. 36		

## Perform 43 or more lessons with the Ultimate System:

### Muscular

A05	Visceral Smooth Muscle
BSL1	Standard & Integrated EMG
BSL2	Motor Unit Recruitment & Fatigue
H06	Finger Twitch
H07	EMG Contractions—Active Learning
H27	Facial EMG
H34	Electrogastrogram (from human)
H36	Muscular Biofeedback

### Cardiovascular

A04	Frog Heart
BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
BSL16	Blood Pressure & Korotkoff
BSL17	Heart Sounds & Cardiac Events
H01	12-Lead ECG

H04	Blood Pressure (Isometric or Straining exercise)
H05	WAnT Wingate Test
H08	Dive Reflex—Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

### Pulmonary Function

BSL8	Respiratory Cycle
BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV
BSL15	Aerobic Exercise Physiology

### Neurophysiology

A01	Frog Pith & Prep
BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL9	GSR and Polygraph


BSL10	EOG Eye Movement, Saccades & Fixation
BSL11	Reaction Time & Learning
BSL14	Biofeedback: Relaxation & Arousal
H03	Nerve Conduction (ulnar nerve)
H10	EEG & Hemispheric Asymmetry
H11	Mirror Test—EDA Sensory motor learning
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image
H16	Reaction Time
H24	Habituation

### Bioengineering

H02	Compartmental Modeling
H20	Filtering
H33	FFT Fast Fourier Transform



# human physiology

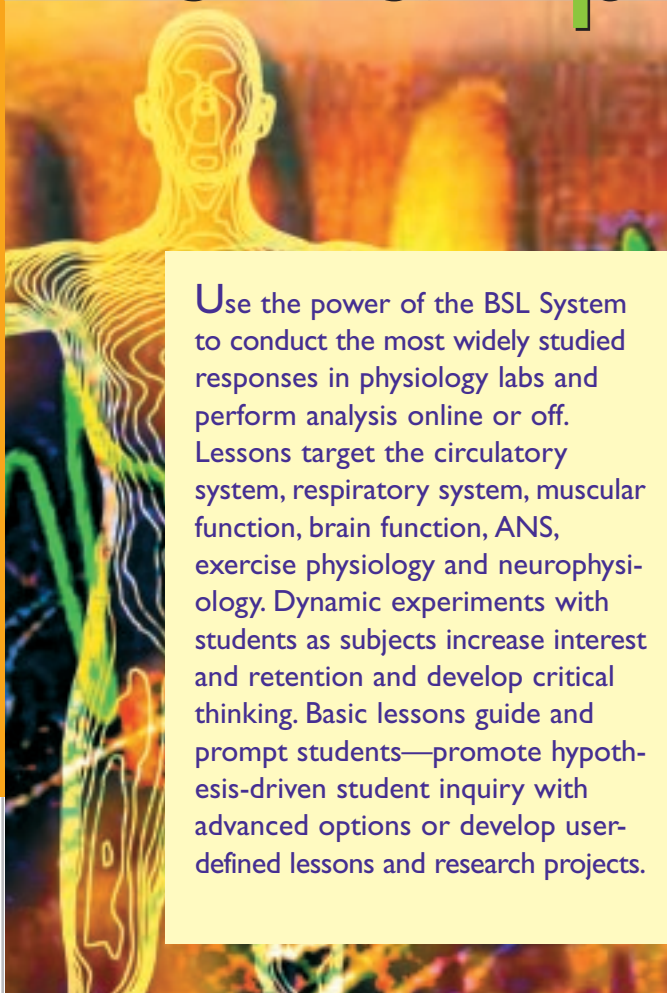

Use the power of the BSL System to conduct the most widely studied responses in physiology labs and perform analysis online or off. Lessons target the circulatory system, respiratory system, muscular function, brain function, ANS, exercise physiology and neurophysiology. Dynamic experiments with students as subjects increase interest and retention and develop critical thinking. Basic lessons guide and prompt students—promote hypothesis-driven student inquiry with advanced options or develop user-defined lessons and research projects.

We have been using the Biopac Student Lab system in our Human Anatomy & Physiology labs at IU South Bend for about 5 years. Students enjoy being able to analyze data that was generated from a member of their lab group. It certainly has more meaning to them than using traditional "canned" lab data.


—Mary Truex, Indiana University South Bend, Biology Department

## Cardiovascular

The BSL System covers a wide range of cardiovascular measures based on the heart's electrical signal. Record standard (LI, LII or LIII) or augmented (aVR, aVL or aVF) leads; add a multi-lead ECG cable to simultaneously record a pre-cordial lead. Students attach Leads I and III and the software uses Einthoven's law to display Lead II. Identify, isolate and measure components of the ECG complex under varying conditions. Use the averaging features to automate ECG analysis. View arrhythmia data obtained from the NIH PhysioBank archive. Calculate R-R interval to display heart rate variability or determine vagal tone. Use the electronic stethoscope to listen to and record heart sounds and then correlate the sounds with the mechanical and electrical events of the cardiac cycle. Record cuff pressure and Korotkoff sounds to measure systolic and diastolic blood pressure. Determine pulse wave velocity. For advanced studies, use the noninvasive cardiac output sensor to record stroke volume and cardiac output.




Overlapped ECG and heart sounds

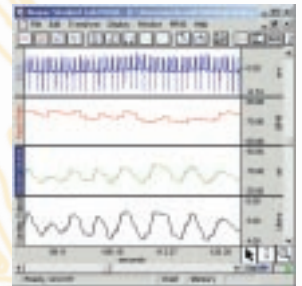

Cardiac output and stroke volume

## Respiratory & Pulmonary Function

The BSL provides an excellent introduction to volumes and capacities, respiratory flow rates, breathing mechanics, and ventilation. Display and/or print a clinical grid on the data for effective interpretation and training. The software guides the student in measuring FVC and in calculating FEV and MVV. The airflow transducer is handheld, lightweight, easy to clean and very easy to use. Take measures before,

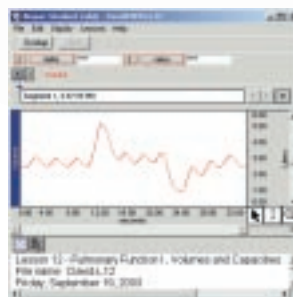

Pulmonary volume

during and after exercise to study ventilation and heat exchange. The new Gas Analysis module provides online measures of CO<sub>2</sub> and O<sub>2</sub> levels for Respiratory Exchange Ratio, O<sub>2</sub> Consumption and Basal/Resting Metabolic Rate lessons. Non-rebreathing T-valves, air chambers, facemasks, and tubing options provide setup options to suit any protocol.

## features

- 25 lessons targeted for Human Physiology
- ECG, EEG, EMG, EOG & EGG
- Pulmonary Function FEV<sub>1,2,3</sub>, MVV, PV Loops
- Airflow
- Blood Pressure
- Heart Sounds & Korotkoff Sounds
- Nerve Conduction
- Electrodermal Activity (GSR)
- Pulse
- Reaction Time
- Temperature
- Stimulation & Response (Somatic Reflexes)
- Gas Analysis (O<sub>2</sub> & CO<sub>2</sub>)
- Colorimetry Tools
- Force & Pressure
- Cardiac Output (via bioimpedance), Stroke Volume
- Auditory, Somatosensory & Visual Evoked Response
- Angle of Movement
- Acceleration

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs

A

10


www.biopac.com

## Muscular

Record EMG data to investigate the properties of skeletal muscle. Record and display raw and integrated EMG signals; overlap the signals for better correlation of the data.

Measure strength and repeat trials for motor unit recruitment, summation and fatigue.

Use the reflex hammer transducer to study reflex response (neural control), or add the dynamometer to study handgrip strength profiles. Use the new Finger Twitch transducer for threshold, summation, tetanus and fatigue analysis. Students can listen to the muscle activity through headphones and note the increase in sound intensity as grip strength is increased through motor unit recruitment.

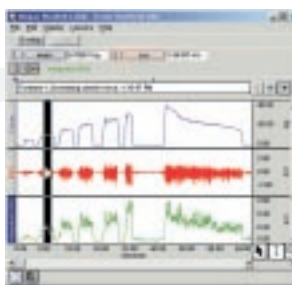

EMG and force

## Nerve Conduction

Combine the human-safe stimulating electrode with the stimulator to record nerve conduction experiments. Stimulate the ulnar nerve at three different points and record nerve conduction time. Measure the distance between the stimulation and recording points and then calculate velocity.

## Neurophysiology


Study EEG under a variety of conditions to explore relaxation and brain rhythms—software will filter and display each rhythm separately: Alpha, Beta, Delta, and Theta. Select from a variety of lessons to study Alpha rhythms in the occipital lobe, reaction times, and hemispheric asymmetry. Add the EOG (occipital signal) to study eye movement, saccades, tracking, angular displacement, or ocular fixations. Use the Stimulator to study evoked response—auditory, visual, or somatosensory. The new high-speed MP35 hardware allows for the recording of spontaneous nerve activity, continuously, at speeds of 100,000 samples a second. Add the SuperLab stimulus presentation package to perform psychophysiology stimulus/response protocols examining high-level brain activity. Conduct microelectrode recordings and study action potentials.

## Autonomic Nervous System

Record a wide variety of ANS-related signals such as EDA/GSR, temperature, ECG, pulse, respiration, airflow, nerve conduction, continuous blood pressure, and continuous non-invasive cardiac output (using bioimpedance technique) to demonstrate changes in the parasympathetic and sympathetic nervous system activity. Use the Polygraph lesson for simultaneous heart rate, electrodermal activity (GSR), and respiration rate. See Psychophysiology & Neurophysiology on page 18 for more details.

## Exercise Physiology & Biomechanics

Combine a wide range of respiratory system and pulmonary function signals with biomechanical data. Use the new Gas Analysis Module for online analysis of expired O<sub>2</sub>

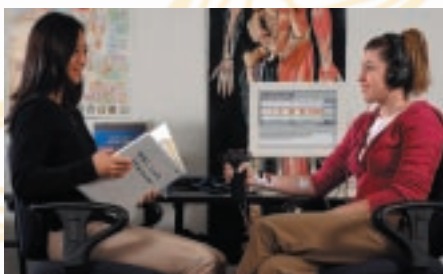

Dynamic lessons engage students

and CO<sub>2</sub> levels. Options are available for continuous blood pressure and cardiac output, and the system easily interfaces with force plates and other instrumentation (more than 18 ready-made connectors available). There are more than 27 specialized lessons—see page 16 for details.

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **human physiology**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

### Human Physiology Core

#### BSLHPY-W (Win) or BSLHPY-M (Mac)

Basic BSL System	BSLBSC, p. 8
Hand Dynamometer	SS25LA, p. 27
BP Cuff Transducer	SS19L, p. 27
Electronic Stethoscope Transducer	SS30L, p. 28
Airflow Transducer	SS11LA, p. 27
Airflow Filters (10/pk)	AFT1, p. 34
Airflow Mouthpieces (10/pk)	AFT2, p. 34
Airflow Nose Clips (10/pk)	AFT3, p. 34
Calibration Syringe (600 ml)	AFT6, p. 34

### Perform 25 or more lessons with this core package:

#### Muscular

BSL1	Standard & Integrated EMG
BSL2	Motor Unit Recruitment & Fatigue
H07	EMG Contractions - Active Learning
H27	Facial EMG
H34	EKG Electrogastrogram
H36	Muscular Biofeedback

#### Cardiovascular

BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL16	Blood Pressure & Korotkoff Sounds
BSL17	Heart Sounds & Cardiac Events
H08	Dive Reflex - Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

#### Pulmonary Function

BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV

#### Neurophysiology

BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL10	Eye Movement, Saccades & Fixation
H10	EEG & Hemispheric Asymmetry
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image

#### Biomedical Engineering

H20	BME Filtering
H33	FFT Fast Fourier Transform

See page 43-45 for a description of all available lessons.

### Increase your lab options with...

Stimulator	BSLSTM, p. 25
Stim. Electrode for humans	HSTM01, p. 28
Finger Twitch Transducer	SS61L, p. 30
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GASSYS2-EA, p. 34
Dissolved O <sub>2</sub> Probe	RXPROBE02, p. 31
pH Probe	RXPROBE01, p. 31
Reflex Hammer Transducer	SS36L, p. 30
Multi-Lead ECG Cable	SS29L, p. 27
Cardiac Output Sensor	SS31L, p. 29
Goniometer	SS21L, p. 29
Respiratory Effort Trans.	SS5LB, p. 26
Temperature Trans.	SS6L, p. 26
Headphones	OUT1, p. 26
EDA (GSR) Lead	SS57L, p. 26
SuperLab System	STP35W, p. 32
Transducer Accessory Pack	BSLHPY-TA, p. 42




# animal physiology & intro human phys

The Biopac Student Lab provides a wide range of options for animal and tissue experiments. Lessons allow students to study animal and human species for comparative physiology programs. Use the new Gas Analysis Module for human and animal experiments for the analysis of expired  $O_2$  and  $CO_2$ . The new Dissolved Oxygen probe allows students to monitor the oxygen consumption of a goldfish. Perform intracellular recording and membrane transport studies.

■ See Human Physiology on pages 10-11 for more details.

## Cardiovascular Hemodynamics

Lessons cover ECG (1-12 lead) and allow students to isolate components of the ECG complex and use the averaging features in the software for further ECG analysis. Use the electronic stethoscope to examine heart sounds and then overlap the data to correlate the sounds with the mechanical and electrical events of the cardiac cycle. Record arterial blood pressure and record systolic, diastolic, mean,  $dp/dt$  max and min, and use the noninvasive Cardiac Output Sensor to record stroke volume and cardiac output. If flow signals are available, record them simultaneously with pressure to examine vascular resistance and compliance. Plot pressure versus flow to obtain P/V Loops. See the Frog Heart and Turtle Heart lessons on page 45 for refractory heartblock and vagal escape.

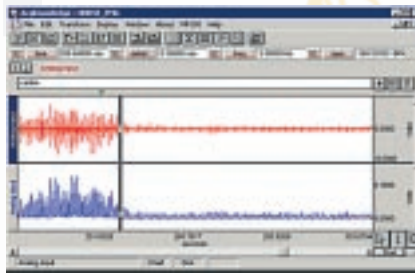

EMG from a horse


## Muscular

Record EMG to investigate the properties of skeletal muscle. Record and display raw and integrated EMG signals; overlap the signals for better correlation of the data. Measure strength and repeat trials for motor unit recruit-


ment/summation and fatigue. Demonstrate the treppe (staircase) phenomenon. Study the contractility of skeletal muscle with the force transducer and stimulator. The Frog Gastrocnemius lesson records threshold, maximal response, summation, tetanus, and fatigue.

## Compound Action Potential

Use the new Nerve Chamber (or an existing chamber) with the Low Voltage Stimulator (SS58L, page 25) to record the compound action potential and nerve conduction from the frog sciatic nerve. Record action potentials from

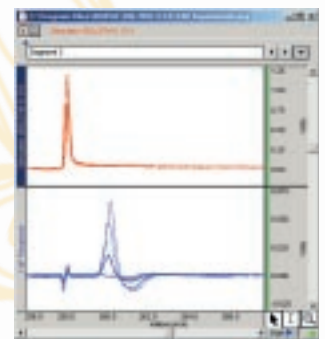

CAP in Overlap Mode

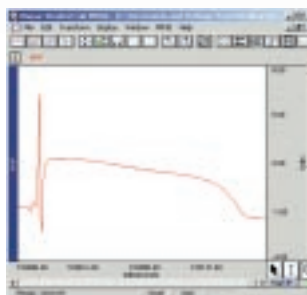

MAP—Guinea Pig

cockroaches, crawfish and earthworms. Add a range of drugs and determine the effect they have on the nerve. The new Nerve Chamber includes a drug delivery chamber (agent well) with lid to maximize the quality of the results and improve experimental repeatability.

## features

- 32 lessons targeted for Animal & Intro. Human Phys.
- ECG, EEG, EMG, EOG & EGG
- Temperature
- Gas Analysis  $CO_2$  &  $O_2$
- $pO_2$
- pH
- Bioimpedance & Cardiac Output
- Force
- Neurophysiology
- Hemodynamics
- Respiratory & Pulmonary Function
- Reaction Time
- Nerve Recordings & Compound Action Potentials
- Membrane Transport (drug delivery)
- Tissue Baths
- Stimulation & Response
- Isolated Heart, Lung, Muscle
- Auditory, Somatosensory & Visual Evoked Response
- Compatible with Crawdad Lab Manual

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs

A

12


www.biopac.com

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **animal physiology**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.


See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

## Neurophysiology

Study EEG under a variety of conditions to explore relaxation and brain rhythms—the software can filter and display each rhythm separately: Alpha, Beta, Delta and Theta. Select from a variety of *PRO* Lessons to study Alpha rhythms in the occipital lobe, reaction times, and hemispheric asymmetry. Add the EOG (ocular signal) to study eye movement, saccades, tracking, angular displacement, or ocular fixations. Use the stimulator to study evoked response — auditory, visual, or somatosensory. The new high-speed MP35 hardware allows the recording of spontaneous nerve activity, continuously, at speeds of 100,000 samples a second. Add the SuperLab stimulus presentation package to perform stimulus response studies investigating higher order neuronal function in humans. Conduct intracellular and extracellular recordings with glass microelectrodes and Ag-AgCl wire electrodes to study action potentials from a variety of subjects.

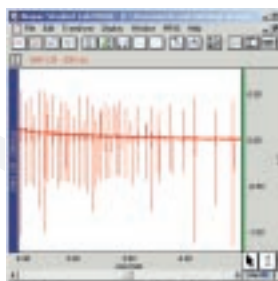

Cockroach nerve

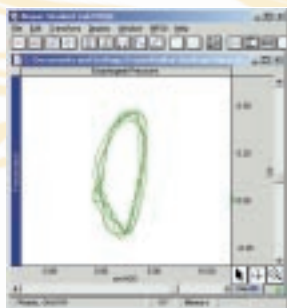

Flow vs. pressure in X/Y mode

## Respiratory & Pulmonary Function

The BSL System provides an excellent introduction to respiratory system and pulmonary function volumes and capacities, respiratory rates, breathing, and ventilation. Display and/or print a clinical grid on the data for effective interpretation and training. A complete range of airflow and pressure transducers is suitable for small, medium and large animals as well as humans.

## Gas Analysis

Use the new gas analysis system for detailed metabolic studies in small and medium sized animals. Chambers, couplers, facemasks, and tubing options provide setup options for any protocol. The system can provide online measures of O<sub>2</sub> and CO<sub>2</sub> for RER, VO<sub>2</sub> and BMR/RMR. Use the Dissolved O<sub>2</sub> probe and lesson to measure the oxygen consumption of a goldfish.

## Intracellular & Membrane Transport

Use the Voltage Drive & Current Monitor Cable (BSLCBL10, page 38) to perform ion transport studies. The cable uses the output from the BSL hardware to drive a set of stimulation electrodes and also monitors stimulation current. Use the new High-Impedance Cable (BSLCBL8/9, page 36) to record from the cockroach ventral nerve and for a variety of intracellular and extracellular recordings.

## In vitro Applications

The new Tissue Bath Stations provide students with research-quality equipment in a modular, flexible configuration. The Visceral Smooth Muscle lesson guides students through the entire recording and analysis process. Students can also electrically stimulate tissue preparations, including field stimulation, with the BSL Stimulator. Interface with Ussing chambers for ion transport studies. Record and analyze data from isolated heart and lung experiments.

## Animal Physiology Core

### BSLAPH-W (Win) or BSLAPH-M (Mac)

Basic BSL System (with BSLCBL8 substituted for SS2L)	BSLBSC, p. 8
Dissolved O <sub>2</sub> Probe Interface	BSL-TCI16, p. 39
Force Transducer (200 g)	SS65L, p. 31
Lead (unshielded) x 2	LEAD110, p. 36
Leads (shielded) x 2	LEAD110S-W/R, p. 36
Low Voltage Stimulator	SS58L, p. 25
Needle Electrodes x 3	EL452, p. 38
Nerve Chamber	NERVE2, p. 38
Recording Nerve Cable	BSLCBL4B, p. 39
Stim. Electrodes for animals	ELSTM2, p. 38
Stimulator Nerve Cable	BSLCBL2A, p. 39

## Perform 32 or more lessons with this core package:

### Muscular

A02	Frog Gastrocnemius
A05	Visceral Smooth Muscle
A11	Resting Potential from Crawdad Manual
A12	Membrane Potential (muscle)
A15	Earthworm Smooth Muscle
BSL1	Standard & Integrated EMG
H07	EMG Contractions - Active Learning
H27	Facial EMG
H34	EKG Electrogastrogram

### Cardiovascular

A04	Frog Heart
A09	Turtle Heart
BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
H08	Dive Reflex - Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

### Pulmonary Function

A07	Dissolved O <sub>2</sub> (goldfish)—with your probe
-----	---

### Neurophysiology

A01	Frog Pith & Prep
A03	Frog Nerve
A06	Cockroach Nerve
A08	Action Potential
A14	CPG Hornworm
BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL10	Eye Movement, Saccades & Fixation
H10	EEG & Hemispheric Asymmetry
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image

### Biomedical Engineering

H20	BME Filtering
H33	FFT Fast Fourier Transform


See page 43-45 for a description of all available lessons.

## Increase your lab options with...

Cardiac Output Sensor	SS31L, p. 29
Dissolved O <sub>2</sub> Probe	RXPROBE02, p. 31
pH Probe	RXPROBE01, p. 31
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GASSYS2-EA, p. 34
Tissue Bath Station	ITBS100, p. 33
Temperature Transducer	SS6L, p. 26
Transducer Accessory Pack	BSLAPH-TA, p. 42



# biology

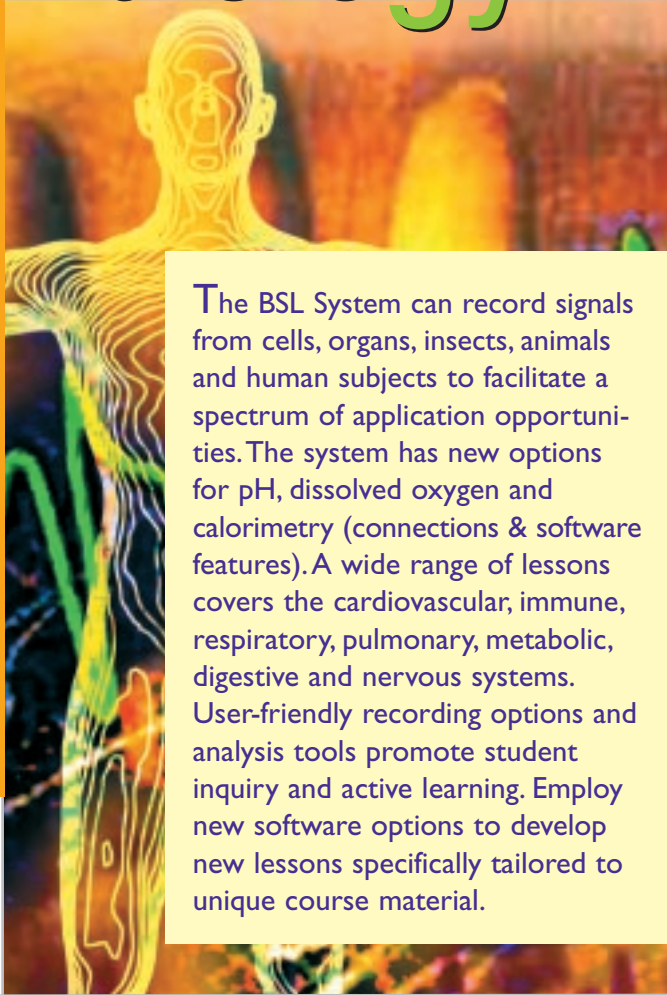

The BSL System can record signals from cells, organs, insects, animals and human subjects to facilitate a spectrum of application opportunities. The system has new options for pH, dissolved oxygen and calorimetry (connections & software features). A wide range of lessons covers the cardiovascular, immune, respiratory, pulmonary, metabolic, digestive and nervous systems. User-friendly recording options and analysis tools promote student inquiry and active learning. Employ new software options to develop new lessons specifically tailored to unique course material.

A

## features

- 53+ lessons targeted for Biology
- ECG, EEG, EMG, EOG & EGG
- Respiratory & Pulmonary Function
- Temperature
- pH
- Dissolved O<sub>2</sub>
- Gas Analysis (O<sub>2</sub> & CO<sub>2</sub>)
- Oxygen uptake
- Blood Pressure
- Cardiac Output (via bioimpedance)
- Stroke Volume
- Tissue Baths
- Force
- Calorimeter Interface
- Pulse
- Autonomic Nervous System
- Nerve Conduction Velocity
- Colorimetry Tools

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs

Biology covers a wide spectrum of applications. To complement those outlined below, see the applications detailed for Human Physiology (page 10) and Animal Physiology (page 12). As with all disciplines, the selected applications only begin to suggest what you can do with the BSL System. Contact a Biopac Student Lab Specialist to discuss your application needs.

### Cardiovascular


Lessons include blood pressure, ECG analysis, heart sounds, and pulse. Students can make single-, three-, six- and 12-lead ECG recordings, familiarizing themselves with Einthoven's triangle and mean electrical axis of the frontal plain. Students perform blood pressure measurements using a cuff, with stethoscope, employing Korotkoff sounds to make determinations of systolic/diastolic pressure. Students can also utilize the latest technology for continuous non-invasive BP recordings. Use the noninvasive Cardiac Output Sensor (SS31L, page 29) on human and animal subjects to record stroke volume and cardiac output. By measuring continuous mean arterial pressure simultaneous with flow (cardiac output), students can examine vascular resistance.



ECG and pulse


### Respiratory & Pulmonary Function

Students can record the pattern of breathing as well as perform comprehensive cardiopulmonary tests. There are lessons for determining tidal volumes and lung capacities, including FEV<sub>1,2,3</sub> and MVV, plus full gas analysis studies.

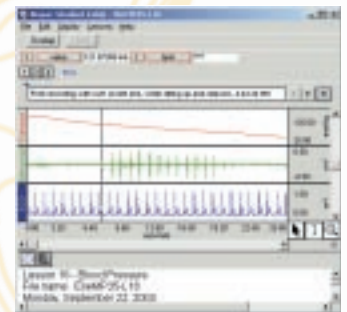

Blood pressure from Lesson 16

### Gas Analysis

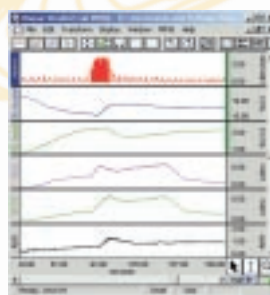

Respiratory exchange ratio


Students can use the new gas analysis system to monitor expired CO<sub>2</sub> and O<sub>2</sub> levels for detailed metabolic studies with human and animal subjects. Take measures to study ventilation and heat exchange. Use the Dissolved O<sub>2</sub> probe and lesson to measure the oxygen consumption of a goldfish. Chambers, facemasks, and tubing accessories provide setup options for any protocol.

### Neurophysiology

Study EEG under a variety of conditions to explore relaxation and brain rhythms—the software can filter and display each rhythm separately: Alpha, Beta, Delta, and Theta. Study Alpha rhythms in the occipital lobe, reaction times, and hemispheric asymmetry. Use the system to demonstrate changes in parasympathetic and sympathetic tone. Add EOG (ocular signal) to study eye movement, saccades, tracking, angular displacement, or ocular fixations. Use the stimulator to study evoked response—auditory, visual, or somatosensory. Record reaction time and measure nerve conduction velocity. The system samples quickly enough to easily record spontaneous nerve activity from small animals and insects.

## Digestive System


Use the system to record gastric signals in humans and animals or monitor isolated intestinal (gut) strips in a tissue bath experiment. Monitor the EGG (Electrogastrogram) before and after food digestion. Examine gastric slow wave propagation, peristaltic (slow wave) propagation, and gastrointestinal motility.

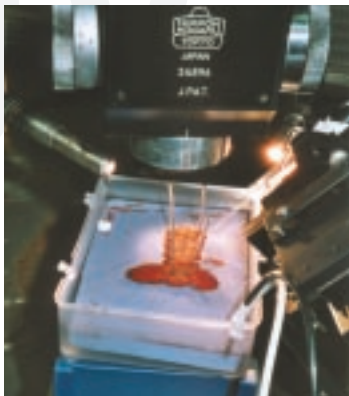

Membrane potential using crawfish and glass microelectrodes

## Cellular Biology

Use the BSL hardware with suitable glass microelectrodes to demonstrate membrane potential with the frog sartorius muscle or the crawfish, and active transport using frog skin. These experiments demonstrate the principle of homeostasis and serve as a good introduction to electrophysiology recording techniques. The system will also work with the popular *Crawdadd* CD-ROM Lab Manual for Neurophysiology by Wytenbach, Johnson, and Hoy (ISBN 0-87893-947-4).

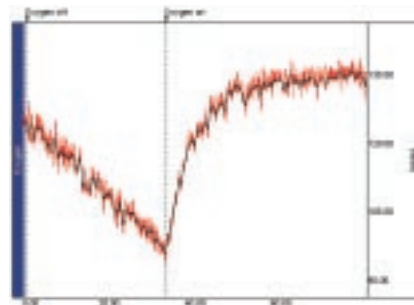

Dissolved oxygen (goldfish)

## Comparative Biology

The Biopac Student Lab system is an excellent tool for comparative biology programs because students can compare data from their own bodies with data recorded from a variety of animals. The curriculum covers most of the major physiological systems, including: brain, muscle, pulmonary, cardiovascular and CNS.

# biology

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **biology**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

### Biology Core

#### BSLBIO-W (Win) or BSLBIO-M (Mac)

Basic BSL System	BSLBSC, p. 8
Airflow Filters (10/pk)	AFT1, p. 34
Airflow Mouthpieces (10/pk)	AFT2, p. 34
Airflow Nose Clips (10/pk)	AFT3, p. 34
Airflow Transducer	SS11LA, p. 27
BP Cuff Transducer	SS19L, p. 27
Calibration Syringe (600 ml)	AFT6, p. 34
Colored Paper (for BSL9)	PAPER1
Dissolved O <sub>2</sub> Probe Interface	BSL-TCI16, p. 39
EDA (GSR) Lead	SS57L, p. 26

EDA (Isotonic) Electrodes (100/pk)	EL507, p. 37
Electrode Gel	GEL1, p. 36
Electronic Stethoscope Transducer	SS30L, p. 28
Force Transducer	SS12LA, p. 27
Hand Dynamometer	SS25LA, p. 27
Hand Switch	SS10L, p. 26
Headphones	OUT1, p. 26
High-Impedance Cable	BSLCBL9, p. 36
Multi-Lead ECG Cable	SS29L, p. 27
Needle Electrodes x 3	EL452, p. 38
Nerve Chamber	NERVE2, p. 38

pH Probe Interface	BSL-TCI21, p. 39
Pressure Transducer	SS13L, p. 27
Pulse Transducer	SS4LA, p. 26
Recording Nerve Cable	BSLCBL4B, p. 39
Respiratory Effort Transducer	SS5LB, p. 26
Stimulator	BSLSTM, p. 25
Stim. Electrode for humans	HSTM01, p. 28
Stim. Electrode for animals	ELSTM2, p. 38
Stimulator Nerve Cable	BSLCBL2A, p. 39
Surgical Tape	TAPE1, p. 36
Temperature Transducer	SS6L, p. 26

### Perform 53 or more lessons with this core package:

#### Muscular

A02	Frog Gastrocnemius
A05	Visceral Smooth Muscle
A11	Resting Potential from <i>Crawdadd</i> Manual
A12	Membrane Potential (muscle)
A15	Earthworm Smooth Muscle
BSL1	Standard & Integrated EMG
BSL2	Motor Unit Recruitment & Fatigue
H06	Finger Twitch
H07	EMG Contractions - Active Learning
H27	Facial EMG
H34	EGG Electrogastrogram
H36	Muscular Biofeedback

H08	Dive Reflex - Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

#### Pulmonary Function

A07	Dissolved O <sub>2</sub> (goldfish)—with your probe
BSL8	Respiratory Cycle
BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV
BSL15	Aerobic Exercise Physiology

#### Neurophysiology

A01	Frog Pith & Prep
A03	Frog Nerve
A06	Cockroach Nerve
A08	Action Potential-Earthworm
A09	Turtle Heart
A14	CPG Hornworm
BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL9	GSR and Polygraph
BSL10	EOG Eye Movement, Saccades & Fixation
BSL11	Reaction Time
BSL14	Biofeedback: Relaxation & Arousal

H03	Nerve Conduction (ulnar nerve)
H10	EEG & Hemispheric Asymmetry
H11	Mirror Test—EDA Sensory motor learning
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image
H16	Reflexes & Reaction Time - Active Learning
H24	Habituation

#### Biomedical Engineering: H02, H20 & H33

See page 43-45 for a description of all available lessons.

#### Cardiovascular

A04	Frog Heart
BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
BSL16	Blood Pressure & Korotkoff Sounds
BSL17	Heart Sounds & Cardiac Events
H01	12-lead ECG
H04	BP Response to Straining
H05	WANT Wingate Test

#### Increase your lab options with...

Cardiac Output Sensor	SS31L, p. 29
Dissolved O <sub>2</sub> Probe	RXPROBE02, p. 31
Finger Twitch Transducer	SS61L, p. 30
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GASSYS2-EA, p. 34
pH Probe	RXPROBE01, p. 31
Reflex Hammer Transducer	SS36L, p. 30




# exercise physiology & biomechanics

With the Biopac Student Lab, it's simple to combine a wide range of respiratory system & pulmonary function signals with biomechanical data. Use the new Gas Analysis Module for online analysis of expired  $O_2$  and  $CO_2$  levels. Simultaneously record continuous, noninvasive blood pressure and cardiac output (bioimpedance method). BIOPAC offers transducers for angle of limb movement, acceleration, heel-toe strike, etc. for gait analysis, range of motion and other related studies. The system easily interfaces with force plates, motion analysis equipment and other instrumentation.

## Respiratory & Pulmonary Function


The BSL includes a range of experiments for pulmonary function and Gas Analysis studies. Detailed lessons guide students through the classical pulmonary function tests, including tidal volume, inspiratory capacity, expiratory capacity, functional residual capacity, vital capacity, total lung capacity, forced vital capacity, forced expiratory volume ( $FEV_{1, 2, 3}$ ) and Maximal Voluntary Ventilation (MVV).

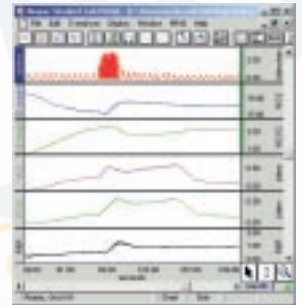

Respiratory exchange ratio

## Gas Analysis

Use the new Gas Analysis Module to create a powerful metabolic analyzer (cart) for students to measure expired  $O_2$  and  $CO_2$  and use lessons for  $VO_2$ , Respiratory Exchange Ratio, and Basal or Resting Metabolic Rate. Combine metabolic and cardiac output measurements to give students a detailed view of a variety of cardiopulmonary responses. If your protocol examines  $CO_2$  above 5% or requires high-speed response, contact BIOPAC to discuss the full range of gas analyzers.

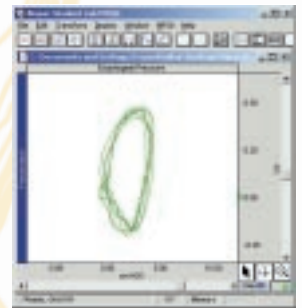


Airflow Pressure/Volume loop

## features

- 27 lessons targeted for Exercise Phys. & Biomechanics
- ECG, EEG, EMG, EOG & EGG
- Respiration
- Temperature
- Airflow & Lung Volume
- Gas Analysis ( $CO_2$  &  $O_2$ ) — Metabolic Cart
- Cardiac Output (via bioimpedance)
- Respiratory Exchange Ratio
- Basal or Resting Metabolic Rate
- Motor Unit Recruitment
- 12-Lead ECG
- Heart Sounds
- Blood Pressure
- Stroke Volume
- Gait Analysis (including Heel-Toe Strike)
- Range of Motion
- Acceleration, Velocity, Distance

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs

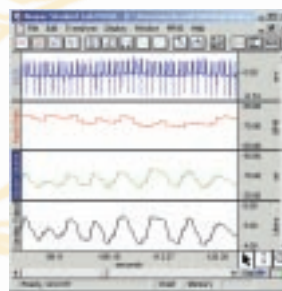

Cardiac output data

## Cardiovascular

The BSL System includes many ECG lessons, including 12-lead ECG, cardiac output, and continuous blood pressure lessons. A Heart Sounds lesson allows students to listen to and record heart sounds, while comparing them to the ECG complex. There are also lessons for the Wingate test and Blood Pressure Response to Isometric Straining Exercise. Record cardiac output from human subjects—connect the bioimpedance electrodes to a subject and record stroke volume and cardiac output in real time pre- and post-exercise. There is also an option for continuous blood pressure monitoring, during exercise.


## Biomechanics

Students can use the new Reflex Hammer in conjunction with EMG recordings to compare responses and reflex actions. For advanced studies,

I purchased the MP30 Advanced System and have found the student labs helpful in explaining and demonstrating physiological concepts and current technology. Personally, I have been using the *PRO* software for my own research purposes. I have instrumented an old cybex isokinetic ergometer with a force transducer to measure force and a potentiometer to measure joint angle. Data acquisition has been going very smoothly. I have been quite impressed with the *PRO* software and its abilities. I have set up a calculation channel that accounts for gravity's affect on limb weight at all joint angles throughout a range of motion. This allows me to accurately estimate muscular resistance. Our area rep took the time to come down to see us and has been helpful with questions.

—Dain LaRoche, Johnson State College, Environmental and Health Sciences

add angle of limb movement as a subject performs a variety of tasks. The system has transducers for recording heel and toe strike and acceleration for gait analysis studies. Use the Hand Dynamometer and Tri-Axial Accelerometer to measure isotonic and isometric performance. Goniometers are available for evaluating one or two degrees of freedom from the same joint (e.g. wrist flexion/extension and radial/ulnar deviations). Record Sit & Reach tests and analyze range of motion. Use the *XY* display mode to monitor motion resulting from two degrees of freedom. The BSL software will determine velocity of motion and calculate acceleration.

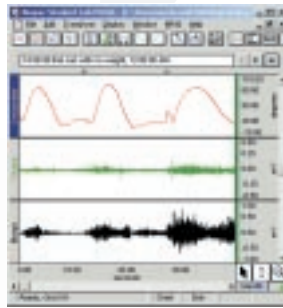

Goniometry data

## Nerve Conduction

Combine the human-safe stimulating electrode with the stimulator to record nerve conduction experiments. Stimulate the ulnar nerve and record nerve conduction time (Lesson H03). The procedure allows students to stimulate a subject at three different points along the ulnar nerve. The distance between the stimulation and recording points is measured and the velocity of signal propagation along the nerve is calculated.

## Muscular

Students can measure EMG and Integrated EMG, including force and angle of limb movement. Students can listen to the sound of the electrical activity coming from muscle as they squeeze a hand dynamometer and equate changes in sound with changes in force. Add the BSL Stimulator and human-safe stimulation electrode for a variety of muscle stimulation experiments. Include visual and auditory feedback with touch for Muscular Biofeedback studies.



Ergometer modified to record WAnT using BIOPAC

Photo courtesy A. Zidermanis, PhD, formerly at Parker College of Chiropractic

## Interface with Existing Equipment

The BSL System offers over 60 industry-standard transducers. Further, the system easily interfaces with other major amplifier and transducer manufacturers encompassing the most commonly used exercise physiology lab products such as force plates, ergometers, motion analysis systems, and metabolic carts. The BSL System can trigger or receive trigger information from other equipment. For interfacing, choose from a variety of ready-made connectors and

cables, or use the custom connector kit (page 39). BIOPAC support staff can help determine the appropriate interface for any complete system.

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **exercise physiology & biomechanics**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

## Exercise Physiology Core

### BSLEXP-W (Win) or BSLEXP-M (Mac)

Basic BSL System	BSLBSC, p. 8
Airflow Filters (10/pk)	AFT1, p. 34
Airflow Mouthpieces (10/pk)	AFT2, p. 34
Airflow Nose Clips (10/pk)	AFT3, p. 34
Airflow Transducer	SS11LA, p. 27
BP Cuff Transducer	SS19L, p. 27
Calibration Syringe (600 ml)	AFT6, p. 34
Electronic Stethoscope Transducer	SS30L, p. 28
Hand Dynamometer	SS25LA, p. 27
Hand Switch	SS10L, p. 26
Headphones	OUT1, p. 26
Pulse Transducer	SS4LA, p. 26
Respiratory Effort Transducer	SS5LB, p. 26
Temperature Transducer	SS6L, p. 26
Surgical Tape	TAPE1, p. 36

## Perform 27 or more lessons with this core package:

### Muscular

BSL1	Standard & Integrated EMG
BSL2	Motor Unit Recruitment & Fatigue
H07	EMG Contractions - Active Learning
H27	Facial EMG
H34	EKG Electrocardiogram
H36	Muscular Biofeedback

### Cardiovascular

BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
BSL16	Blood Pressure & Korotkoff Sounds
BSL17	Heart Sounds & Cardiac Events
H05	WAnT Wingate Test
H08	Dive Reflex - Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

### Pulmonary Function

BSL8	Respiratory Cycle
BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV
BSL15	Aerobic Exercise Physiology

### Neurophysiology

BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL11	Reaction Time
H10	EEG & Hemispheric Asymmetry
H16	Reflexes & Reaction Time - Active Learning

### Biomedical Engineering

H02	Compartmental Modeling
H20	BME Filtering
H33	FFT Fast Fourier Transform

See page 43-45 for a description of all available lessons.

## Increase your lab options with...

Airflow Transducer for Ex. Phys.	SS52L, p. 33
Calibration Syringe (2 L)	AFT26, p. 35
Cardiac Output Sensor	SS31L, p. 29
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GAS-SYSTEM2-EA, p. 34
Finger Twitch Transducer	SS61L, p. 30
Goniometer	SS21L, p. 29
Reflex Hammer	SS36L, p. 30
Stimulator	BSLSTM, p. 25
Transducer Accessory Pack	BSLEXP-TA, p. 42




# psychophysiology & neurophysiology

Present a wide array of psychophysiology experimental techniques with the BSL System. Lessons guide students through recording and analysis modalities to provide building blocks that empower students to perform increasingly advanced studies. Acquire signals for ECG, EDA (GSR), EEG, EMG, EOG and EGG. Combine physiological data with trigger information time-synched with advanced stimulus presentation paradigms. Record noninvasive cardiac output and blood pressure while students perform tasks or respond to a presentation.

## Autonomic Nervous System

Use the BSL system to record changes in the parasympathetic and sympathetic nervous system activity. Acquire a wide variety of ANS-related signals such as EDA (GSR), skin temperature, ECG, pulse, respiration, airflow, nerve conduction, continuous blood pressure, and continuous noninvasive cardiac output (bioimpedance method). Simultaneous graphing shows heart rate, electrodermal activity (GSR) and respiration rates. After the experiment, evaluate the data using the powerful and user-friendly analysis functions.

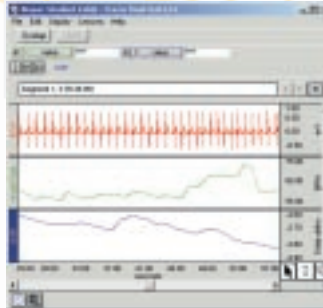

Electrodermal activity (GSR) from Lesson 14

## Event Related Potentials

Combine the BSL with a visual presentation system such as SuperLab to explore a wide variety of experiments like Stroop, Oddball, Habituation, and Startle Response. SuperLab is user-friendly and greatly simplifies development of a range of visual and auditory presentations. As each stimulus is presented, SuperLab sends a digital pulse to the BSL System to mark the onset of the stimulus. The stimuli are classified into different groups with each group assigned a unique digital channel. The BSL software will identify the digital pulse and automatically measure the associated response, and provide the average or perform measurements on each response in the stimulus classification.



Facial EMG electrodes

## Evoked Response

Lessons guide students through evoked response recordings. Use the system to trigger a stimulus and derive the average response. Use with headphones, stroboscope, or electrical or mechanical stimulation for auditory, visual or somatosensory response.



EEG filtered from Lesson 3

## features


- 28 lessons targeted for Psychophys. & Neurophys.
- ECG, EEG, EMG, EOG & EGG
- Autonomic Nervous System
- Auditory & Visual Evoked Response
- Event Related Potential
- Startle Eye Blink Experiments
- Nerve Conduction
- Habituation
- Stroop
- Sensory Motor Learning
- Electrodermal Activity (GSR)
- Heart Rate Variability
- Visual Presentation System (SuperLab & E-Prime)
- Cardiac Output (via bioimpedance)
- Pre-ejection Period (PEP)
- Automatic Continuous Noninvasive Blood Pressure

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs.

Powerful set-up and analysis tools make the BSL ideal for graduate level studies and personal research.

## EEG

Introductory lessons allow students to record EEG and look at the differences between Alpha, Beta, Delta and Theta activity levels. Students can record EEG from both hemispheres while listening to music, reading and performing mental tasks. The software guides them through the recording and then shows them how to analyze the frequency components of the signal. The Fast Fourier Transformation is used to show the frequency levels of the activity on each hemisphere.

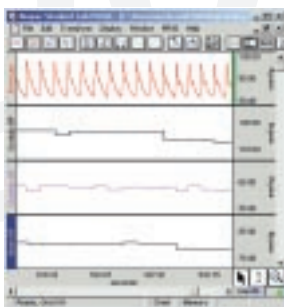


Noninvasive blood pressure data

## Cardiovascular

Students can record the changes in pre-ejection period and cardiac output relative to a variety of different stimuli. The new noninvasive Cardiac Output Sensor greatly simplifies previously complex procedures to allow students to record stroke volume and cardiac output. Combine these signals with ECG and blood pressure to demonstrate cardiovascular regulation by the Autonomic Nervous System. The system can also perform detailed Heart Rate Variability tests and Averaged ECG recordings.

## Muscular

Students can measure EMG and Integrated EMG, including force and angle of limb movement. They can listen to the sound of the electrical activity coming from the muscle as they squeeze a hand dynamometer and equate changes in sound form and intensity with changes in force. Add the BSL Stimulator and human-safe stimulation electrode for a variety of muscle stimulation experiments. Use the Tri-Axial Accelerometer to examine movement of limbs, head and torso in three-dimensional space. Couple the resultant acceleration, velocity and distance data with EMG recordings to obtain a comprehensive picture of skeletal muscle performance.

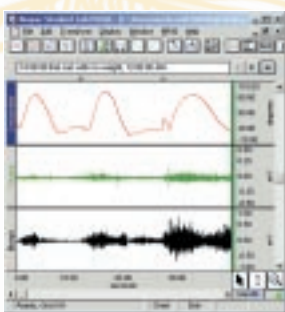

EMG data: Goniometer, Triceps, Biceps

## EOG

Students can record horizontal and vertical eye movements while observing fixation and tracking. The XY display mode will plot horizontal vs. vertical eye movement to track eye position relative to an image or object. By using the zoom and measurement tools, students can measure duration of saccades and fixation. A simple light fixture becomes an excellent tool for measuring angular displacement.



EDA mirror test (Lesson H11)

## Nerve Conduction

Combine the human-safe stimulating electrode with the stimulator to record nerve conduction experiments. Stimulate the ulnar nerve at three different points and record nerve conduction time. Measure the distance between the stimulation and recording points and then calculate the velocity of signal propagation along the nerve (motor response). Use the Finger Twitch transducer coupled with nerve stimulation to examine the relationship between applied (external) nerve stimulation and associated motor recruitment.

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **psychophysiology & neurophysiology**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

### Psychophysiology Core

#### BSLPSY-W (Win) or BSLPSY-M (Mac)

Basic BSL System	BSLBSC, p. 8
BP Cuff Transducer	SS19L, p. 27
Colored Paper (for BSL9)	PAPER1
EDA (GSR) Lead	SS57L, p. 26
EDA (Isotonic) Electrodes (100/pk)	EL507, p. 37
Electronic Stethoscope Transducer	SS30L, p. 28
Hand Switch	SS10L, p. 26
Headphones	OUT1, p. 26
Pulse Transducer	SS4LA, p. 26
Respiratory Effort Transducer	SS5LB, p. 26

### Perform 28 or more lessons with this core package:

#### Muscular

BSL1	Standard & Integrated EMG
H07	EMG Contractions - Active Learning
H27	Facial EMG
H34	EGG Electrogastrogram
H36	Muscular Biofeedback

#### Cardiovascular

BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
BSL16	Blood Pressure & Korotkoff Sounds
BSL17	Heart Sounds & Cardiac Events
H08	Dive Reflex - Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

#### Neurophysiology

BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL9	GSR and Polygraph
BSL10	EOG Eye Movement, Saccades & Fixation
BSL11	Reaction Time
BSL14	Biofeedback: Relaxation & Arousal
H10	EEG & Hemispheric Asymmetry
H11	Mirror Test—EDA Sensory motor learning
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image
H16	Reflexes & Reaction Time - Active Learning
H24	Habituation

#### Biomedical Engineering

H02	Compartmental Modeling
H20	BME Filtering
H33	FFT Fast Fourier Transform


See page 43-45 for a description of all available lessons.

### Increase your lab options with...

Cardiac Output Sensor	SS31L, p. 29
Noninvasive Blood Pressure Monitor	NIBP100A, p. 32
Psych. Variable Assessment Trans.	SS43L, p. 30
Stimulator	BSLSTM, p. 25
SuperLab Package	STP35, p. 32
Finger Twitch Transducer	SS61L, p. 30
Stim. Electrode for Humans	HSTM01, p. 32
Transducer Accessory Pack	BSLPSY-TA, p. 42



# biomed engineering

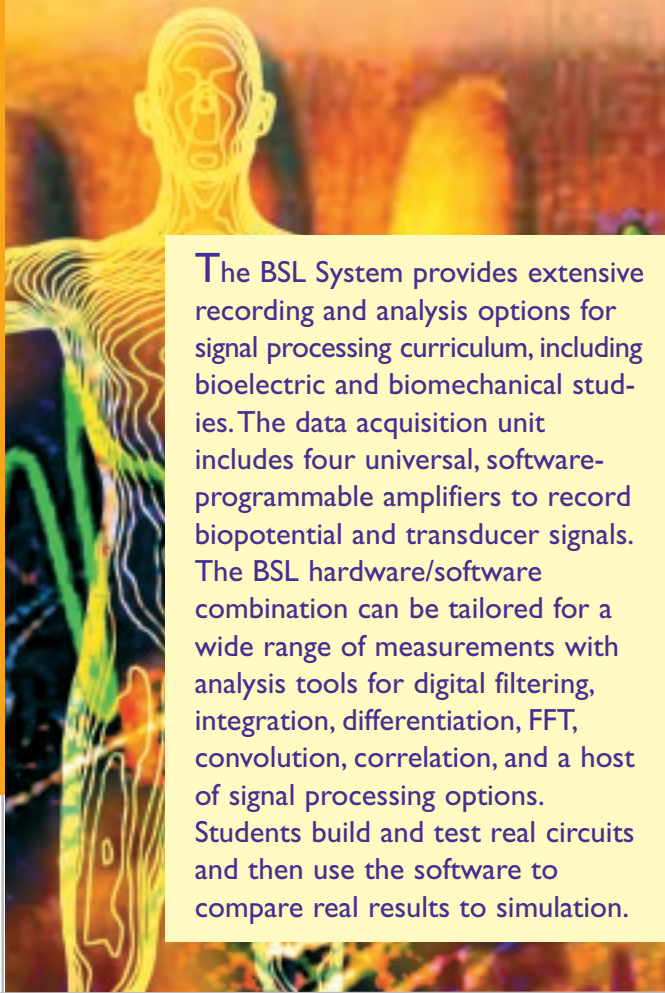


The BSL System provides extensive recording and analysis options for signal processing curriculum, including bioelectric and biomechanical studies. The data acquisition unit includes four universal, software-programmable amplifiers to record biopotential and transducer signals. The BSL hardware/software combination can be tailored for a wide range of measurements with analysis tools for digital filtering, integration, differentiation, FFT, convolution, correlation, and a host of signal processing options. Students build and test real circuits and then use the software to compare real results to simulation.

*“During each [BIOPAC] laboratory exercise, students follow detailed procedures that are designed to demonstrate principles of biophysics and biological measurement. Student feedback on the laboratories has been phenomenal, with the hands-on experiences motivating the students in a way that no lecture-only course can.” — IEEE Engineering in Medicine and Biology, July/August 2003 (Vol. 22, No. 4, pg. 106)*

## Signal Analysis & Processing


The BSL software has an extensive library of signal processing functions permitting graphical insight to analytical methods. The software can demonstrate the procedure and consequences associated with simple to complex signal processing methodologies. For example, students can view data before and after IIR or FIR filter processing, build a complex waveform from periodic signals (i.e., create a square wave from multiple sine waves) and decompose the result, or apply non-linear processing methods to data. Use the X/Y display mode to generate Lissajous patterns and investigate chaotic phenomena and demonstrate phase relationships between two variables. Signals can be correlated and convolved. Use the histogram function to focus on distribution of specific signal measures.

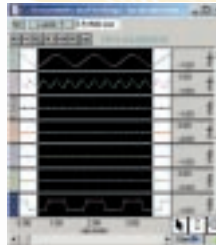

BME Square

## Transducers & Calibration


The BSL System employs a wide array of transducers that transform physical measures into electrical signals. The generic input design of the MP35 acquisition unit allows it to interface a huge variety of third-party or completely unique transducers. Students can use the BSL software to linearize and calibrate transducers and then compare results to expected values. Relate fundamental physical standards to more complicated measures. For example, calibrate the Airflow Transducer with a syringe, and then use the Airflow Transducer to calibrate a respiration sensor designed to monitor thoracic circumference.

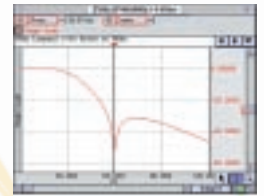

FFT of Biopotential filter frequency response

## Human & Animal Physiology

The wide range of human and animal physiology experiments provide a powerful tool for teaching students

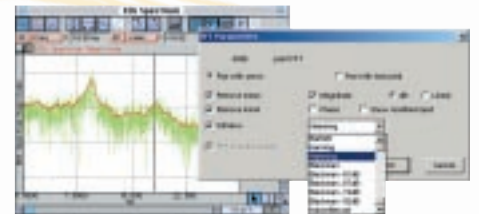

FFT dialog and EEG spectrum

the best technique and methodology for making a measurement. Each experiment demonstrates fundamental physiological concepts and educates students in the setup, recording and analysis process. Physiology basics are clearly explained. See pages 10-13 for details.

## features

- 60+ lessons targeted for Human & Animal Physiology
- Signal Analysis & Processing
- ECG, EDA (GSR), EEG, EGG, EMG & EOG
- Force, Pressure, Strain, Flow, Temperature, Sound, Light
- Filters (FIR & IIR)
- Instrumentation Design
- Respiratory System & Pulmonary Function
- Bioimpedance (Cardiac Output & Blood Flow)
- Biomechanics—Angle, Acceleration, Distance, Velocity
- Transducers & Calibration
- Physiological Control Systems
- Compartmental Modeling
- Blood Pressure & Heart Sounds
- Gait Analysis
- Chart, Overlap, Scope & X/Y Displays
- Spectral Analysis & Histograms
- Export to MatLab®, LabVIEW®, and MS Excel/Word®

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs




### Programming Options

Students can create their own programs to control the MP35 hardware with the BHAPI hardware application program interface. Students can also develop their own analysis programs to read the BIOPAC file format with the ACKAPI software application program interface. See page 24 for details.

### Physiological Control Systems & Compartmental Analysis


Implement simple experiments illustrating physiological control systems and compartmental analysis with the BSL System. The students can observe signal changes and then effect a change to observe a particular response. Investigate linear and nonlinear control paradigms. Create simple to intricate feedback loops where students perform a specific role in the loop operation. For instance, students can explore Westheimer's saccadic eye movement model—which represents the eye as a 2nd order system—then record eye motion via EOG set up, and then compare the real results to the modeled results to validate or adjust the model.

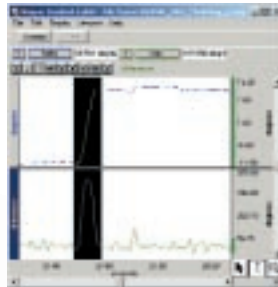

Compartmental Modeling (EOG)


### Biomechanics

The Student Lab System has a comprehensive ability to monitor gait and other mechanical responses. The system works with Goniometers, Accelerometers, Heel-Toe Strike transducers and Tri-Axial Accelerometers. Biopotential signals such as EMG can be synchronously recorded. Use the Hand Dynamometer and Tri-Axial Accelerometer to measure isotonic and isometric performance. Goniometers are available for evaluating one or two degrees of freedom from the same joint (e.g. wrist flexion/extension and radial/ulnar deviations). Use the XY display mode to monitor motion resulting from two degrees of freedom. Model mechanical systems, demonstrate principles of biomechanical resonance or inertial navigation (acceleration, velocity and position), or convert gravity vectors (from Tri-Axial Accelerometers) into associated “tilt” angles for use in ergonomic evaluations. See page 16 for details.

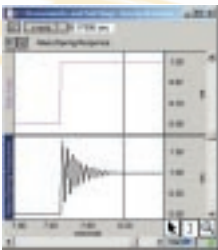

BME Step

### Instrumentation Design

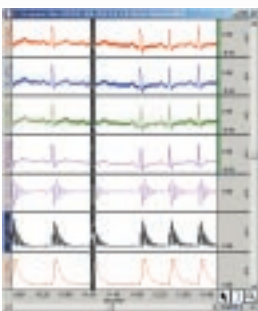

Circuit simulation

The new signal processing breadboard allows students to build and test real-world signal processing circuit modules and then verify their performance against mathematical simulation using graphical comparisons. Students can combine circuit modules, collect physiological signals and then analyze the results. Each circuit module constitutes an important subset of circuit design when recording and processing physiological signals. The BSL system is used like an oscilloscope to make measurements for circuit module evaluation.

### Interface with Existing Equipment

The BSL System offers over 60 industry-standard transducers. Further, the BSL System interfaces with other major amplifier and transducer manufacturers encompassing the most commonly used biomedical engineering instruments and sensors by using a wide variety of interface connectors and cables. Choose from 18 ready-made interface connectors, or build your own with the custom interface kit.

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **biomedical engineering**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23) for all available transducers, electrodes and accessories.

### Biomedical Engineering Core

#### BSLBME-W (Win) or BSLBME-M (Mac)

Basic BSL System	BSLBSC, p. 8
Airflow Filters (10/pk)	AFT1, p. 34
Airflow Mouthpieces (10/pk)	AFT2, p. 34
Airflow Nose Clips (10/pk)	AFT3, p. 34
Airflow Transducer	SS11LA, p. 27
BP Cuff Transducer	SS19L, p. 27
Calibration Syringe (600 ml)	AFT6, p. 34
Electronic Stethoscope Transducer	SS30L, p. 28
Hand Dynamometer	SS25LA, p. 27
Hand Switch	SS10L, p. 26
Headphones	OUT1, p. 26
Pulse Transducer	SS4LA, p. 26
Signal Processing Breadboard Lab	SS39L, p. 30

### Perform 29 or more lessons with this core package:

#### Muscular

BSL1	Standard & Integrated EMG
BSL2	Motor Unit Recruitment & Fatigue
H07	EMG Contractions—Active Learning
H27	Facial EMG
H34	EKG Electrogastragram
H36	Muscular Biofeedback

#### Cardiovascular

BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
BSL7	ECG & Pulse
BSL16	Blood Pressure & Korotkoff Sounds
BSL17	Heart Sounds & Cardiac Events
H08	Dive Reflex—Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

#### Pulmonary Function

BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV

#### Neurophysiology

BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL10	EOG Eye Movement, Saccades & Fixation
BSL11	Reaction Time
H10	EEG & Hemispheric Asymmetry
H12	EOG Saccades & Displacement
H13	EOG Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image
H16	Reflexes & Reaction Time - Active Learning

#### Bioengineering

H02	Compartmental Modeling
H20	BME Filtering
H25	BME Signal Processing (8 modules)
H26	ECG R-wave Detector
H33	FFT Fast Fourier Transformation

See page 42-44 for a description of all available lessons.

### Increase your lab options with...

Cardiac Output Sensor	SS31L, p. 29
Volt. Drive/Current Monitor Cable	BSLCBL10, p. 38
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GASSYS2-EA, p. 34
Finger Twitch Transducer	SS61L, p. 30
Heel/Toe Strike Transducer	SS28L, p. 29
Stimulator	BSLSTM, p. 25
Tri-Axial Accelerometer	SS26L, p. 29
Transducer Accessory Pack	BSLBME-TA, p. 42




# pharmacology & toxicology


The BSL System combines a range of experiment options for cardiovascular hemodynamics, respiratory system & pulmonary function, *in vitro* tissue and cellular studies. The new range of tissue bath stations provide an extra level of functionality for *in vitro* tissue experiments. Powerful, real-time analysis functions for pressure recordings include systolic, diastolic, mean BP, and dP/dt max and min. Similar tools are available for smooth muscle experiments, including peak, area, and derivative measurements. Add the noninvasive Cardiac Output Sensor to record stroke volume and cardiac output.

## Respiratory & Pulmonary Function

The BSL lessons include measurements of tidal volume, inspiratory capacity, expiratory capacity, functional residual capacity, vital capacity, total lung capacity, forced expiratory volume and maximal voluntary ventilation. The new Gas Analysis Module provides a powerful tool for metabolic studies. Measure expired O<sub>2</sub> and CO<sub>2</sub> with lessons for VO<sub>2</sub> max, respiratory exchange ratio, and basal or resting metabolic rate. Perform metabolic, respiratory and pulmonary measurements on a variety of species.

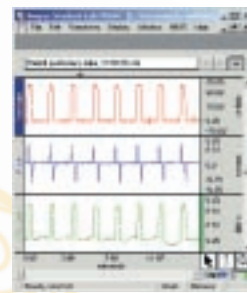

Rabbit pulmonary function

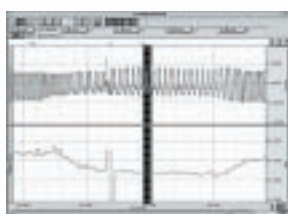

Frog heart rate response to drugs

## Cardiovascular Hemodynamics

Lessons cover ECG (1- through 12-lead), cardiac output, and continuous blood pressure. A heart sounds lesson allows students to listen to and record heart sounds, while comparing them to the ECG complex. The frog heart lesson explores cardiac rate and contractile response using a range of drug doses. Students can analyze blood pressure signals in real time.

## In vitro & Cellular Pharmacology

The new Tissue Bath Stations provide students with research-quality equipment in a modular, flexible configuration. The Visceral Smooth



Muscle lesson Drug effect and marker summary guides students through the entire recording and analysis process. Students can also electrically stimulate tissue preparations, including field stimulation, with the BSL Stimulator. Interface with Ussing chambers for ion transport studies. Record and also analyze data from isolated heart and lung experiments. Use the Nerve Chambers (page 38) for compound action potential studies. The system can record monophasic action potentials and spontaneous nerve activity.



New Tissue Bath

## features

- 32 lessons targeted for Pharmacology & Toxicology
- ECG, EEG, EGG, EOG & EMG
- Temperature
- Tissue Bath Station with integrated heating circulator
- Stimulator (direct or field)
- Cardiac Output (via bioimpedance)
- Dose Response Studies
  - Smooth Muscle
  - Cardiac Muscle
  - Skeletal Muscle
- Epithelial Transport
- Ion Transport/Ussing Chamber Measurements
- Gas Analysis Module
- Blood Pressure
- Isolated Heart/Lung
- Isolated Muscle
- Nerve Activity

Suitable for inquiry-based, active learning in 2-yr. & 4-yr. programs, medical schools, and nursing programs

A

22

www.biopac.com

The following hardware suggestions will enable you to perform a wide variety of applications targeted for **pharmacology & toxicology**. Use BIOPAC lessons or easily create your own experiments with the BSL *PRO* software included with each system. Order the core package or select items à la carte.

See **BSL Hardware** (page 23-40) for all available transducers, electrodes and accessories.

**Pharmacology & Toxicology Core**

**BSLPHA-W (Win) or BSLPHA-M (Mac)**

Basic BSL System (with BSLCBL8 substituted for SS2L)	BSLBSC, p. 8
Airflow Filters (10/pk)	AFT1, p. 34
Airflow Mouthpieces (10/pk)	AFT2, p. 34
Airflow Nose Clips (10/pk)	AFT3, p. 34
Airflow Transducer	SS11LA, p. 27
Pressure Transducer	SS13L, p. 27
Calibration Syringe (600 ml)	AFT6, p. 34
Force Transducer (200g)	SS65L, p. 31
Lead (unshielded) x 2	LEAD110, p. 36
Leads (shielded) x 2	LEAD110S-W/R, p. 36
Needle Electrodes x 3	EL452, p. 38
Nerve Chamber	NERVE2, p. 38
Recording Nerve Cable	BSLCBL4B, p. 39
Stimulator	BSLSTM, p. 25
Stimulator Nerve Cable	BSLCBL2A, p. 39
Stim. Electrodes for animals	ELSTM2, p. 38

**Perform 31 or more lessons with this core package:**

**Muscular**

A03	Frog Gastrocnemius
A05	Visceral Smooth Muscle
A11	Resting Potential from Crawdad Manual
A12	Membrane Potential (muscle)
A15	Earthworm Smooth Muscle
BSL1	Standard & Integrated EMG
H07	EMG Contractions—Active Learning
H27	Facial EMG
H34	EGG Electrogastrogram

**Cardiovascular**

A04	Frog Heart
A09	Turtle Heart
BSL5	Components of the ECG (Lead II)
BSL6	Leads I, II, III & Einthoven's Law
H08	Dive Reflex—Active Learning
H23	Signal Averaged ECG
H32	Heart Rate Variability

**Pulmonary Function**

BSL12	Pulmonary Function: Vol. & Capacities
BSL13	Pulmonary Flow Rates: FEV and MVV

**Neurophysiology**

A01	Frog Pith & Prep
A02	Frog Nerve
A06	Cockroach Nerve
A08	Earthworm Action Potential
A14	CPG Hornworm
BSL3	EEG Relaxation & Brain Rhythms
BSL4	Alpha Rhythms in the Occipital Lobe
BSL10	EOG Eye Movement, Saccades & Fixation
H10	EEG & Hemispheric Asymmetry
H12	EOG Saccades & Displacement
H13	Visual Tracking vs. Imagination
H14	Ocular Fixation while reading
H15	Ocular Fixation while viewing an image

See page 43 for a description of all available lessons.

**Increase your lab options with...**

Cardiac Output Sensor	SS31L, p. 29
O <sub>2</sub> & CO <sub>2</sub> Analysis Module	GASSYS2-EA, p. 34
Temperature Transducer—Immersible	SS8L, p. 28
Transducer Accessory Pack	BSLPHA-TA, p. 42

# intro to bsl hardware

The Biopac Student Lab offers a comprehensive line of electrodes, transducers and accessories that work with the MP35 high performance data acquisition unit for a wide variety of applications on human, animal, organ or tissue preparations.

Electrodes are easy to use and transducers are durable and simple to handle. “Simple Sensor” connectors detect which electrode or transducer is plugged into each channel, and generate an on-screen prompt if students mistakenly plug in the wrong transducer for a lesson.

The BSL software contains presets that simplify setup for each transducer and signal type offered (i.e., Temperature, Airflow, Dissolved O<sub>2</sub> and ECG, EEG, EMG). Presets keep the hardware student-friendly...just connect and collect. You can easily customize presets to match your protocol and existing transducers.

If you have legacy transducers that you'd like to continue using, you can select from more than 20 interface options for manufacturers like Narco, Gilson, Lafayette, Harvard, and Vernier (see page 39). The BSLs interfacing flexibility can save a great deal of money and increases the system's usefulness.

BSL hardware can be ordered individually or as a system package.

- Packages contain everything—including consumable items—necessary to run a specified number of lessons, specific to that particular discipline.
- The BASIC, ADVANCED and ULTIMATE Systems have been offered since the Biopac Student Lab was launched in 1995, and are now in use in thousands of labs worldwide.
- The new CORE packages are offered to meet the diverging requirements of specific departments.
- Any package can be used as a starting point to develop your own custom system.

Please contact a **Biopac Student Lab Specialist** to request a formal quotation and to learn more about our quantity discounts. If you prefer, you can build a quote online at [www.biopac.com](http://www.biopac.com).

**BSL Hardware**

Data Acquisition Unit	page 24
Stimulators	page 25
Transducers (“SS Series”)	page 26-31
Stimulus Presentation	page 32
Noninvasive Blood Pressure	page 32
Tissue Bath Stations	page 33
Airflow & Gas Analysis	page 33-35
Electrodes	page 36-38
Interface Options	page 39
Core Package Summary	page 41



**IMPORTANT USAGE NOTICE**

BIOPAC Systems, Inc. instruments, components, and accessories are designed for educational and research oriented life science applications and investigations. BIOPAC Systems, Inc. does not condone the use of its instruments for clinical medical applications. Instruments, components, and accessories provided by BIOPAC Systems, Inc. are not intended for the diagnosis, cure, mitigation, treatment, or prevention of disease.

- See **Compliance & Ratings** on page 24 for details.





## Data Acquisition Unit - MP35

The MP35 data acquisition unit is the heart of the Biopac Student Lab System and is included in all BSL System orders. The unit connects to your computer using a USB port. The four universal, analog input channels are human safe (certified to IEC60601-1) and can record data from a wide range of signal sources, including: biopotential signals such as ECG, EEG, EMG, and EOG; transducer signals such as force, pressure, temperature, pH, and bioimpedance; and microelectrode signals from intra- and extracellular sources. Built-in control features support low- and high-voltage stimulation, trigger pulses, sound output and digital I/O options. Use the BSLSTM Stimulator for human subjects, or the new Low Voltage Stimulator (SS58L) for animal studies. The system can also record from and send digital trigger information to other software and devices such as SuperLab® and E-Prime®, switches, and relays. The MP35 fast processor, coupled with a variety of triggering functions, allows the system to be used like an oscilloscope, chart recorder or X/Y monitor.

The BSL System software controls the MP35 hardware to create a comprehensive teaching system. The teaching system, centered around dozens of lesson experiments, can be applied to complement a wide range of curriculum from human physiology to biomedical engineering. Powerful measurement tools help extract meaningful data. The BSL System software can automatically calculate many measurements from raw data, such as heart rate, systolic and diastolic blood pressure, lung volume, etc. See the BSL Software overview (pages 2-7) to learn more about the power and flexibility of the MP35 and the BSL System.

### BSL System Safety Compliance & Ratings

The BIOPAC MP35 System and accessories are safe for use on human, animal, organ and tissue preparations.



The MP35 System was tested by an accredited product safety testing and certification agency and classified as Class II Type BF equipment.

The MP35 System complies with applicable requirements for the following product and safety standards:

- CE mark
  - IEC 60601-1
  - EN 60601
  - EMC: IEC 60601-1-2
- UL 60601-1
- CAN/CSA-C22.2 No. 601.1-M90
- JIS T 1001



## Application Programming Interface

Student software developers can use BIOPAC API options to design and execute their own programs to control BIOPAC hardware or analyze data in other programs.

### BIOPAC Hardware API - BHAPI

Control the BIOPAC MP35 acquisition unit. Use API functions to: acquire data; acquire at different sample rates; set triggers; get the MP35 status; use the Analog Output channels or the Digital I/O.

- Reference manual and sample programs available for C/C++, C#, LabVIEW, MATLAB, and VB.NET.

### BIOPAC File Format API - ACKAPI

Use the base functions of this software library in a variety of combinations to use and parse specific data from BIOPAC's binary file format in other analysis programs. Retrieval options include: channel information; samples by segment of a specified channel; all samples of a specified channel; a particular sample of a specified channel; samples by time slice of a specified channel; marker information; text of a specified marker.

- Designed to be compatible with Microsoft C++ and Microsoft Visual Basic.

## MP35 Specifications

<b>Analog Inputs</b>	Front panel DSUB 9f labeled "CH1"-"CH4"
Number of channels:	4 isolated, human-safe universal biopotential/transducer input amplifiers
A/D resolution:	24-bit
Gain ranges:	10x to 50,000x in 12 steps
Input voltage range:	Gain dependent 400 microvolts-2 V peak-peak ( $\pm 10$ V with SS70L, page 39.)
Input Noise:	0.222 $\mu\text{V}_{\text{RMS}}$ (0.5 Hz-35 Hz)
CMRR:	85 dB minimum
Filters:	Programmable analog and digital (IIR) filters; automatic or user-adjustable
<b>Analog Output</b>	Back panel DSUB 9m labeled "Analog Out"
Voltage Output:	Range 0-4 V; Resolution 12-bit
Pulse Output:	Width: variable Repetition: variable
Pulse Level:	TTL-compatible with BSLSTM Stimulator: 0-10 V or 0-100 V with SS58L Stimulator: -10 V to +10 V
<b>Input Triggering Options</b>	
External Trigger:	Back panel BNC labeled "Trigger" TTL positive or negative edge
Analog Trigger:	Any of four input channels (front panel "CH1"-"CH4"); Threshold going positive or negative
Digital trigger:	Any of 8 input lines (back panel DSUB 25m)
<b>Sample Rate</b>	400 k samples/second aggregate across four channels - 100 k s/s on all four channels for short durations - 100 k s/s on three channels continuously - 50 k s/s on all four channels continuously
<b>Electrode Check</b>	
Impedance Range:	0-100 K $\Omega$ ; Checks electrode impedance from Vin+ and Vin- leads with respect to GND lead.
<b>Serial Interface Type</b>	USB
<b>Certification</b>	Complies with IEC60601-1 EMC complies with IEC60601-1-2 CE Marked
<b>Dimensions</b>	7 cm x 29 cm x 25 cm
<b>Weight</b>	1.4 kg

(See [www.biopac.com](http://www.biopac.com) for detailed specifications)

## Biopac Student Lab Stimulator - BSLSTM



The BSL stimulator can be used on Human, Animal, Organ or Tissue Preparations for the following physiological measurements:

- Twitch sub-threshold
- Tetanic contraction
- Twitch threshold
- Muscle tension/length versus force
- Maximum twitch response
- Fatigue
- Single twitch, summation
- Nerve conduction velocity
- Field Stimulation

### Stimulator Features

- Safety locking key switch to establish the operating range
- Output via the front panel BNC (female) connector
- Set the stimulus voltage level and view it on the digital display
- Output pulse trains (1-254 pulses), continuous or individual pulses
- Lock pulse width and frequency safety limits for your protocol
- Red warning LED flashes when a stimulus pulse is output
- Connects to the Analog Out on the back of the MP35
- Reference Output Cable: Stimulus marker output cable connects to any of the four analog input channels on the MP35 to record the stimulator marker pulse
- Reference Switch controls the output signal pulse width to Actual or Fixed (15 ms)
- Diagnostic test switch sends a fixed pulse width of 2.5 ms

The BSL Stimulator plugs directly into the MP35 and is controlled by the BSL PRO software. The PRO software is used to set up the stimulation pulse width, frequency and output (single pulse or series) options, and automatically notes any change to the stimulator parameters (e.g. pulse, width, and frequency) by inserting an event marker with descriptive text at the point of change. The stimulator Preset automatically scales the display for pulse amplitude. The BSL Stimulator can be set to start automatically with the recording, or manually at any point during the recording. Stimulus data can be displayed and monitored independently of the sample rate or stimulus frequency. Use with the HSTM01 (page 28) for human subjects.

### BSLSTM Specifications

<b>Pulse level amplitude:</b>	10 V Range or 100 V Range
<b>Range (selectable):</b>	.025-10 V or .12-100 V
<b>Pulse width:</b>	0.05-100 milliseconds
<b>Pulse repetition:</b>	5 seconds-0.5 milliseconds (0.2-2,000 Hz)
<b>Stimulator isolation:</b>	2,000 V <sub>RMS</sub> DC (HI POT test)
<b>Capacitance coupling:</b>	60 pF
<b>Power requirements:</b>	BSLSTMB for MP35: no additional power required BSLSTMA for MP30: 12 V DC 1 Amp adapter (included)
<b>Fuse:</b>	250 V fast blow 2 Amp
<b>Module dimensions:</b>	16 cm x 16 cm x 5 cm (610 grams)
<b>Compliance:</b>	IEC-60601-2-10 section 51.104

(See [www.biopac.com](http://www.biopac.com) for detailed specifications)

*BSLSTMB - For MP35 Systems*

*BSLSTMA - For MP30 Systems*

See Lessons A02, A03, A08, A09, H03, H06, H09, page 43.

## Low Voltage Stimulator - SS58L



SS58L



SS58L connected to MP35 with BSLCBL2A.

Use the Low Voltage Stimulator with any electrode or lead with a BNC connector (such as needle electrodes or clip leads, page 38)

for direct stimulation of animal or tissue preps. Interface with nerve chambers via BSLCBL3A or BSLCBL4B, page 39. Use with headphones or speakers for sound output. Control the stimulus with the Output Control option of the BSL PRO software. You can monitor the output directly on the computer without any external cable.

### SS58L Specifications

<b>Interface:</b>	MP35 Analog Out port (unisolated)
<b>Pulse level:</b>	-10 V to +10 V, software adjustable in 5 mV increments
<b>Pulse width:</b>	0.05-100 milliseconds
<b>Pulse repetition:</b>	5 seconds-0.1 millisecond (0.2-10,000 Hz)
<b>Power:</b>	No additional power required

See Lessons A02, A03, A08, A09, page 45.

### See these other Output Options:

- OUT1 Headphones, page 26
- HSTM01, Stim. Electrode for humans, page 28
- TSD122 Stroboscope, page 32
- STP35W Stimulus Presentation System, page 32
- ELSTM2 Stim. Electrodes, page 38
- BSLCBL7/11/12 Clip Leads, page 38



## Shielded Lead Set - SS2L

**General-purpose electrode lead—two SS2Ls are included with every BSL System\***

The SS2L cable connects disposable electrodes to the MP35 unit to measure biopotential signals. Each lead set has three pinch leads that snap directly onto standard disposable electrodes (such as the EL500 series electrodes): Red (positive), white (negative), and black (ground). Each 1-meter pinch lead terminates in a yoke connected to a 2-meter cable (total length 3 meters). Used in over 30 Lessons to record ECG, EEG, EGG, EMG, EOG, etc., see page 46. \*Core Animal and Core Pharmacology, substitute BSLCBL8 for SS2L. See also SS1LA, BSLCBL8 and BSLCBL9 (page 36).





## EDA Lead Transducer for Disposable Setups - SS57L

The SS57L EDA Lead Transducer snaps to two disposable EDA (isotonic gel) electrodes (EL507 on page 37). See the EDA (GSR) Transducer (SS3LA on page 28) as an alternative reusable option.

Range: .1-100  $\mu$ Mho (normal human range is 1-20  $\mu$ Mho)  
 Excitation: 0.5 V DC (constant)  
 Pinch Leads: Red (+), Black (GND)

See Lessons BSL9, BSL14, H11, H24, page 44.



## Pulse Photoplethysmogram - SS4LA

Measure blood density changes in the fingertip or other body locations, caused by varying blood pressure. Record the pulse pressure waveform. Attach to the finger by the Velcro® strap or tape to other body parts.

Sensor type: IR  
 Dimensions: 16 mm (long) x 17 mm (wide) x 8 mm (high)

See Lessons BSL7, H05, pages 43-44.



## Respiratory Effort Transducer - SS5LB

Record respiration via chest or abdominal expansion and contraction. The strap presents minimal resistance to movement and is extremely unobtrusive. The novel (non-Piezo) design permits the recording of arbitrarily slow respiratory activity.

Response: True DC  
 Circumference Range: 10 cm - 140 cm (increase with a longer nylon strap)  
 Dimensions: 95 mm (long) x 47 mm (wide) x 15 mm (thick)

See Lessons BSL8, BSL9, pages 43-44.



## Fast-response Thermistor - SS6L

Measure small variations in temperature, either on the skin surface or in an air stream. Record temperature changes in airflow during breathing to indicate respiration rate. Attach to the skin surface with surgical tape (TAPE1).

Response time: 0.6 sec  
 Max operating temp: 100°C  
 Accuracy & Interchangeability:  $\pm 0.2^\circ\text{C}$   
 Compatibility: YSI® series 400  
 Dimensions: 5 m (long) x 1.7 m (diameter)

See Lessons BSL8, BSL15, pages 43-44.

\*See other Temperature Sensors on page 28 (SS7L and SS8L).




## Hand Switch - SS10L

Use this hand switch for remote event marking or for psychophysiological response tests. Monitor switch data as an input channel. Connects to any analog input channel. Switch type: momentary push-button.

See Lessons BSL11, H11, H16, H24, H27, H30, pages 43-44.

\*See digital switch options for the MP35 unit on page 30 (SS53L-SS55L).

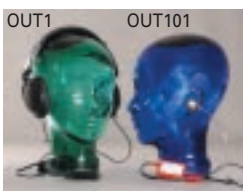

## Headphones - OUT1

Use these headphones to listen to real-time physiological signals (like EMG). Connect to the Analog Out port on the MP35. Weight: 85 grams

See Lessons BSL1, BSL2, BSL11, H09, H16, H31, pages 43-44.

\*See the TubePhone (OUT101) on page 31.

## connecting to the mp system



Unless otherwise specified, all transducers connect directly to the MP35 acquisition unit and have a 3-meter cable. All BIOPAC "Simple Sensor" transducer connectors are DSUB 9m for analog input and DSUB 9f for output. There is only one way to plug Simple Sensor connectors into the MP35, so you

don't have to worry about plugging things in upside down or into the wrong socket. We offer a wide range of transducer interface connectors so that you can connect your existing equipment, or you can build your own interface by using one of our custom interface kits (see page 39).

## Airflow Transducer (Med. Flow) - SS11LA



Airflow Transducer for resting human & light exercise. Use the SS11LA to perform a variety of tests relating to airflow and lung volume; integrate the airflow signal to obtain volume measurements. The pneumotach can be mounted on a camera tripod and has a clear, removable head for sterilization (using Cidex or a dishwasher) and replacement. Connects to industry standard bacteriological filters (AFT1) and disposable mouthpieces (AFT2). For hygiene, do not share disposable mouthpieces and disposable filters.

Flow Rate:  $\pm 300$  liters/min  
 Dead space: 93 ml  
 Dimensions: Head: 82.5 mm diameter x 101.5 mm length  
 Handle: 127 mm (long) x 23 mm (thick) x 35 mm (wide)  
 Ports: 22 mm ID/30 mm OD

The following consumable items are required for the SS11LA Airflow transducer to perform the pulmonary function lessons: AFT1 Disposable Filter, AFT2 Disposable Mouthpiece, AFT3 Disposable Noseclip, AFT6 Calibration Syringe. See page 34 for Airflow Accessories. Available in bulk, low-cost quantities for student accessory packs.

### Replacement Airflow Head - RX117

The RX117 is a sterilizable replacement transducer head for the SS11LA Airflow Transducer.

See Lessons BSL12, BSL13, BSL15, BSL18, H19, pages 43-44.

\*See High Performance Air Flow options on page 35.

## Variable Range Force Transducer - SS12LA



Use the SS12LA force transducer for tissue bath and isolated organ experiments, frog gastrocnemius and human finger twitch experiments. The SS12LA will operate in five ranges (50 g, 100 g, 200 g, 500 g or 1,000 g),

is extremely stable and incorporates impact and drop shock protection to insure against rough laboratory handling. The SS12LA mounting rod can be attached in three different locations, two on the top and one on the end surfaces of the transducer, to provide a variety of mounting options. The SS12LA includes two aluminum S-hooks: one has a 1.26 mm (.032") wire diameter; the other is 2 mm (.051").

Sensitivity:

Range	Noise (1 Hz LP)
50 g	1.0 mg
100 g	2.0 mg
200 g	4.0 mg
500 g	10.0 mg
1,000 g	20.0 mg

Temperature Range: -10°C to 70°C

Dimensions: 19 mm x 25 mm x 190 mm

See Lessons A02, A04, A05, A09, A15, H06, pages 43-45.

\*See the Tension Adjuster (HDW100A) on page 31.

See the Fixed-Range Force Transducers (SS63L-SS66L) on page 31.

## Pressure Transducer - SS13L



Measure direct arterial or venous blood pressure in animals or record pressure changes within a closed system (such as an organ or tissue bath system). Connect tubing via male Luer-lock fittings. The disposable transducer has a 30 cm cable that attaches to a reusable cable. Supplied non-sterile but can be cold sterilized.

Range: -50 to 300 mm Hg  
 Operating temperature: 10°C to 40°C  
 Transducer dimensions: 67 mm long X 25 mm wide

### Replacement Element - RX104A

The RX104A replacement element for the SS13L Pressure Transducer does not include the Simple Sensor connector and cable.

See Lessons A10, H04, pages 44-45.

## Blood Pressure Cuff - SS19L



Use to measure blood pressure via oscillometric or auscultatory methods. Use with the SS30L to record BP and Korotkoff sounds. The BSL System records the pressure of the cuff during BP measurements. This assembly

includes an adult blood pressure cuff (RX120D), pump bulb, mechanical indicator, and pressure transducer. Alternate cuff sizes are available (order as RX120A-F).

Manometer accuracy:	$\pm 3$ mmHg		
Pressure range:	20 mmHg to 300 mmHg		
Cuff	Circumference	Width	Length
<b>RX120A</b>	9.5-13.5 cm	5.2 cm	18.5 cm
<b>RX120B</b>	13.0-19.0 cm	7.5 cm	26.1 cm
<b>RX120C</b>	18.4-26.7 cm	10.5 cm	34.2 cm
<b>RX120D</b>	25.4-40.6 cm	14.5 cm	54.0 cm
<b>RX120E</b>	34.3-50.8 cm	17.6 cm	63.3 cm
<b>RX120F</b>	40.6-66.0 cm	21.0 cm	82.5 cm

See Lessons BSL16, H04, pages 43-44.

\*See BSLUPG16&17 (page 40) to order BP Cuff with Stethoscope.



## Hand Dynamometer - SS25LA

Use this in isolation or combine with EMG recordings for in-depth studies of muscular activity. The lightweight, ergonomically designed transducer provides direct readings in kilograms or pounds. Use to measure grip force.

Dimensions: 17.78 cm x 5.59 cm x 2.54 cm  
 Weight: .323 kg  
 Range: 0-90 kg

See Lesson BSL2, page 43.




## Multi-Lead ECG - SS29L

The multi-lead ECG cable connects to standard snap-connector disposable electrodes (EL503 series) to simultaneously record Leads I, II, III, aVR, aVL, aVF, plus one alternating precordial chest lead V(1-6) for 12-lead studies. Incorporates a Wilson terminal. Each 1-meter pinch lead terminates in a yoke connected to a 2-meter cable that has three Simple Sensors (three input channels required).

See Lesson H01, page 43.



Connect directly to the MP35 Data Acquisition Unit

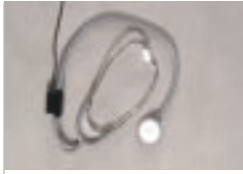

## Electronic Stethoscope Transducer - SS30L

This is a standard clinical stethoscope with a built-in electronic microphone to simultaneously capture sound. Listen to heart sounds and Korotkoff sounds, and simultaneously record the sound data (a variety of acoustical signals can be recorded). When recording ECG, you can correlate the timing of the heart sounds with the cardiac cycle. Use with the SS19L Blood Pressure Cuff to record Korotkoff sounds for easy determination of systolic and diastolic blood pressure. Microphone Bandwidth: 20-100 Hz (does not interfere with stethoscope earphones).

See Lessons BSL16, BSL17, H04, H21, pages 43-44.

\*See also the Physiological Sounds Microphone (SS17L, below) for a higher bandwidth alternative and the Speech Frequency Microphone (SS62L, page 30). Order BSLUPG16&17 Upgrade Pack (page 40) to get BP Cuff (SS19L, page 27) and Stethoscope.



## Human-safe Stimulation Electrode - HSTM01

Provides a superior degree of safety and comfort when using the BSLSTM Stimulator for human stimulation. The ergonomic design allows the user to focus on electrode placement when locating a subcutaneous nerve. Stimulus presentation is allowed by pressing the red safety switch and stopped by releasing it. Complies with IEC 60601-2-10 for limitation of stimulation signal and interfaces with the BSLSTM stimulator (BNC connector).

Isolation Voltage: 1500 V<sub>RMS</sub>  
Isolation Capacitance: 120 pF

See Lessons H03, H06, pages 43-44.




## EDA (GSR) Transducer - SS3LA

Measure electrodermal activity and response (galvanic conductance/GSR). The built-in, reusable electrodes fit around the tip of a person's finger and attach via Velcro® straps or can be taped to any other body part.

Electrode Type: Ag-AgCl, shielded  
Excitation: 0.5 V DC (constant)  
Range: 0.1-100 μMho (normal human range is 1-20 μMho)  
Surface Area: 6 mm diameter contact area  
Gel Cavity Depth: 1.66 mm  
Dimensions: 16 mm (long) x 17 mm (wide) x 8 mm (high) [each]

See Lessons BSL9, BSL14, H11, H24, page 44.

\*See the EDA (GSR) Disposable Option (SS57L) on page 26.

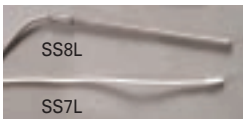

## Waterproof Temp. Probe - SS7L

Use this vinyl probe for core (oral/rectal) temperature recordings.

Response time: 1.1 sec  
Max operating temp: 100°C  
Accuracy & Interchangeability: ±0.2°C  
Compatibility: YSI® series 400  
Dimensions: 3 mm (dia) x 9.8 mm (long)

\*See the Fast-response Thermistor (SS6L) on page 26.

## Liquid Immersion Temp. Probe - SS8L

Use this stainless steel probe for dry or wet bath temperature measurements.

Response time: 3.6 sec  
Max operating temp: 100°C  
Accuracy & Interchangeability: ±0.2°C  
Compatibility: YSI® series 400  
Dimensions: 4 mm (dia) X 115 mm (long)



## Displacement Transducer - SS14L

Record very slight movements (up to 100 mm) in a range of physiological preparations. Incorporates a semi-isotonic strain gauge (500 ohm silicon) and a nickel-plated cantilever beam device (27 cm) for holding the transducer. Features high linearity.

Sensitivity range: Up to 100 mm  
Dimensions (mm): Body: 95 x 25 x 25; Blade: 305 x 1.27 x 0.6; Support rod: 127 x 9.5 (dia)



## Physiological Sounds Microphone - SS17L

Use to record a variety of acoustical signals, including heart sounds and sounds associated with rubbing or grinding (e.g., Bruxism). Use with the SS19L noninvasive Blood Pressure Cuff to record Korotkoff sounds for easy determination of systolic and diastolic blood pressure. Use as a higher bandwidth alternative to the electronic stethoscope (SS30L, above).

Microphone Bandwidth: 35 Hz to 3,500 Hz  
Housing: Stainless Steel  
Transducer Weight: 9 grams

See Lesson H04, page 44.

\*See also Speech Frequency Microphone SS62L, page 30.

## Goniometers & Torsiometers - SS20L-SS24L



Goniometers measure bending strain along or around a particular axis to transform angular position into a proportional electrical signal. All goniometers have a telescopic endblock that compensates for changes in distance between the two mounting points as the limb moves. Goniometers can be attached to the body surface using tape (TAPE1 or TAPE2, page 36).

- Twin-axis goniometers (SS20L/21L) are dual output devices and measure angular rotation about two orthogonal planes simultaneously (e.g. wrist flexion/extension and radial/ulnar deviations).
- Torsiometers (SS22L/23L) measure angular twisting (as on the torso, spine or neck) as opposed to bending and measure rotation about a single axis (e.g. forearm pronation/supination).
- The single axis goniometer (SS24L) measures finger, thumb or toe joint movement and will measure the angle in one plane only.

	<b>SS20L</b>	<b>SS21L</b>	<b>SS22L</b>	<b>SS23L</b>	<b>SS24L</b>
Type:	Goniometer	Goniometer	Torsiometer	Torsiometer	Goniometer
Channels:	2	2	1	1	1
Max Length:	110 mm	150 mm	110 mm	150 mm	35 mm
Min Length:	75 mm	130 mm	75 mm	130 mm	30 mm
Range:	±180°	±180°	±90°	±90°	±180°
Typical Placement:	wrist or ankle	elbow, knee or shoulder	neck	torso or spine	fingers, thumb or toes
Weight:	17 g	19 g	17 g	19 g	8 g

See Lessons H17, H28, H35, pages 44-45.

## Accelerometers - SS26L-SS27L



These tri-axial accelerometers measure accelerations in the X, Y, and Z directions. For simultaneous measurement, three input channels are required. The design conforms to body contours and includes a Velcro® strap for easy attachment.

Range (Output): SS26L: ±5 g, SS27L: ±50 g

Bandwidth: DC - 500 Hz (-3 dB)

Dimensions: 33 mm (long) x 28 mm (wide, at base) x 19 mm (high)

**SS26L:** Optimal for measuring slow movements, such as walking.

**SS27L:** Optimal for measuring quick movements, such as swinging a tennis racket.

## Heel/Toe Strike - SS28L



Use this transducer to record heel and toe strike activity as the subject walks. The heel/toe strike data is recorded as a single channel; the heel strike generates a negative deflection and the toe strike results in a positive deflection. Two force sensitive resistors (FSR) attach to the sole of a shoe; use two SS28L transducers to simultaneously record from both feet.

Nominal Contact Force: 200 g to indicate heel/toe strike

Attachment: TAPE1, TAPE2

FSR Dimensions: 18.3 mm wide x 0.36 mm (thick) x 30 cm long

FSR Active Area: 12.7 mm (dia)

Cable Length: 7.6 meters

## Cardiac Output Sensor - SS31L



Measure cardiac output noninvasively using electrical bioimpedance techniques; the sensor is suitable for human and animal measurements. Use the sensor to record stroke volume and CO before and after exercise, or during a psychophysiology test. The sensor provides a noninvasive, powerful cardiovascular hemodynamic demonstration that allows students to see the real-time changes in stroke volume and cardiac output. Typically used with disposable bioimpedance strip electrodes (EL506 on page 37), but can function with spot or band electrodes, reusable electrodes, or needle electrodes.

Number of Channels: 2 - Impedance ( $Z_0$ ) and  $dZ/dt$

Operational Frequency: 100 kHz


Outputs: Impedance  $Z_0$  0-100 ohms  
 $dZ/dt$  ±20 ohms/sec

Configuration: Tetrapolar—4 Electrodes

Excitation: 400  $\mu$ A RMS constant current

See Lesson H21, page 43.



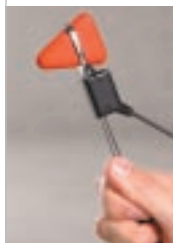

## Reflex Hammer - SS36L

This is a classic reflex hammer with a transducer attached to perform reflex measurements. It uses a Taylor Hammer®—the most common type of reflex hammer used by doctors—and incorporates electronics to record the time and the relative strength of the impact. This allows students to measure how much of an impact is needed to elicit a response.

See Lesson H28, page 44.

## Signal Processing Breadboard - SS39L

The Bioengineering Breadboard Lab consists of circuitry hardware and lessons (with schematics and design notes) that demonstrate a very important subset of circuit design for recording and processing physiological signals. Includes electrode signal lead interface. Use with SS60L cables to record circuit performance on more than one channel. Students will use the MP35 and BSL PRO software to evaluate their designs.



Included Hardware:

- 1 - Breadboard
- 1 - Power/Signal Cable (includes fuses with built-in, automatic reset)
- 1 - Parts kit including op-amps, capacitors, diodes, resistors, and jumper wires as required to complete projects

Lesson Modules:

- |  |   |
|--|---|
| Lab 1: Square Wave Oscillator                | Lab 5: Notch Filter for 60 Hz Rejection                       |
| Lab 2: Instrumentation Amplifier             | Lab 6: QRS Detection: Band Pass Filter                        |
| Lab 3: High Pass Active Filter               | Lab 7: QRS Detection: Absolute Value Circuit                  |
| Lab 4: Active Gain Block and Low Pass Filter | Lab 8: QRS Detection: Low Pass Filter and Overall System Test |
- Combine modules to build a complete ECG Signal Processor.

## Optional Signal Processing Cable - SS60L

Use this signal cable to record additional channels with the SS39L Signal Processing Breadboard.

See Lessons H25, H26, page 45.

## Differential Pressure - SS40L-SS42L



These transducers interface a variety of small animal breathing circuits to the MP35 for air pressure monitoring. The transducers are extremely sensitive and come in three ranges to suit a number of different applications. Included with each SS45L-SS52L (page 33).

Range:

- SS40L:** ±2.5 cm H<sub>2</sub>O
- SS41L:** ±12.5 cm H<sub>2</sub>O
- SS42L:** ±25.0 cm H<sub>2</sub>O

Dynamic Response:

100 Hz

Connection ports/ID:

3 mm to 4.5 mm tubing accepted

Dimensions:

8.3 cm (high) x 3.8 cm (wide) x 3.2 cm (deep)

Weight:

76 grams

## Variable Assessment Transducer - SS43L



Use this handheld, slide control transducer to record subjective responses for a variety of different stimuli. Use multiple transducers to allow several people to simultaneously answer the same question or otherwise respond to stimuli. Easily customize the response scale by inserting parameters into the scale sleeve on the front of the unit. 7.6 meter cable.

## Digital Switches - SS53L-SS55L



Use for remote event marking or to externally trigger data acquisition for psychophysiological response tests. Monitor switch data as a digital input channel. The following switches inter-

face with the I/O port on the rear of the MP35 unit for digital input.

### Hand Switch, Digital - SS53L

Dimensions: 19 mm (dia) x 63 mm (long)  
Cable Length: 2 meters  
Connector Type: DSUB 25f

### Foot Switch, Digital - SS54L

Dimensions: 69 mm (wide), 90 mm (long), 26 mm (high)  
Cable Length: 1.8 meters  
Connector Type: DSUB 25f

### Eight-channel Marker Box, Digital - SS55L

Independently mark events, or provide responses, on up to eight channels simultaneously. Assign separate digital channels as event markers for individual analog input channels. Easily customize the switch indicators by inserting parameters into the label sleeve on the front of the unit.

Dimensions: 19 cm (wide), 11 cm (deep), 4 cm (high)  
Cable Length: 2 meters  
Connector Type: DSUB 25f

See Lessons H11, H16, H24, H27, H30, pages 43-44.

## Finger Twitch Transducer - SS61L



Use this transducer to record finger twitch responses from human subjects receiving electrical stimulation (using the HSTM01, page 28). The transducer conforms to the shape of the finger and attaches via a Velcro® strap and tape.

Transducer Dimensions: 14.6 cm (long), 0.50 cm (wide)  
Weight: 6 g

See Lesson H06, page 43.

## Speech Frequency Microphone - SS62L



Use this precision microphone for speech frequency analysis and other acoustic studies. Requires continuous high-speed sample rate - use with the MP35 only.

Frequency Range: 60-12,000 Hz  
Impedance: 600 Ohms  
Type: Cardioid  
Cable: 6 meters  
On/Off Switch: none

\* See also Physiological Sounds Microphone (SS17L) and Electronic Stethoscope Transducer (SS30L) on page 28.

## Fixed-Range Force Transducers - SS63L-SS66L



Use force transducers for *in vitro* tissue bath studies, *in vivo* force measurements and other applications where low noise, accuracy and repeatability are critical. The most sensitive units are ideal for use with small preparations such as aortic rings. When fine tension and position

adjustments are required, use the Tension Adjuster (HDW100A, page 31).

Part #	Full Scale Range (FSR)	Noise with 10 Hz LP Filter	Noise with 1 Hz LP Filter
SS63L	50 g	2.5 mg	1.0 mg
SS64L	100 g	5.0 mg	2.0 mg
SS65L	200 g	10.0 mg	4.0 mg
SS66L	500 g	25.0 mg	10.0 mg

Temperature Range:	-10°C to 70°C
Weight:	250 grams
Mounting Rod:	9.5 mm (dia)- variable orientation
Dimensions:	100 mm (long) x 19 mm (wide) x 25 mm (thick)

See Lessons A02, A04, A05, A09, A15, H06, pages 43-45.

See also the Variable Range Force Transducer (SS12LA, page 27).

## Pneumogram Transducer - SS67L



Capture respiratory signals from small, unconscious animals—just lay them on top of the sensor pad. Ideal for MRI-applications. Affix to

subjects with single-sided adhesive (TAPE1, page 36). This unobtrusive, multipurpose pneumogram transducer can:

1. Noninvasively measure respiration—from a small mouse to large rodent.
2. Measure small pressing forces (like pinching fingers together) for Parkinson's evaluations.
3. Measure human smiling (with the sensor on the cheekbone).
4. Measure pulse when placed close to the heart.
5. Measure spacing and pressure between teeth coming together.

SS67L components:

- 1- Differential Pressure Transducer (SS41L, page 30)
- 1- Pneumogram Sensor (RX110, below)
- 1- Tubing (1.6m)

### RX110 - Replacement Sensor

The RX110 is a self-inflating pressure pad connected to tubing terminating in a Luer male connector. The RX110 sensor is included with the SS67L Pneumogram Transducer.



## pH Probe - SS68L

This double-junction pH probe measures pH within the range of 0-14. Use the pH Probe Adapter (BSL-TCI21, page 39) to use a third party pH probe with the MP35. Order probe only as RXPROBE01.

SS68L Components:	RXPROBE01 and BSL-TCI21
Output:	5 mV/pH (0 mV @ pH=7)
Weight:	99.22 g
Length:	3.25 cm
Diameter:	1.2 cm

## Dissolved O<sub>2</sub> Probe - SS69L



Measures dissolved oxygen levels. Includes electrolyte solution, replacement membrane cap and replacement O-ring.

Use the Vernier Adapter (BSL-TCI16, page 39) if you already have a Vernier Dissolved O<sub>2</sub> Probe. Order probe only as RXPROBE02.

SS69L Components:	RXPROBE02 and BSL-TCI16
Type:	Polarographic
Oxygen Range:	0-40 ppm
Electrodes:	Platinum and silver
Minimum sample velocity:	20 cm/sec.

See Lesson A07, page 45.



## Tubeophone - OUT101

Use the OUT101 with a stimulator module to deliver clicks in auditory evoked response applications (i.e., ABR). The tubeophone design consists of a monaural acoustic transducer

attached to a short, flexible, plastic tube, which fits into the subject's ear with the aid of a foam tip. Compares to TDH-39, 49 or 50 audiometric headphones.

Dimensions:	3.8 cm (wide) x 5 cm (high) x 1 cm (thick)
Cable termination:	6.3 mm (1/4") phone plug
Cable length:	1.8 meter

See Lesson H09, page 44.

\*See the Headphones (OUT1) on page 26

## Tension Adjuster - HDW100A



Use the HDW100A to adjust the tension between the preparation and a Force transducer (SS12LA on page 27 or SS63L-SS66L on page 31). The position adjuster is located on the top for easy access and smooth operation and mounts on typical lab stands. Vertical scales are provided for both metric and standard units. Works with most commercial force transducers.

Resolution:	0.0025 mm per degree rotation
Stand Clamp:	12.80 mm ID
Transducer Clamp:	14.60 mm ID
Dimensions:	93 mm (high) X 19 mm (thick) x 74 mm (deep)

See Lessons A02, A04, A05, A09, A15, page 45.

## Tension Adjuster Adapter - HDW200

This adapter allows third-party tension adjusters with an arm diameter of 6.35 mm (1/4") or less to work with BIOPAC force transducers.



# visual and auditory stimulus presentation

## SuperLab System - STP35W



The SuperLab package presents visual or auditory stimuli on a second computer screen, or auditory stimuli via headphones or speakers, and simultaneously

(1 ms resolution) sends trigger signals to the MP35 for data synchronization and collection purposes. Images and sounds are grouped and have a digital synch signal assigned to each group. As each image or sound is presented, a synch pulse is sent to the MP35. The BSL software identifies the pulses and provides the average response to the group stimulus and takes measurements over the response data such as max, min, mean, time of max, etc. The STP35W includes an interface cable, permitting up to eight synchronization signals (input or output) between the STP35W and the BSL System. Different trigger channels can be paired to different visual or auditory stimuli to perform sophisticated evoked response averaging tests (e.g., P300). The STP35W includes a six-pushbutton response box for performing accurate reaction time measurements. **Note: Second PC required.**

### The STP35W System includes:

- SuperLab Software (available for Windows or Mac)
- Digital I/O Card and Support Pack software (PCI slot required)
- Six-Pushbutton Response Box
- Pushbutton Keycap Color Change Kit
- Interface cable for MP35 System

See Lessons H24, H27, H30, H31, pages 43-44.

## Interface Options



### SuperLab Interface Cable for MP35 - STP35

If you already have SuperLab and an MP35 unit, you can use the STP35 Interface Cable to connect the two systems. The STP35 Cable interfaces with the I/O Port on the rear of the unit.

### Parallel Port Interface Cable - STP35A

Use to connect with systems that use the parallel (printer) port, such as E-Prime, MediaLab, DirectRT, and Inquisit.



### Stroboscope - TSD122

The stroboscope connects directly to the MP35 Input or Output channels for visual evoked response applications, and includes BSLCBL5 (page 39) to interface with the BSL System. The stroboscope operates from zero to 12,000 flashes a minute and can generate or respond to a

TTL synchronizing signal.

I/O Ports:	TTL (Sync input and output)-3.5mm phone jacks
Flash Duration:	30 $\mu$ sec
Power:	Battery, built-in, rechargeable
Battery Life:	60 hours at 100 strobes/sec
Interface Cable:	BSLCBL5 (included) connects to an analog input (CH1-CH4) to record the strobe flash as a pulse and trigger the MP unit.
Body Dimensions:	9.3 cm (wide) x 9 cm (high) x 23 cm (long)
Weight:	1.1 kg

TSD122C 120 V/60 Hz

TSD122D 220 V/50 Hz

See Lesson H22, page 44.

## 32 noninvasive blood pressure



### Noninvasive Blood Pressure Monitoring - NIBP100A

The NIBP100A is a noninvasive solution for automatic blood pressure monitoring, with accuracy comparable to an indwelling radial artery catheter. The system provides a continuous reading of BP (updates every 10-15 heartbeats) and will function while subjects undertake light exercise or psych analysis

tasks. The BSL System calculates systolic, diastolic and mean pressures in real time. The NIBP100A is easy to use—no complicated setup or calibration requirements!

### NIBP100A Specifications


Performance:	BP: 20-240 mmHg Min/Max; SD 8 mmHg Accuracy Pulse: 40-200 bpm Min/Max; $\pm$ 5 bpm or 10% Accuracy
Interface:	BSL-TCI5 (page 39; supplied)
Size (cm):	12.70 (h) x 11.43 (w) x 21.59 (l)
Weight:	2.04 kg with power cord and wrist module
Compliance:	FDA clinical standards and UL60601-1 and IEC 60601-2-30

### RXNIBPA - Replacement sensor

The sensor in the NIBP100A requires replacement every six (6) months. An internal processor monitors the age of the sensor and then automatically stops on the 6-month anniversary.

See Lesson H18, page 44.

# noninvasive blood pressure for animals

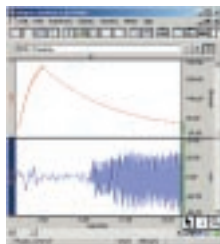

## Noninvasive Small Animal Tail Blood Pressure System

The NIBP200A incorporates a built-in pump that automatically inflates the blood pressure cuff to occlude the vessel. Once the pump reaches the inflation point it slowly deflates the cuff, providing a linear drop in pressure. A single pushbutton controls both the inflation and deflation cycles, making the system very operator friendly. The amplifier provides two analog outputs for pressure and pulse waveforms. The interface cables provide simple connection to your BIOPAC or third-party A/D hardware. An optional tail heater is available.

The NIBP200A includes:

- 1 - amplifier unit (110 V or 220 V)
- 1 - sensor (7.5 mm is included, but can be specified as 6 mm, 6.5 mm, 7 mm, 8 mm, or 8.5 mm when ordering)
- 1 - cuff (size 11 cm is included, but can be specified as 9.5 cm or 13 cm when ordering)
- 1 - restrainer (size medium is included, but can be specified as small, large, or mice when ordering)
- 1 - Pulse cable to BSL System
- 1 - Pressure cable to BSL System
- 1 - User's Manual

NIBP200A1 Noninvasive BP-Sm. Animal Tail 110 V

NIBP200A2 Noninvasive BP-Sm. Animal Tail 220 V

Optional accessories: TAILHEATA Tail Heating Unit 110 V / 60 Hz

TAILHEATB Tail Heating Unit 220 V / 50 Hz

## tissue baths



## Integrated Tissue Bath & Heater System - ITBS100

A modular, durable solution for your lab

Features include:

- Jacketed bath and reservoir in a range of volumes
- Integrated, heating circulator, programmable temperature range 20° - 44° C
- 500 ml/min circulation flow
- Movable micrometer-transducer assembly
- User-friendly display and controls
- One-switch control of fill/drain cycle
- Microprocessor control
- Low-level alarm for water temperature
- Acrylic, robust bath
- Small, lightweight setup

Components:

- 1 x Bath (specify 20 ml, 30 ml, or 50 ml)
- 1 x Reservoir (800 ml)
- 1 x Integrated Heater (1600 ml volume, 20° - 44° C)
- 1 x Circulator Pump (15 W; 500 ml/min)
- 1 x Micrometer (fits BIOPAC transducers)
- 2 x Triangle Tissue Clip (Stainless Steel; reorder as RXCLIP-TRI)
- 2 x Tissue Clip (Stainless Steel; reorder as RXCLIP)
- 1 x Tissue Holder (Stainless Steel; reorder as RXHOLDER-S)
- 1 x 3-way Rotary Valve
- 1 x Power Supply (specify 110 V/60 Hz. or 220 V/50 Hz)

See Lessons A05, A15, page 45.

\*Contact BIOPAC for modular, multi-channel tissue bath stations and independent heating circulator.

## Field Stimulation Electrode - BSLSTMHLD



This tissue holder with built-in field stimulation electrodes works with Tissue Bath Stations and the BSLSTM Stimulator.

Electrodes: Platinum rings  
 Ring size: 5 mm  
 Ring spacing: 2.0 cm  
 Connector: BNC

## airflow & gas analysis

### Airflow Transducers- SS45L-SS52L

\*See also the Medium Flow Airflow Transducer (SS11LA) on page 27.



Connect these Airflow Transducers directly to a breathing circuit or plethysmogram chamber to perform a variety of pulmonary measurements relating to airflow, lung volume and expired gas analysis of small animals or exercising humans. Each Airflow Transducer consists of a flow head coupled to a precision, highly sensitive, differential pressure transducer (SS40L) with 1.8 meters of tubing. Each flow transducer has an internal heating element that can be connected to an optional 6 V power supply (AC137, page 40). For airflow and lung volume measurements, connect a short airflow cannula to the flow head. For switchable or replacement head options, use the RX137 Series replacement part number cited.

Part	Replacement	Max. Range (ml/sec)	Dead Space (cc)	Output (µV/[ml/sec])	Ports ID; OD (mm)	Length (mm)	Animal Size; Weight
SS45L	RX137A	±12	0.1	25.700	1.35; 7.00	75	Small Mouse; 30 gm
SS46L	RX137B	±20	0.8	15.400	6.00; 7.00	75	Mouse; 50 gm
SS47L	RX137C	±60	0.9	5.780	6.00; 7.00	75	Rat/Guinea Pig; 350.0 gm
SS48L	RX137D	±150	2.0	2.100	9.00; 10.00	75	Cat/Rabbit; 750 gm
SS49L	RX137E	±350	4.0	0.924	10.00; 11.00	60	Small Dog; 5.5 kg
SS50L	RX137F	±1200	14.0	0.231	17.00; 19.00	60	Med. Dog; 15 kg
SS51L	RX137G	±3000	35.0	0.0963	28.00; 30.00	60	Large Dog; 25 kg
SS52L*	RX137H	±8000	80.0	0.0385	43.00; 45.00	60	Exercising Human

\*Requires one coupler (AFT11F, page 35) to interface with the GASSYS2 (page 34) and other airflow accessories.



# airflow & gas analysis

SPECIAL

## Gas Analysis Module - GASSYS2



O<sub>2</sub> and CO<sub>2</sub> Gas Analysis Module measures expired O<sub>2</sub> and CO<sub>2</sub> concentrations. Obtain real-time Oxygen Consumption (VO<sub>2</sub>), Basal Metabolic Rate

(BMR) and Respiratory Exchange Ratio (RER) measurements using the MP35 System with the GASSYS2 module. Available with



a 5-liter chamber or a 1-liter chamber for small children/medium sized animals. Chambers are modular so you can interchange them with one base module and

they are easy to clean. The GASSYS2 includes tubing (AFT7), coupler (AFT11E), non-rebreathing "T" valve (AFT22), and a separate 12 V power supply (AC100A), creating a low-cost solution for BSL users already using the SS11LA Airflow Transducer and accessories. Flap valve prevents expired air from returning to chamber, and gas calibration port allows O<sub>2</sub>/CO<sub>2</sub> sensor calibration.

CO<sub>2</sub> Range: 0-5%

O<sub>2</sub> Range: 0-25%

Note: Two couplers (2 x AFT11F) required when interfacing with the Airflow Transducer (SS52L).

GASSYS2-EA - Module with 5-liter chamber

GASSYS2-EB - Module with 1-liter chamber


RX-GASA - 5-liter chamber/screw fixture

RX-GASB - 1-liter chamber/screw fixture

See Lessons BSL18 and H19, page 43.

\*See gas calibration options on page 35 (GASCAL and GASREG). The GASSYS2 can measure both O<sub>2</sub> consumption and CO<sub>2</sub> production in real time. CO<sub>2</sub> production is measured to a 5% maximum for RER or RQ calculations, and real-time VO<sub>2</sub> measurements can be performed on subjects at rest or exercising. If your protocol examines CO<sub>2</sub> above 5% or requires high-speed response, contact BIOPAC to discuss the full range of gas analyzers.

## SS11LA accessories

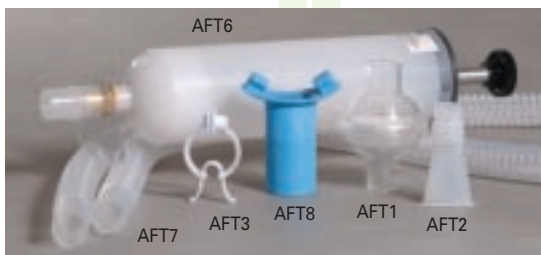

\*See SS11LA Airflow Transducer on page 27.

### Disposable Bacterial Filters (22 mm ID/OD) - AFT1

Use between the Airflow Transducer (SS11LA) and the 22mm disposable Mouthpiece (AFT2) to remove airborne bacteria.

Ports: 22 mm ID/OD.

AFT1 Pack of 10

AFT1-250 Pack of 250

See Lessons BSL12, BSL13, BSL15, BSL18, H19, pages 43-44.

### Disposable Mouthpieces (22 mm OD) - AFT2

Connects to the Airflow Transducer (SS11LA) via the disposable Bacterial Filter (AFT1).

AFT2 Pack of 10

AFT2-250 Pack of 250

See Lessons BSL12, BSL13, BSL15, BSL18, H19, pages 43-44.

### Disposable Nose Clips - AFT3

Plastic clip gently squeezes the nostrils shut. Use with the Airflow Transducer (SS11LA).

AFT3 Pack of 10

AFT3-250 Pack of 250

See Lessons BSL12, BSL13, BSL15, BSL18, H19, pages 43-44.

### BSL Accessory Pack - BSL-ACCPACK



Have students purchase their own disposable accessories for hygienic purposes and reduce the burden on department budgets—sell these student accessory packs through your school bookstore. Each pack includes

disposable (one-use only) items required to run BSL Lessons:

60 x EL503 Disposable Electrodes

1 x AFT3 Nose Clip

1 x AFT2 Disposable Mouthpiece

8 x ELPAD Abrasion Pads

1 x AFT1 Disposable Bacterial Filter

8 x EL507 EDA Electrodes

### Standard Syringe - AFT6

Use the 0.6 liter calibration syringe with the SS11LA Airflow Transducer or the GASSYS2 (incorporates a 22 mm OD, 15 mm ID coupler); use the AFT26 for increased calibration precision.

See Lessons BSL12, BSL13, BSL18 and H19, page 43.

### T-Valve - AFT22

Interfaces with the SS11LA Airflow Transducer and the AFT7 tubing via the AFT11E coupler. Suitable for static applications with low to medium flow rates. Requires AFT1 and AFT2 for proper operation. All ports are 22 mm OD; includes 22 mm OD coupler.

See Lessons BSL18 and H19, page 43.

### Reusable Mouthpiece - AFT8

Autoclavable—Use with the SS11LA Airflow Transducer to reduce the cost of disposable parts. 30 mm ID.


AFT8 Pack of 1

AFT8-10 Pack of 10

### Disposable Adult Facemask - AFT10

This disposable facemask connects to 22mm breathing circuits (22mm ID, 25mm OD).

Connects directly to the AFT1, AFT22 non-rebreathing T-valve or SS11LA Airflow transducer via tubing (AFT7) and two couplers (AFT11B and AFT11E). Includes hook-ring to secure adjustable head strap (AFT10S).

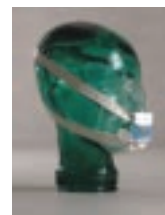

### Strap - AFT10S

This fully adjustable non-latex reusable head strap holds the AFT10 disposable facemask securely to the subject's head.

Use one strap to securely fasten the mask. (1 per pkg.)

## High Performance Airflow Accessories

The following items interface with the Airflow Transducer SS52L, (page 33), which is intended for exercise physiology and other high volume pulmonary function studies. To use with the SS11LA Airflow Transducer (page 27), use AFT7 tubing and AFT11E coupler.

### T-Valve - AFT21

High-performance, very low dead space, low airflow resistance valve; suitable for high airflow applications. Interface with the Airflow Transducer (SS52L) via the Tubing (2 x AFT7) and Coupler (AFT11F) and the GASSYS2. All ports are 35 mm OD and 30 mm ID; AFT11D coupler included to fit AFT4 filter.

### Head Support - AFT24



Use the head support when breathing directly into the AFT21 non-rebreathing T-valve for exercise physiology measurements. The AFT24 secures the AFT21 directly in front of the subject and minimizes the strain associated with valves and tubing.

See Lessons BSL18 and H19, page 43.


### Facemask with Valve - AFT25



Use the AFT25 facemask when performing gas analysis tests on an exercising subject or for other high airflow applications. The adult facemask has an integral high-performance non-rebreathing T-valve, and very low dead space and low flow resistance. Interface with the Airflow Transducer (SS52L) via tubing (2 x AFT7) and coupler (AFT11E) and the GASSYS2.

All ports are 35 mm OD and 30 mm ID.

See Lessons BSL18 and H19, page 43.

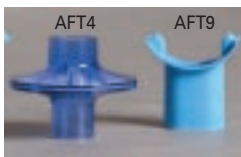

### Disposable Bacterial Filter - AFT4

Use with the Airflow Transducer (SS52L) and the 35 mm reusable mouthpiece (AFT9) to remove airborne bacteria when measuring bi-directional airflow.

### Reusable Mouthpiece - AFT9

Use with the Airflow Transducer (SS52L) or the 35mm bacterial filter (AFT4) to connect to the non-rebreathing T-valve (AFT21).

35 mm ID.

AFT9 Pack of 1 AFT9-10 Pack of 10

## Airflow Guide

Item	Applications	BIOPAC Part
Airflow Transducers (Pneumotachs)	Human/Animal respiration, pulmonary function, exercise physiology. See Lessons BSL11, BSL12, BSL18, H21, BSL15, H19 (page 43-44).	SS11LA (med), page 27 SS45L-SS52L, page 33
Calibration Syringes	Use the AFT26 for increased precision. See Lessons BSL11, BSL12, BSL18, H19 (page 43-44).	AFT6, page 34 • AFT26, page 35
Calibration Gas	Use to verify calibration settings.	GASCAL, page 35
Couplers	Interface tubes and ports. Fit to specified ID/OD sizes—more port sizes available upon request.	AFT11B, AFT11E, AFT11F, page 35
Facemask & Strap	Exercise physiology.	AFT10 and AFT10-S, page 34
Facemask with Valve	Exercise physiology, pulmonary function, metabolic rate (BMR). See BSL18 and H19 (page 43).	AFT25, page 35
Filters, Disposable	AFT1 for SS11LA. AFT4 for SS52L.	AFT1, page 34 • AFT4, page 35
Head Support	Exercise Physiology. See Lessons BSL18 and H19 (page 43).	AFT24, page 35
Mouthpieces	Disposable and reusable options.	AFT2, page 34 • AFT8, page 34 AFT9, page 35
Nose Clips	Reduce leakage errors in mouth-breathing circuits.	AFT3, page 34
Student Pack	Multi-purpose airflow accessories for use with the SS11LA. Used in BSL Lessons.	BSL-ACCPACK, page 34
T-valves, non-rebreathing	Standard and High-performance options. See Lessons BSL18 and H19 (page 43).	AFT21, page 35 • AFT22, page 34
Tubing	Interface the Airflow Transducers (SS11LA and SS52L) with the GASSYS2. See Lessons BSL18, H21, H19 (page 43).	AFT7 and AFT7-L, page 35

## General Airflow Accessories

### Premium Calibration Syringe - AFT26



This 2.0 liter syringe provides a larger volume calibration value and is recommended for exercise physiology applications and higher quality, more consistent

calibrations where greater precision is required. Aluminum casing improves durability. Works with Airflow Transducers or the GASSYS2, and fits AFT7 tubing (35 mm OD coupler).

See Lessons BSL12, BSL13, BSL18, H19, page 43.

### Tubing - AFT7

Smooth bore tubing for use with Airflow Transducers (SS11LA or SS52L) and other large bore accessories such as the AFT25 Facemask with Valve or the AFT21 T-valve for gas analysis studies; included with the GASSYS2. 35 mm ID, 38 mm OD, 1 meter.

AFT7-L: 3 meters.

See Lessons BSL18 and H19, page 43.

### Couplers - AFT11 Series



#### AFT11B

Rigid coupler. 15 mm OD - 22 mm ID.

#### AFT11E

Flexible coupler, included with GASSYS2. Use to connect SS11LA to AFT7 tubing.

Couples 22-25 mm OD to 22-25 mm OD or 35-38 mm ID, or 35 mm OD to 38 mm ID.

#### AFT11F

Flexible coupler. Use two couplers to interface the SS52L Airflow Transducer with the GASSYS2 and other airflow accessories. Couples 35 mm ID to 45 mm OD.

### Gas Cylinder - GASCAL

Calibration Gas Cylinder: 4% CO<sub>2</sub>, 16% O<sub>2</sub>, Bal. N<sub>2</sub> (not for drug use, not for inhalation). Contains 560 Liters @ 2200 psig, 70°F (21°C). Please allow time for drop shipment.



#### Regulator - GASREG

Regulator output connects to GASSYS2. Calibration port via tubing.



# electrodes & recording accessories

## input adapters

Recording cables require multiple leads and electrodes and accept Touchproof connections.



### Shielded Electrode Adapter - SS1LA

Interfaces the MP35 with shielded or unshielded reusable surface and needle electrodes (such as the EL250 and EL450

series). Accepts Touchproof 1.5 mm socket electrode leads. Can be used as an alternative to the SS2L for reusable electrodes.

**Adapter - CBL201:** If you have an older model SS1L, use CBL201 to convert the 2 mm pin connection to the Touchproof 1.5 mm connection.

## High-Impedance Cables



The fully-shielded high-impedance electrode interface cable permits high resolution recording of extracellular signals from small preparations using reusable electrodes (EL250, EL350 or EL450 series, pages 37-38), your own

custom Ag-AgCl electrodes, or glass microelectrodes. The 2-meter cable terminates with standard Touchproof connectors.

**Input Impedance (Common-Mode):** 5e11 Ohms (500 GigaOhm)

**Input Bias Current:** 3 pA Typical

**Voltage Noise:** 1.3  $\mu$ V p-p

**Voltage Noise Density:** 36 nV /  $\sqrt{\text{Hz}}$

**Current Noise Density:** 0.01 pA /  $\sqrt{\text{Hz}}$

**BSLCBL8 Input Range:** MP35  $\pm$ 1 V, MP30  $\pm$ 70 mV

**BSLCBL9 Input Range:** MP35  $\pm$ 10 V, MP30  $\pm$ 700 mV

See Lessons A06, A11, A12, A13, A14, page 45.

## recording electrode leads

### Pinch Leads - LEAD110 Series



Lead110 series electrode pinch leads terminate in standard Touchproof connectors for interfacing to the SS1LA, BSLCBL8, or BSLCBL9. These leads have no ferrous parts. Use shielded leads for mini-

mal noise interference. Generally, for biopotential recordings, one each of LEAD110S-W, LEAD110S-R and LEAD110 are required.

LEAD110 works best as a ground or reference lead.

**LEAD110 Unshielded Electrode Lead - Black (1 m)**

**LEAD110A Unshielded Electrode Lead - Black (3 m)**

**LEAD110S-R Shielded Electrode Lead - Red (1 m)**

**LEAD110S-W Shielded Electrode Lead - White (1 m)**

See these lead options:

- General-purpose Electrode Lead Set (SS2L), page 25
- Multi-lead ECG Cable (SS29L), page 27
- Nerve Chamber, page 38
- Nerve Chamber Leads, page 39
- EDA Snap Lead, see SS57L, page 26
- Cardiac Output Sensor, (SS31L), page 29

## electrode accessories


See also: BSL-ACCPACK (page 34)—includes all consumable items required to run BSL Lessons.

### Radiotranslucent Lead - LEAD108



This 1.8-meter Touchproof lead works with EL508 radiotranslucent electrodes (page 37) for MRI applications.

### Contact Post Lead - LEAD120

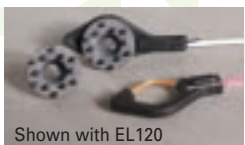

Shown with EL120

This 1-meter lead with Touchproof connector works with the reusable EL120 electrode (page 38). Snap the electrode into place and then plug the lead in via the SS1LA adapter.

**LEAD120-R Red Cable**

**LEAD120-W White Cable**

### Clip Leads - LEAD140 Series

Use these leads to interface with silver wire/platinum recording



electrodes, or to connect directly with an animal preparation. Choose from three styles: electronic test clip, alligator clip



and toothless alligator clip. Leads terminate with a Touchproof connector to interface with the SS1LA shielded electrode adapter, or BSLCBL8 or BSLCBL9.

**LEAD140 Alligator Clip Lead**

**LEAD141 Toothless Alligator Clip Lead**

**LEAD142 Electronic Test Clip Lead (spring-loaded)**



### Hypoallergenic Gel

Conductant for EL250 series reusable electrodes and the EDA (SS3LA) Transducer.

**GEL1 30 g**

**GEL100 250 g**

### Isotonic Gel - GEL101

GEL101 is specially formulated with 0.5% saline in a neutral base and is primarily used as a conductant for the electrodermal activity (GSR) studies. Recommended for research studies only. 125 g.

### Abrasive Pads - ELPAD

Use to abrade the skin to remove non-conductive skin cells for improved signal quality.

### Tape

Attach goniometers and other devices to the body.

**TAPE1** Single-sided adhesive, 9.1 meters

**TAPE2** Double-sided adhesive, 25.6 meters

### Adhesive Disks

Two-sided adhesive collars hold reusable electrodes in place.

**ADD204** For EL254 and EL254S; 4 mm ID; 19 mm OD

**ADD208** For EL258 and EL258S; 8 mm ID; 22 mm OD

## electrodes - disposable Ag-AgCl “snap” series

For the complete range of BIOPAC electrodes, including radiotranslucent MRI-compatible options, visit [www.biopac.com](http://www.biopac.com) or [electrodesales.com](http://electrodesales.com)



Use for added convenience and hygiene; get the same signal transmission as reusable electrodes. Each peel-and-stick electrode is pre-gelled with hypoallergenic gel and intended for one use only. Snap fastener attaches to BIOPAC's SS2L electrode leads.

**SPECIAL**

### General-purpose Electrodes - EL503

Included in every BSL System (pack of 100). Used in over 25 Lessons. These unique electrodes provide excellent signals and are suitable for all biopotential and nerve conduction recordings. 35 mm diameter vinyl tape, 10 mm contact area.

EL503 100/pack

EL503-10 1,000/pack

See the *Student Accessory Pack (BSL-ACCPACK, page 34)* as a packaged alternative for consumables.

### Paired Electrodes - EL500

Use for general-purpose EMG measurements, nerve conduction measurements, and cardiac output. Spacing (center to center): 41 mm x 82 mm x 1.5 mm foam; 25 pairs/pack.

See Lesson H21, page 43.

### Stress Test Electrodes - EL501

Use for short-term recordings where the subject may be in motion or when electrodes should be closely placed, such as multi-channel ECG, EGG, EMG or EOG recordings. 38 mm diameter, mounted on 1.5 mm thick foam with strong adhesive; 50/pack.

### Long-term Electrodes - EL502

Moisture-resistant backing; solid gel adheres well to skin. 41 mm diameter tape, 10 mm contact area.

EL502 50/pack

EL502-10 500/pack

### Cloth Base Electrodes - EL504



Use these compliant, trimmable and flexible electrodes in difficult locations, such as the face for EMG or fingers for nerve conduction. 2.5 cm squares.

EL504 30 pack

EL504-10 300/pack

See Lessons H27, H31, page 43-44.

### Bioimpedance Strip Electrodes - EL506

25 cm (trimmable), Ag-AgCl—replace band electrodes! Use with the noninvasive Cardiac Output Sensor (SS31L, page 29).

EL506 8/pack

EL506-10 80/pack

See Lesson H21, page 43.

### EDA (GSR) Electrodes - EL507

Use with the EDA Lead (SS57L, page 26) for skin conductance and resistance setups. Electrodes have an increased contact area and isotonic wet gel.

EL507 100/pack

EL507-10 1,000/pack

See Lesson BSL9, BSL14, H11, H24, page 44.

### Radiotranslucent Electrodes - EL508



These disposable, radiotranslucent electrodes are MRI-compatible and pre-gelled. Use with LEAD108, page 36.

EL508 100/pack

## electrodes - reusable

### Ag-AgCl Electrodes - EL250 Series



Silver-silver chloride reusable electrodes. Use EL250 series electrodes as an alternative to the SS2L and EL500 series disposable electrodes. Typically, for one biopotential input, you will

need two shielded electrodes for signal inputs and one unshielded electrode for ground. The 1-meter leads terminate in Touchproof 1.5 mm socket connectors for direct connection to the SS1LA, BSLCBL8 or BSLCBL9 lead adapters (page 36). All EL250 Series electrodes require adhesive disks (ADD200 series) and recording gel (GEL1 or any preferred recording gel); see page 36. Available shielded or unshielded. When shielded, the gray lead is for electrode contact and the black lead is for the lead shield.

EL254 4 mm contact, 7.2 mm housing, unshielded Ag-AgCl

EL254S 4 mm contact, 7.2 mm housing, shielded Ag-AgCl

EL258 8 mm contact, 12.5 mm housing, unshielded Ag-AgCl

EL258S 8 mm contact, 12.5 mm housing, shielded Ag-AgCl

See Lesson H27, page 43.



### Bar Electrodes for Animals - EL350 Series

Use on animals to record a signal during nerve conduction, somatosensory or muscle twitch studies. The bar configuration permits easy electrode placement

without disturbing electrode-to-electrode spacing. Two tin electrode disks are placed 30 mm apart in a watertight acrylic bar; non-ferrous. When using bar electrodes for signal recording, a single ground lead (LEAD110 with EL503) is required. The leads (61cm long) terminate in standard Touchproof 1.5 mm socket connectors, which connect to the SS1LA, BSLCBL8 or BSLCBL9 lead adapters.

EL350 Unshielded, concave bar lead electrode

EL350S Shielded, concave bar lead electrode

EL351 Unshielded, convex bar lead electrode

See the *HSTM01* on page 28 for use on human subjects.



## electrodes - reusable

### Needle Electrodes - EL450 Series



Use for recording in animal subjects and tissue preparations. Shipped non-sterile, so pre-sterilization is required. 28-gauge stainless steel needles, with a flexible cable terminating in standard Touchproof connectors, which connect to the SS1LA shielded electrode lead adapter, or BSLCBL8 or BSLCBL9.

#### EL450 - Unipolar, Teflon®-coated

Teflon® coating covers the entire length of the electrode, exposing just the tip for effective isolation of subcutaneous recording points.

Dimensions: 300 µm (dia), 3.7 cm (long), 1.2 mm x 61 mm lead

#### EL451 - Bipolar, Teflon®-coated

Use when recording from a single site, as in studies of single muscle fibers.

Dimensions: 460 µm (dia), 3.0 cm (long), 1.2 mm x 91 mm lead

#### EL452 - Unipolar, Uncoated

Use for small animal preparations.

Dimensions: 300 µm (dia), 1.5 cm (long), 1.2 mm x 61 mm lead



### Contact Posts - EL120

The EL120 Ag-AgCl electrode has contact posts designed to improve contact through fur or hair. Twelve posts create a 10 mm contact area. The posts are 2 mm deep to

push through fur/hair to provide good contact with the skin surface. Requires LEAD120, page 36. Shipped 10/pack.

### Micromanipulator - MANIPULATOR

This manual micromanipulator is a reliable, durable, and economical solution for high-precision experiments. Vernier scales allow readings to 0.1 mm. X-axis fine control allows readings to 10 µm. Includes a tilting base and ships with a standard 12 mm clamp

and electrode holder (14 cm long). All control knobs project to the rear, so units can be tightly grouped. Specify left- or right-handed unit when ordering.

Travel Range:	Resolution:
X-axis Fine	10 mm 0.01 mm
X-axis	35 mm 0.1 mm
Y-axis	25 mm 0.1 mm
Z-axis	25 mm 0.1 mm



## stimulator clip leads

Use these clip leads to interface stimulating electrodes, or to connect directly with an animal preparation. Each 1 meter cable has two clips and terminates with one BNC connector to interface with the BSLSTM or SS58L Stimulator (page 25) and silver or platinum wire electrodes.



BSLCBL7



BSLCBL11



BSLCBL12

BSLCBL7 - BNC to 2x Alligator

BSLCBL11 - BNC to 2x Electronic Test Clip (spring-loaded)

BSLCBL12 - BNC to 2x Toothless Alligator

## stimulation options

### Animal Stim. Needle Electrode - ELSTM2



Use when applying a stimulus to animal and tissue preparations (not for use on humans); connect to the BSLSTM or SS58L stimulators (page 25). The dual stainless steel needles

are 2.5 cm long x 0.3 mm diameter and are Teflon® coated.

Needle electrodes are shipped non-sterile, so pre-sterilization is required. 2.5 meter cable with BNC input.

See Lessons A02, A09, page 45.

See also Human-safe Stimulation Electrode (HSTM01, page 28) and Field stimulation electrode (BSLSTIMHLD, page 33)

### Output Adapter - BSLCBL6

Each stimulator lead connects to the BNC connector on the front panel of the BSLSTM stimulator. BSLCBL6 outputs signals from



the BSLSTM with a cable that terminates in a 3.5 mm phone plug socket with a 1.3 m cable and can be used to connect headphones (OUT100) for auditory evoked response.

### Voltage Drive & Current Monitor Cable - BSLCBL10



Use this cable with an MP35 to perform ion transport experiments. Output from the MP35 drives a set of voltage electrodes and also monitors the current. The maximum recordable current is 50 micro amps. Built-in circuitry allows the MP35 to perform in a voltage clamping mode; the voltage clamp cable terminates in a standard connector. Use with the SS1LA, BSLCBL8 or BSLCBL9 lead adapters to interface with Ag-AgCl electrodes.

## nerve chamber & accessories



The acrylic, desktop Nerve Chambers incorporate 15 stainless steel electrodes for recording and stimulating a variety of different nerve preparations. Each stainless steel electrode is spaced 5mm apart to provide a variety of recording and stimulating configurations. Chambers have a 35 ml reservoir and include a drain with valve and hose. Interface with the BSLSTM or SS58L stimulator via BSLCBL2A and BSLCBL4B nerve conduction cables. The NERVE1 chamber includes a convenient agent well for adding compounds (ether or dry ice, etc.) and a lid to enclose the preparation when protocol requires it.

The NERVE1 chamber includes a convenient agent well for adding compounds (ether or dry ice, etc.) and a lid to enclose the preparation when protocol requires it.


Dimensions: (cm): 4.5 (h) x 7 (w) x 14 (l)

NERVE1 - with agent well and lid

NERVE2 - standard chamber

See Lesson A03, page 45.

## Nerve Chamber Recording

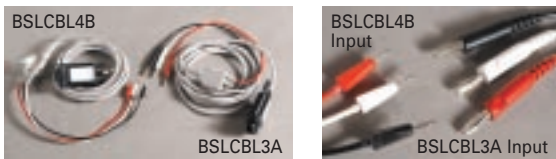

**BSLCBL3A** - banana plugs x 3; for third-party nerve chambers  
**BSLCBL4B** - 2 mm pin plugs x 3; for BIOPAC NERVE1 & 2

### BSLCBL3A and BSLCBL4B Specifications

Gain: 1/10 (divide by 10)  
 Input Impedance (Common-Mode): 500 GigaOhm  
 Common-Mode Rejection: 90 dB  
 Input Bias Current: 3 pA  
 Noise Voltage: 1.3  $\mu$ V p-p  
 Cable: 1.2 meters

## Nerve Chamber Stimulation



**BSLCBL1** - BNC to 2 x banana.  
 Use to connect the BSLSTM or SS58L stimulator to third-party nerve chambers; 1.2-meter cable.

**BSLCBL2A** - BNC to 2 x 2 mm banana pin with ground jumper. Use

to connect the BSLSTM or SS58L stimulator to BIOPAC nerve chambers; 1.2-meter cable.

## interface cables

### Analog Input

#### BNC Input

Connect to the analog input ports on the front panel of the MP35 or MP30 ("CH1" – "CH4") to record signals from other devices, such as amplifiers, third-party chart recorders, and signal generators.

SS70L Isolated Female BNC ( $\pm 10$  V max.; attenuation x 10)

SS9L Unisolated Male BNC ( $\pm 50$  V max.; attenuation x 1,000)

**Safety Note:** Never connect the SS9L to an MP35 or MP30 unit if electrodes from other channels are attached to human subjects—this may void the isolation.

### 3.5 mm phone plug adapter - BSLCBL5

Use this 3.5 mm phone plug adapter to interface the MP35 or MP30 with equipment that outputs high-level voltage signals. The built-in attenuation of 1/200 translates 10 V to 50 mV. Included with the TSD122.



### Input Adapter for Research Amplifiers - BSLCBL14

Use this 3.5 mm phone plug cable to interface with BIOPAC research amplifiers via the IPS100C Isolated Power Supply. Contact BIOPAC for details.



### Touchproof "Y" Adapter - CBL204

This "Y" electrode lead adapter (25 cm long) provides two Touchproof sockets and one Touchproof plug. Use to connect multiple electrode sites (such as Vin- and

GND from BSLCBL8/9, page 36) to a glass microelectrode. Connect multiple CBL204s to reference three or more electrode leads to the same input or output.

## Analog Out Cables

### BNC Output - OUT2

Analog output cable connects to the Analog Out port on the rear of the MP35. DSUB 9m connector. The cable terminates in a male BNC connector.

\*See OUT1 Headphones on page 26.

### Custom Output Interface - SS-KIT-OUT



The kit comes with a Simple Sensor connector, cable, and components to properly interface with speakers, stimulators and other equipment requiring a trigger or analog output signal. See the BSL-TCI series for existing transducer interfaces.

## transducer interface options

### Transducer Connector Interfaces - BSL-TCI SERIES



Save money and interface with your existing transducers. BSL-TCI Series connectors for common transducer manufacturers are listed below—if you don't see the part you need, call BIOPAC Support or use the SS-KIT to build a custom interface.

Part #	Connector Make	Works with Transducers from...
<b>BSL-TCI0</b>	6-pin	Grass and Gilson
<b>BSL-TCI1</b>	5-pin	Beckman
<b>BSL-TCI2</b>	8-pin	AD Instruments, iWorx, and WPI Transducers
<b>BSL-TCI3</b>	9-pin hexagonal	Lafayette and Narco
<b>BSL-TCI4</b>	6-pin	Honeywell
<b>BSL-TCI5</b>	4-pin phone jack	BIOPAC NIBP100A BP module
<b>BSL-TCI6</b>	12-pin	Beckman
<b>BSL-TCI7</b>	5-pin	Nihon Koden
<b>BSL-TCI8</b>	7-pin	Narco
<b>BSL-TCI9</b>	8-pin	Fukada
<b>BSL-TCI10</b>	12-pin	Gould
<b>BSL-TCI11</b>	6-pin male	Hugo Sachs and Harvard Apparatus
<b>BSL-TCI12</b>	5-pin, 240 degrees	Thornton
<b>BSL-TCI13</b>	Piezo (BNC)	iWorx PT100
<b>BSL-TCI14</b>	1/4" phono socket	Lafayette Force
<b>BSL-TCI15</b>	5-pin DIN	Vernier
<b>BSL-TCI16</b>	BT Connector	BIOPAC RXPROBE02 and Vernier
<b>BSL-TCI17</b>	5-pin DIN	Intelitool
<b>BSL-TCI18</b>	2 x 2 mm sockets	Mercury Strain Gauge or Indium Gallium
<b>BSL-TCI19</b>	6-pin mini DIN	Intelitool
<b>BSL-TCI20</b>	3.5 mm phono jack	Intelitool
<b>BSL-TCI21</b>	BNC	BIOPAC RXPROBE01 and 3rd-party pH probes
<b>BSL-TCI22</b>	Breadboard/Electrode	BIOPAC Breadboard SS39L to LEAD SS2L

Note: All company names listed above are ®.

### Custom Input Interface - SS-KIT-IN

Adapt third-party transducers to the MP35. The kit comes with a Simple Sensor connector, cable, and components to properly interface with quarter, half or full bridge transducers (pressure, force, strain, acceleration, sound, etc.). See the BSL-TCI series for existing transducer interfaces.







## USB Cable - CBLUSB

2.5-meter replacement USB cable connects the MP35 to a USB Port. Includes and provides EMI protection to maintain BSL System certified safety rating (CE, EMC).

## power transformers



### IEC 60601-1 Certified - AC300A

+12V, 1.25 amp—for MP35 to mains wall outlet. Included with each BSL System. Specify power cord: ACCORD-HUS (Hospital grade, USA) or ACCORD-EURO (Europe).

### +12V, 1.00 amp - AC100A

For GASSYS2 or MP30 to mains wall outlet. Specify power cord: ACCORD-US (USA) or ACCORD-EURO (Europe).

### +6V, 1.50 amp - AC137A

For heating elements to mains wall outlet. Powers the heating element for any of the SS45L-SS52L pneumotachs (or RX137 replacement heads). Specify power cord: ACCORD-US (USA) or ACCORD-EURO (Europe).

## BSL Laboratory Manual - MANBSL3S

BIOPAC lessons provide online instructions and real-time prompts to guide students through setup, calibration and recording. This Laboratory Manual provides further details for recording (such as hardware and subject setup), background material and a student Lab Report with guided analysis for BSL Lessons 1-17. One manual is included with each system—order additional copies through your bookstore for students. Please specify a language when ordering.

MANBSL3S - English

MANBSL3S-ES - Spanish

MANBSL3S-FR - French

MANBSL3S-J - Japanese

\*See also published Lab Manuals, page 8.

## Lessons Upgrade - BSLUPG16&17

This lesson upgrade package includes a Blood Pressure Cuff Transducer (SS19L, page 27) and an Electronic Stethoscope Transducer (SS30L, page 28) to run BSL Lessons 16 & 17.

## Battery Pack - BAT100



For portability, use the BAT100 rechargeable battery pack and charger with the BSL System. The maintenance-free battery pack is built into a carrying case with a shoulder strap. The fully-charged battery will operate a

BSL System for a minimum of 24 hours. Includes battery pack, universal recharger and all necessary cables. Specify power cord: ACCORD-US (USA) or ACCORD-EURO (Europe).

Output Capacity:	12 V @ 13 amp-hours
Operating Time for MP35 with 4 transducers:	24-36 hours nominal
Charge Time:	15 hours
Dimensions:	22 cm (high) x 8 cm (wide) x 24 cm (deep)
Weight:	5.6 kg
Recharger:	
Output Capacity:	12 V @ 1.0 amps
Input:	120/240 VAC @ 50/60 Hz
Dimensions:	8 cm (high) x 13 cm (wide) x 15 cm (long)
Weight:	1.8 kg

## Platform Change MP30 Platform Change

If you have an existing MP30 System that interfaces with your computer via an ISA100B card, use a Platform Change to update your system. Contact a Biopac Student Lab Specialist or visit the BIOPAC web site ([www.biopac.com](http://www.biopac.com)) for details and upgrade options for your computer, hardware, and operating system.

PLAT30W1 for Windows

PLAT30M1 for Macintosh

*Packages, specifications and lessons subject to change without notice.*

# core packages

BIOPAC offers a comprehensive line of electrodes, transducers and accessories that work with the MP35 System for a wide variety of applications. Core packages are just one ordering option, based on typical systems sold to users from a particular discipline.

- The BASIC System is the foundation for all packages.
- The BASIC, ADVANCED and ULTIMATE Systems provide a general system for cross-disciplinary users.
- The BASIC, ADVANCED and ULTIMATE Systems have been offered since the Biopac Student Lab was launched in 1995, and are now in use in thousands of labs worldwide.
- The new CORE packages are offered to meet the diverging requirements of specific departments.
- Any system can be used as a starting point to develop your own custom package.

- Review the Core Package Components table to determine which package meets your requirements, or talk to a Biopac Student Lab Specialist to create your own custom package.
- Packages contain everything—including consumable items—necessary to run a specified number of lessons, specific to that particular discipline.

Each package is discounted to reflect the number of transducers and accessories included.

If you have existing transducers that you'd like to continue using, you can select from more than 20 interface options (page 39) for manufacturers like Grass, Gould, Gilson, Harvard and Vernier.

BIOPAC is continually adding and improving products and lessons. Visit [www.biopac.com](http://www.biopac.com) for the latest offerings.

# core packages

- Packages are offered as a comprehensive building block for discipline-specific applications.
- Increase application potential by adding more hardware.

## Core Package Components

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
				Basic	Advanced	Ultimate	Human	Animal	Biology	Exercise Phys.	Psychophys.	Biomed. Eng.	Pharm. & Tox.
<b>BIOPAC Hardware</b>	<b>Part #</b>	<b>Page</b>											
MP35 Data Acquisition Unit & Cables	ships with system	24	x	x	x	x	x	x	x	x	x	x	x
BSL Software—Lessons & PRO	ships with system	2-7	x	x	x	x	x	x	x	x	x	x	x
Lab Manual—17 BSL lessons	MANBSL3S	40	x	x	x	x	x	x	x	x	x	x	x
PRO Lessons—40+ experiments	on web	43-45	x	x	x	x	x	x	x	x	x	x	x
Electrode Lead (x2)	SS2L	25	x	x	x	x			x	x	x	x	
High-impedance Cable ±1V Input (x2)	BSLCBL8	36						x					x
High-impedance Cable ±5V Input	BSLCBL9	36							x				
Electrodes, Disposable (100/pk)	EL503	37	x	x	x	x	x	x	x	x	x	x	x
Abrasive Pads (10/pk)	ELPAD	36	x	x	x	x	x	x	x	x	x	x	x
Airflow Transducer	SS11LA	27		x	x	x			x	x		x	x
Calibration Syringe—600ml	AFT6	34		x	x	x			x	x		x	x
Bacterial Filters, Disposable (10/pk)	AFT1	34		x	x	x			x	x		x	x
Mouthpieces, Disposable (10/pk)	AFT2	34		x	x	x			x	x		x	x
Nose Clips, Disposable (10/pk)	AFT3	34		x	x	x			x	x		x	x
Electrodermal Activity (GSR) Lead	SS57L	26		x	x				x		x		
EDA (GSR) Electrodes (100/pk)	EL507	37		x	x				x		x		
Hand Switch—Pushbutton	SS10L	26		x	x				x	x	x	x	
Headphones	OUT1	26		x	x				x	x	x	x	
Pulse Plethysmograph Transducer	SS4LA	26		x	x				x	x	x	x	
Respiratory Effort Transducer	SS5LB	26		x	x				x	x	x		
Temperature Transducer	SS6L	26		x	x				x	x			
Colored Paper—for L09 Polygraph	PAPER1	-		x	x				x		x		
Electrode Gel	GEL1	36			x				x				
Tape—Single-sided	TAPE1	36		x	x				x	x			
Blood Pressure Cuff Transducer	SS19L	27			x	x			x	x	x	x	
Electronic Stethoscope Transducer	SS30L	28			x	x			x	x	x	x	
Force Transducer— Variable range	SS12LA	27			x				x				
Hand Dynamometer	SS25LA	27			x	x			x	x		x	
Multi-Lead ECG Cable	SS29L	27			x				x				
Pressure Transducer	SS13L	27			x				x				x
Stimulator—BSL	BSLSTM	25			x				x				x
Stim. Electrode—human-safe	HSTM01	28			x				x				
Stimulator—Low Voltage	SS58L	25						x					
Stim. Electrode—animal	ELSTM2	38						x	x				x
Dissolved O <sub>2</sub> Probe	SS69L	31											
Dissolved O <sub>2</sub> Interface	BSL-TCI16	39						x	x				
Signal Processing Breadboard	SS39L	30										x	
Electrode Lead (x2)	SS1LA	36											
Electrode Lead—unshielded (x2)	LEAD110	36						x					x
Electrode Leads—shielded (x2)	LEAD110S-W & -R	36						x					x
Force Transducer—200g	SS65L	31						x					x
Needle Electrodes (x3)	EL452	38						x	x				x
Nerve Chamber	NERVE2	38						x	x				x
Nerve Cable—Recording	BSLCBL4B	39						x	x				x
Nerve Cable—Stimulator	BSLCBL2A	39						x	x				x
pH Interface (SS68L or other pH probe)	BSL-TCI21	39							x				
				Basic	Advanced	Ultimate	Human	Animal	Biology	Exercise Phys.	Psychophys.	Biomed. Eng.	Pharm. & Tox.
				Cross-discipline			10	12	14	16	18	20	22

### Hardware Options

- This table only lists parts included in CORE PACKAGES.
- For a full list of hardware options, see the BSL Hardware section (pages 24-40) or check the Index (page 47).
- All parts can be ordered individually—you can add hardware options, or create your own package to suit your curriculum.

Visit [www.biopac.com](http://www.biopac.com) or contact a Biopac Student Lab Specialist to discuss your specific application needs.

See the page indicated for an overview of each discipline, including lessons the Core Package supports and suggested hardware options and potential applications.





# transducer accessory “TA” packs

The following “TA” Packs are suggested transducer accessories for any core package to increase the application potential of your lab. Each “TA” Pack allows you to perform the additional lessons listed—or you can create your own experiments. All items can be ordered individually.

- See the Core Packages summary (p41).
- For a full list of hardware options, see the BSL Hardware section (page 23-40).
- For an overview of each lesson, see the Lessons Summary (page 43-46).

---

## Human Physiology TA — BSLHPY-TA

### *This hardware*

Stimulator BSLSTM, page 25  
Stimulation Electrode for Humans HSTM01, page 28  
Finger Twitch Transducer SS61L, page 30  
Reflex Hammer SS36L, page 30

*See Core Human Phys. on page 11*

### *Adds these lessons*

H03 Nerve Conduction Velocity, page 44  
H06 Finger Twitch, page 43  
H28 Reflex Response (patellar tendon), page 44

---

## Animal Physiology TA — BSLAPH-TA

### *This hardware*

Cardiac Output Sensor SS31L, page 29  
Voltage Drive & Current Monitor Cable BSLCBL10, page 38  
Bioimpedance Strip Electrodes EL506-10 (80/pk), page 37

*See Core Animal Phys. on page 13*

### *Adds these lessons*

H21 Impedance Cardiography (CO), page 43  
A10 Regulation of Circ. & Resp., page 45  
A13 Active Transport (frog skin), page 45

---

## Exercise Physiology TA — BSLEXP-TA

### *This hardware*

Multi-lead ECG Cable (12-lead) SS29L, page 27  
Cardiac Output Sensor SS31L, page 29  
CO<sub>2</sub> & O<sub>2</sub> Analysis Module GAS-SYSTEM2, page 34  
Bioimpedance Strip Electrodes EL506-10 (80/pk), page 37

*See Core Exercise Phys. on page 17*

### *Adds these lessons*

BSL18 Metabolic Rate, page 43  
H01 12-lead ECG, page 43  
H19 VO<sub>2</sub> & RER, page 43  
H21 Impedance Cardiography (CO), page 43

---

## Biomechanics TA — BSLEXM-TA

### *This hardware*

Reflex Hammer Transducer SS36L, page 30  
Goniometer (Twin axis) SS21L, page 29  
Stimulator BSLSTM, page 25  
Stimulation Electrode for Humans HSTM01, page 28  
Heel/Toe Strike Transducer SS28L, page 29

*See Core Exercise Phys. on page 17*

### *Adds these lessons*

H03 Nerve Conduction Velocity, page 44  
H17 Biomechanics (Goniometry & EMG), page 44  
H28 Reflex Response (patellar tendon), page 44  
H35 Range of Motion: Sit & Reach, page 45

---

## Psychophysiology TA — BSLPSY-TA

### *This hardware*

Cardiac Output Sensor SS31L, page 29  
SuperLab Stimulus Presentation Pkg. STP35W, page 32  
Bioimpedance Strip Electrodes EL506-10 (80/pk), page 37

*See Core Psychophys. on page 19*

### *Adds these lessons*

H21 Impedance Cardiography (CO), page 43  
H30 Stroop Effect, page 44  
H31 Prepulse Inhibition, page 44

---

## Biomedical Engineering TA — BSLBME-TA

### *This hardware*

Cardiac Output Sensor SS31L, page 29  
Stimulator BSLSTM, page 25  
Stimulation Electrode for Humans HSTM01, page 28  
Voltage Drive & Current Monitor Cable BSLCBL10, page 38  
Bioimpedance Strip Electrodes EL506-10 (80/pk), page 37

*See Core Biomed. on page 21*

### *Adds these lessons*

H03 Nerve Conduction Velocity, page 44  
H21 Impedance Cardiography (CO), page 43  
A13 Active Transport (frog skin), page 45

---

## Pharmacology & Toxicology TA — BSLPHA-TA

### *This hardware*

Tissue Bath Station ITBS100, page 33

*See Core Pharm. & Tox. on page 23*

### *Adds this lesson*

A05 Visceral Smooth Muscle, page 45

---

## CARDIOVASCULAR

### BSL LESSON 5: ECG I

Record Lead II ECG and examine components of the ECG complex as an introduction to the electrocardiograph and the recording of the heart's electrical signal.

### BSL LESSON 6: ECG II

Record ECG using bipolar Leads I and III; the software calculates Lead II to demonstrate Einthoven's law.

### BSL LESSON 7: ECG & PULSE

Use a pulse plethysmogram transducer and Lead II ECG to examine the mechanical action of the heart and peripheral pulse pressure to learn how the heart pumps blood throughout the body.

### BSL LESSON 16: SYSTEMIC BLOOD PRESSURE

Record arterial blood pressure using the auscultatory (cuff) technique, Korotkoff sounds using an amplified stethoscope, and ECG using Lead II.

### BSL LESSON 17: HEART SOUNDS

Record ECG Lead II and place an amplified stethoscope at four different locations to listen to the sounds of the heart's valves and correlate the sounds with the cardiac cycle.

### H01 12-LEAD ECG

Record a 12-lead ECG and observe changes in the frontal plane vectors throughout a cardiac cycle.

### H08 ECG DIVE REFLEX—ACTIVE LEARNING

Subjects immerse their face in cold water and record the change in heart rate that occurs to investigate the physiological reason for the observed response.

### H21 IMPEDANCE CARDIOGRAPHY


Noninvasively record and measure stroke volume and heart rate data and correlate with cardiac output.

### H23 SIGNAL AVERAGED ECG

Record ECG data under different experimental conditions and perform a Signal Averaged ECG recording for each segment of data.

## MUSCULAR

### BSL LESSON 1: EMG I

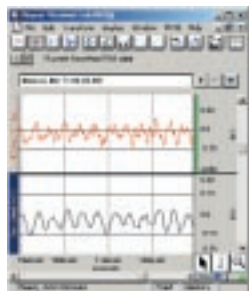

Record maximum grip clench for the dominant and non-dominant hand to investigate the properties of skeletal muscle.

### BSL LESSON 2: EMG II

Use a hand dynamometer to record maximum grip strength for both hands and explore the role of skeletal muscle in performing mechanical tasks.

### H06 FINGER TWITCH—HUMAN

Record the force generated from a finger twitch and measure the stimulus frequency required to induce fatigue. (Alternative to Frog Gastroc.)

## H07 EMG—ACTIVE LEARNING


Investigate the electrical activity of different muscles as they contract with varying degrees of force and design experiments by selecting muscles to record from and creating activities those muscles will perform.

### H27 FACIAL EMG

Record EMG response on the corrugator supercilii & zygomaticus major muscles.

### H34 ELECTROGASTROGRAM

Record electrical activity through stomach muscles (EGG) and note the power and frequency of contractions at rest and after eating.

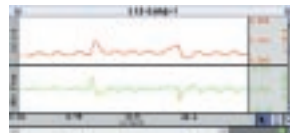

## RESPIRATORY & PULMONARY FUNCTION

### BSL LESSON 8: RESPIRATORY CYCLE I

Record chest contraction & expansion and ventilation, then correlate respiration changes with ventilation to examine the effects of cerebral influence and chemoreceptor influence on the medullary control centers.

### BSL LESSON 12: PULMONARY FUNCTION I

Perform a variety of pulmonary measurements: Tidal volume, Inspiratory capacity, Expiratory capacity, Functional residual capacity, Vital capacity and Total lung capacity.

### BSL LESSON 13: PULMONARY FUNCTION II

Record and analyze Forced Vital Capacity, Forced Expiratory Volume ( $FEV_{1,2,3}$ ) and Maximal Voluntary Ventilation (MVV) to build on the principles established in Lesson 12.

### BSL LESSON 18: METABOLIC RATE

Record indirect basal metabolic rate (BMR) and post-exercise metabolic rate.

### H19 $VO_2$ & RER

Record and measure oxygen consumption (absolute  $VO_2$ ) and respiratory exchange ratio (RER) under a variety of conditions and observe the relationship between  $VO_2$  and RER.

## NEUROPHYSIOLOGY

### BSL LESSON 3: EEG I

Record EEG from the occipital lobe while performing a variety of tasks to demonstrate how the brain's electrical activity varies dependent upon the task being performed.

### BSL LESSON 4: EEG II

Discover how the brain constantly receives sensory input and integrates the information before processing it. The system records and displays raw EEG, alpha wave and alpha-RMS activity.

### Lesson Update

See [www.biopac.com](http://www.biopac.com) for lesson additions and product updates.

## BSL LESSON 9: GSR & POLYGRAPH (EDA)

Record changes in respiratory rate, heart rate and electrodermal activity (skin conductance) to become familiar with the standard physiological measures recorded by a polygraph and study the effects of cognitive behavior and emotion.

## BSL LESSON 10: ELECTROOCULOGRAM (EOG) I

Record horizontal and vertical eye movement to demonstrate eye fixation and tracking. Students perform a number of tasks that allow them to record the duration of saccades and fixation.

## BSL LESSON 11: REACTION TIME I

Subject hears two schedules of clicks through headphones and reacts by pressing a pushbutton hand switch as quickly as possible to demonstrate the effect of learning and physiological processes on reaction times.

## BSL LESSON 14: BIOFEEDBACK

Record ECG, heart rate and electrodermal activity, and try to influence heart rate and EDA (GSR) to control the position of a bar graph to demonstrate the principles of biofeedback training for relaxation purposes.

## H03 NERVE CONDUCTION VELOCITY

Record responses along the ulnar nerve of a human subject to observe the Threshold, Maximal and Supra-Maximal response levels and determine nerve conduction velocity along the ulnar nerve.

## H09 AUDITORY EVOKED POTENTIAL (AEP)

Present an auditory stimulus to a human subject and record Auditory Evoked Potential.

## H10 HEMISPHERIC EEG

Record EEG and study effects of sensory stimulation or change in attitude/attention on alpha rhythm, beta rhythm, and hemispheric asymmetry.

## H11 MIRROR TEST: SENSORY MOTOR LEARNING & EDA

Correlate efficiency in a task requiring movement and attention focus with reticular tone (which indirectly indicates emotional fluctuations) and analyze performance over repeated trials.

## H12 SACCADES: EOG


Explore applications of electrooculography and observe the constant saccade durations for a variety of given angular displacements.

## H13 TRACKING: EOG

Observe tracking movements used while watching a moving object and demonstrate the difference between eye movement based on actual visual stimulation and imagined recreations.

## H14 FIXATION I: EOG

Record horizontal EOG and observe Ocular Fixation while reading.

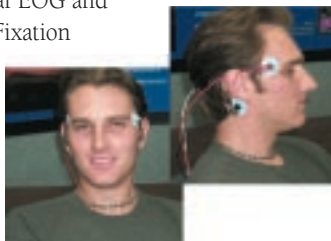

## H15 FIXATION II: EOG

Record a horizontal and vertical EOG and observe spontaneous gaze changes produced when viewing an image, and then correlate results from the plot with the subject's attitude or level of interest.

## H16 REFLEXES & REACTION TIME

Measure basic reflex and reaction time (visual stimulus) exercises and record reaction time to auditory stimulus. Compare reaction times from fixed interval and pseudo-random presentation to study the effects of learning and physiological processes on reaction times.

## H22 VISUAL EVOKED RESPONSE

Present a visual stimulus to a human subject and record Visual Evoked Potentials (P100 test).

## H24 HABITUATION

Record EDA (GSR) and Heart Rate response to repeated stimulus to demonstrate habituation and its probabilistic trend toward decreased response.

## H28 REFLEX RESPONSE

Record knee and ankle reflex response with the SS36L Reflex Hammer transducer. Option: Use the SS20L Goniometer to measure angular movement in response to varying strike force.

## H30 STROOP EFFECT

Record strength of interference between two associative tasks: naming and reading.

## H31 PREPULSE INHIBITION

Record the startle response with and without a prepulse inhibition stimulus.

## H32 HEART RATE VARIABILITY

Explore statistical measures, geometric measures, and spectral analysis in heart rate variability.

## EXERCISE PHYSIOLOGY

### BSL LESSON 15: AEROBIC EXERCISE PHYSIOLOGY

Record ECG, heart rate, airflow and skin temperature under a variety of conditions to see how the body responds to changing metabolic demands.

## H04 BLOOD PRESSURE

Record (noninvasive) blood pressure response to isometric or straining exercise.

## H05 WINGATE TEST

Record the Wingate Anaerobic Test (on a modified, plate loaded, Monark 818E work ergometer) and complete WAnT calculations.

## H17 BIOMECHANICS (Goniometry & EMG)

Record muscle activity from the triceps and biceps while recording angle of limb movements.

## H18 EXERCISE PHYSIOLOGY (Blood Pressure)

Record Automatic Noninvasive Blood Pressure in pre- and post-exercise conditions and observe the change between the conditions.



### H35 RANGE OF MOTION: SIT & REACH

Students use a goniometer to record angle of joint movement (i.e., hip, ankle etc.) during a Sit & Reach test.

### H36 MUSCULAR BIOFEEDBACK

Students record EMG and use auditory and visual (bar graph) biofeedback and touch to increase muscle performance.

## BME - BIOMEDICAL ENGINEERING

### H02 COMPARTMENTAL MODELING

Explore Westheimer's saccadic eye movement model, which represents the eye as a 2nd order system. Record eye motion via EOG setup and compare to modeled results.

### H20 FILTERING

Design and develop software-based digital filters to perform a variety of physiological signal filtering tasks. Cascade 2nd order biquads to create high order filters.

### H25 SIGNAL PROCESSING BREADBOARD I (8 Circuits)

Schematic and design notes for Square Wave Oscillator, Instrumentation Amplifier, High Pass Active Filter, Active Gain Block and Low Pass Filter, Notch Filter for 60 Hz Rejection, QRS Detection: Band Pass Filter, QRS Detection: Absolute Value Circuit; QRS Detection: Low Pass Filter.


### H26 SIGNAL PROCESSING BREADBOARD II (System)

Block diagram to build ECG Signal Processor with SS39L.

### H33 FFT FAST FOURIER TRANSFORM

Build up a square wave from cosine components and use the FFT function to analyze the composite response.

## ANIMAL

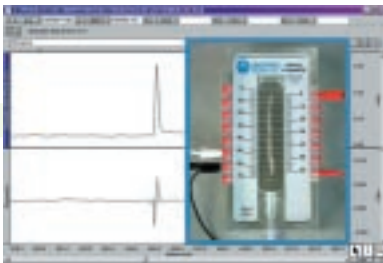

### A01 FROG PITH & PREPARATION

Explanation of how to pith and prepare a frog for experiments A02, A03 and A04.

### A02 FROG GASTROCNEMIUS

Directly stimulate the frog gastrocnemius muscle (or stimulate the muscle via the sciatic nerve) and record threshold voltage and contractile responses.

### A03 FROG SCIATIC NERVE

Record compound action potentials of the dissected sciatic (somatic motor and sensory) nerve.

### A04 FROG HEART

Record cardiac rate and contractile responses of the surgically exposed frog heart. Option: Study the effects of chronotropic and inotropic agents on the heart.

### A05 VISCERAL SMOOTH MUSCLE

Study the effects of media ionic composition, temperature, and various pharmacological agents on the contraction of the visceral smooth muscle of the rabbit ileum.

### A06 COCKROACH VENTRAL NERVE

Record nerve activity from the ventral nerve cord while stimulating the cerci with puffs of air.

### A07 Q<sub>10</sub> PRINCIPLE (Dissolved O<sub>2</sub>—Goldfish)

Demonstrate the Q<sub>10</sub> principle by measuring the metabolic rate of goldfish at two temperatures: 22° C (acclimation temperature) and 32° C (acute exposure temperature).

### A08 ACTION POTENTIALS IN EARTHWORM

Use extracellular recording techniques to stimulate and record action potentials from an earthworm's nerve cord. Measure conduction velocity and refractory period, and plot a strength versus duration curve.

### A09 PROPERTIES OF CARDIAC MUSCLE (Turtle Heart)

Measure the duration of systole and diastole and observe the effects of diastolic loading. Monitor the effect of vagal stimulation, temperature changes and spontaneous rhythmicity of the heart.

### A10 REGULATION OF CIRCULATION & RESPIRATION

Record aortic and femoral blood pressure, ECG and cardiac output from an anesthetized pig (or dog) to study homeostasis via cardiovascular physiology and short term regulation of blood pressure.

### A11 RESTING POTENTIAL (Crayfish Muscle)

Follow BSL setup and use "Lab 4: Crayfish Muscle Resting Potential" from the Crawdad CD-ROM Lab Manual for Neurophysiology (ISBN 0-87893-947-4) to record and alter resting potential by changing external ion concentration.

### A12 MEMBRANE POTENTIAL (Muscle)

Use a glass microelectrode to record the membrane potential from the frog sartorius muscle. Record the potential change when the electrode penetrates a muscle cell.

### A13 ACTIVE TRANSPORT


Use skin from a frog in an Ussing chamber to record active transport. Measure the skin potential and use Ohm's law to calculate the NaCl transport across the skin.

### A14 CENTRAL PATTERN GENERATORS

Perform extracellular recording on tobacco hornworm pupae to study central pattern generators (CPGs) and neural mechanisms.

### A15 EARTHWORM SMOOTH MUSCLE

Setup earthworm gut with a force transducer and tissue bath to measure contractions and the effect of drugs.

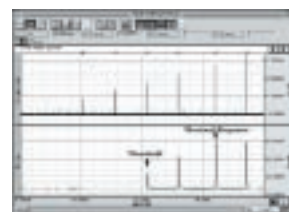

# Lesson hardware guide

## using Core Package hardware

	Page	Basic	Adv.	Ult.	Human	Animal	Biology	Psych	Ex. Phys.	BME	Pharm.	Hardware Used (optional)
BSL1: EMG I	43	x	x	x	x	x	x	x	x	x	x	SS2L p25, (OUT1 p26)
BSL2: EMG II	43			x	x		x		x	x		SS2L p25, SS25LA p27, (OUT1 p26)
BSL3: EEG I	43	x	x	x	x	x	x	x	x	x	x	SS2L p25
BSL4: EEG II	43	x	x	x	x	x	x	x	x	x	x	SS2L p25
BSL5: ECG I	43	x	x	x	x	x	x	x	x	x	x	SS2L p25
BSL6: ECG II	43	x	x	x	x	x	x	x	x	x	x	SS2L x2 p25
BSL7: ECG & Pulse	43		x	x			x	x	x	x		SS2L p25, SS4LA p26
BSL8: Respiratory Cycle I	43		x	x			x		x			SS5LB and SS6L p26
BSL9: GSR & Polygraph	44		x	x			x	x				SS2L p25, SS5LB and SS57L p26
BSL10: Electrooculogram (EOG) I	44	x	x	x	x	x	x	x	x	x	x	SS2L x2 p25
BSL11: Reaction Time I	44		x	x			x	x	x	x		SS10L and OUT1 p26
BSL12: Pulmonary Function I	43		x	x	x		x		x	x	x	SS11LA p27, AFT6 p34
BSL13: Pulmonary Function II	43		x	x	x		x		x	x	x	SS11LA p27, AFT6 p34
BSL14: Biofeedback	44		x	x			x	x				SS2L p25, SS57L p26
BSL15: Aerobic Exercise Physiology	44		x	x			x		x			SS2L p25, SS6L p26, SS11LA p27
BSL16: Systemic Blood Pressure	43			x	x		x	x	x	x		SS2L p25, SS19L p27, SS30L p28
BSL17: Heart Sounds	43			x	x		x	x	x	x		SS2L p25, SS30L p28
BSL18: Metabolic Rate	43											SS11LA p27, GASSYS2 and AFT6 p34
H01 12-Lead ECG	45			x			x					SS29L p27
H02 BME Compartmental Modeling	44		x	x			x	x	x	x		SS2L p25
H03 Nerve Conduction Velocity	44			x			x					SS2L p25, BSLSTM p25, HSTM01 p28
H04 Blood Pressure Response to Straining Exercise	44			x	x		x		x			SS19L p27, SS30L p28
H05 Wingate Test (WAnT)	43		x	x			x		x			SS4LA p26
H06 Finger Twitch	43			x			x					SS61L p30, BSLSTM p25, HSTM01 p28
H07 EMG - Active Learning	43	x	x	x	x	x	x	x	x	x	x	SS2L p25
H08 ECG Dive Reflex - Active Learning	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H09 Auditory Evoked Potential (AEP)	44											BSLSTM p25, OUT101 p31, BSLCBL6 p38
H10 Hemispheric EEG	44	x	x	x	x	x	x	x	x	x	x	SS2L x2 p25
H11 Mirror Test: Sensory Motor Learning & EDA	44		x	x			x	x				SS10L and SS57L p26
H12 Saccades: EOG	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H13 Tracking: EOG	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H14 Fixation I: EOG (3-lead)	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H15 Fixation II: EOG (6-lead)	44	x	x	x	x	x	x	x	x	x	x	SS2L x2 p25
H16 Reflexes & Reaction Time - Active Learning	44		x	x			x	x	x	x		SS10L and OUT1 p26
H17 Biomechanics	44											SS2L x2 p25, SS21L p29
H18 Exercise Phys. - Continuous Noninvasive BP	43											NIBP100A p32
H19 VO <sub>2</sub> & RER	45											SS11LA p27, GASSYS2 and AFT6 p34
H20 BME Filtering	43	x	x	x	x	x	x	x	x	x	x	no transducers required
H21 Impedance Cardiography (Cardiac Output)	44											SS2L p25, SS30L p28, SS31L p29
H22 Visual Evoked Response (VER)	43											SS2L p25, TSD122 p32
H23 Signal Averaged ECG	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H24 Habituation	45		x	x			x	x				SS2L p25, SS10L and SS57L p26
H25 BME Signal Processing I (8 circuits)	45									x		SS39L and (SS60L) p30
H26 BME Signal Processing II (ECG R-Wave Detector)	43									x		SS39L and (SS60L) p30
H27 Facial EMG	44	x	x	x	x	x	x	x	x	x	x	SS2L x2 p25, (SS10L p26)
H28 Reflex Response (patellar tendon)	43											SS2L p25, SS36L p30, (SS20L p29)
H30 Stroop Effect	44											SS10L p26, STP35W p32
H31 Prepulse Inhibition	44											SS2L p25, STP35W p32, (OUT100 web)
H32 Heart Rate Variability Analysis	44	x	x	x	x	x	x	x	x	x	x	SS2L p25
H33 BME FFT Fast Fourier Transform	45	x	x	x	x	x	x	x	x	x	x	no transducers required
H34 Electrogastrogram from human (stomach)	43	x	x	x	x	x	x	x	x	x	x	SS2L p25
H35 Range of Motion: Sit & Reach	45											SS21L p29
H36 Muscular Biofeedback (auditory, visual, touch)	45		x	x			x	x	x	x		SS2L p25, OUT1 p26
A01 Frog Pith & Prep	45	x	x	x	x	x	x	x	x	x	x	no transducers required
A02 Frog Gastrocnemius	45					x	x				x	SS12LA p27, STIM p25, ELSTM2 p38
A03 Frog Sciatic Nerve	45					x	x				x	STIM p25, NERVE1/2 p38, BSLCBL p39
A04 Frog Heart	45			x		x	x				x	SS12LA p27, (HDW p31)
A05 Visceral Smooth Muscle	45			x+		x+	x+				x+	SS12LA p27, (HDW p31), (ITBS100 p33)
A06 Cockroach Ventral Nerve	45					x	x				x	BSLCBL8/9 p36, EL452 x3 p38
A07 Q <sub>10</sub> Principle (Dissolved O <sub>2</sub> -Goldfish)	45					x*	x*					SS69L p31 or BSL-TCI16 p39
A08 Action Potential in Earthworm	45					x	x				x	STIM p25, NERVE1/2 p38, BSLCBL p39
A09 Properties of Turtle Heart Cardiac Muscle	45					x	x				x	SS12LA p27, STIM p25, STM2+452 p38, CBL8 p36
A10 Regulation of Circ. & Resp: Homeostasis	45											SS2L p25, SS13L p27, SS31L p29
A11 Resting Potential - Crayfish Muscle	45					x†	x†				x†	BSLCBL8 p36, (CBL204 x2 p39)
A12 Membrane Potential (muscle)	45					x-	x-				x-	BSLCBL8 p36, (CBL204 x2 p39)
A13 Active Transport (frog skin)	45											BSLCBL8 p36, BSLCBL10 p38
A14 Central Pattern Generator (hornworm pupa)	45					x	x				x	BSLCBL8 p36, EL452 x3 p38
A15 Earthworm Smooth Muscle	45			x		x	x				x	SS12LA p27, (HDW p31), (ITBS100 p33)

\*interface included, probe required

+tissue bath required

†requires Crawdad CD-ROM Lab Manual

-glass microelectrode required

- Advanced System, 9**
- Airflow & Respiration**
- Accessories (Mouthpieces, Nose Clips, Filters)
    - Student Consumables BSL-ACCPACK, 34
    - SSI 1LA Accessories, 34
    - High-Performance Accessories, 35
  - Calibration Syringe
    - 600 ml AFT6, 34
    - 2 liter AFT26, 35
  - Couplers AFT11 series, 35
  - Differential Pressure Trans. SS40L-SS42L, 30
  - Facemask with T-valve AFT25, 35
  - Pneumogram Trans. (sm. animal) SS67L, 31
  - Pneumotach Trans. (medium flow) SSI 1LA, 27
  - Pneumotach Trans. Series (wide flow range),
    - SS45L-52L, 33
  - Tubing AFT7, 35
  - T-valve AFT21, 35
- Animal Physiology, 12**
- Basic System, 8**
- Bioimpedance (Cardiac Output) Sensor SS31L, 29**
- Biology, 14**
- Biomechanics, 16**
- Biomedical Engineering (BME), 20**
- Biopac Student Lab Overview, 2-7**
- Biopotential Electrodes, 36-38**
- For ECG, EDA (GSR), EEG, EGG, EMG, EOG
- Blood Pressure**
- Animal or *In Vitro* SS13L, 27; NIBP200A, 33
  - Cuff Trans. SS19L, 27
  - Noninvasive Monitoring NIBP100A, 32
  - See also Sound Detection
- Breadboard (Signal Processing) SS39L, 30**
- BSL Biopac Student Lab Overview, 2-7**
- Cables**
- High-impedance BSLCBL8/9, 36
  - Interface, 39
  - Nerve Recording BSLCBL3A/4B, 39
  - Nerve Stimulating BSLCBL1/2A, 39
  - USB CBLUSB, 40
  - Voltage Drive & Current Monitor BSLCBL10, 38
- Cardiac Output Sensor SS31L, 29**
- Consumables AFT series, 34**
- Conversion (Mac↔PC) PLAT30 series, 40**
- Core Packages, 8-23, 41**
- Customer Loyalty Program (Upgrades), contact BIOPAC**
- Data Acquisition Unit MP35, 24**
- Differential Pressure Trans. SS40L-SS42L, 30**
- Displacement Trans. SS14L, 28**
- Dissolved O<sub>2</sub> Probe Trans. SS69L, 31**
- Electrodes, 36-38**
- Accessories (Gel, Tape, Abrasive Pads), 36
  - Adhesive Disks/Collars ADD200, 36
  - Bar Electrode Series EL350, 37
  - Disposable Series EL500, 37
- Leads**
- Clip Leads, Recording LEAD140/141/142, 36
  - Clip Leads, Stimulation BSLCBL7/11/12, 38
  - Multi-lead (12-lead ECG) Cable SS29L, 27
  - Pinch Leads LEAD110 series, 36
  - Shielded Adapter SSI 1LA, 36
  - Shielded General-purpose Lead Set SS2L, 25
  - Needle Series EL450, 38
  - Reusable Series EL250, 37
  - Stimulating, for animals ELSTM2, 38
  - Stimulating, human-safe HSTM01, 28
- EDA (GSR) Electrodermal Activity**
- Disposable Setup SS57L, 26
  - Reusable Trans. SS3LA, 28
  - Electrodes with Isotonic Gel EL507, 37
- Exercise Physiology, 16**
- Finger Twitch Trans. SS61L, 30**
- Force**
- Fixed-range Trans. SS63L-SS66L, 31
  - Hand Dynamometer SS25LA, 27
  - Tension Adjuster HDW100A, 31
  - Variable Range Trans. SS12LA, 27
- Gas Analysis**
- CO<sub>2</sub> and O<sub>2</sub> Analysis GASSYS2, 34
  - Calibration Gas, 35
  - Regulator, 35
- Goniometers SS20L-SS24L, 29**
- GSR Galvanic Skin Response, see EDA**
- Hardware, 23-42**
- Table: Hardware Used in Lessons, 46
- Headphones OUT1, 26**
- Heel/Toe Strike SS28L, 29**
- Human Physiology, 10**
- Interface**
- Input, BNC/DSub 9, SS9L and SS70L, 39
  - Input, 3.5 mm/DSub 9, BSLCBL5, 39
  - Output, BNC/3.5 mm, BSLCBL6, 38
  - Output, DSub 9/BNC, OUT2, 39
  - Custom Adapters SS-KIT, 39
  - Research Amplifier Adapter BSLCBL14, 39
  - Trans. Connector Interface BSL-TCI series, 39
  - Y-adaptor CBL204, 39
- Lab Manual BIOPAC MANBSL3S, 40**
- Lessons Summary, 43-45**
- Micromanipulator, MANIPULATOR, 38**
- Manuals**
- BIOPAC MANBSL3S, 40
  - Published 3rd-party Manuals, 8
- Movement**
- Accelerometers SS26L-SS27L, 29
  - Finger Twitch Trans. SS61L, 30
  - Goniometers SS20L-SS24L, 29
  - Heel/Toe Strike Trans. SS28L, 29
  - Reflex Hammer Trans. SS36L, 30
- MP35 Data Acquisition Unit, 24**
- Multi-lead ECG SS29L, 27**
- Nerve Chambers NERVE1/2, 38**
- Neurophysiology, 18**
- Noninvasive Blood Pressure, 32-33**
- Nursing Programs, see Human Physiology, 10**
- pH Probe Trans. SS68L, 31**
- Pharmacology, 22**
- Platform Change PLAT30, 40**
- Pneumogram Trans. (sm. animal) SS67L, 31**
- Power, AC series, 40**
- Pressure Trans. SS13L, 27**
- Pulse Plethysmogram Trans. SS4LA, 26**
- Psychophysiology, 18**
- Reflex Hammer Trans. SS36L, 30**
- Respiratory Effort Trans. SS5LB, 26**
- Safety Notice, 23, 24**
- Signal Processing (Breadboard) SS39L, 30**
- Simple Sensor Transducers, 26-40**
- Software Overview, 2-7**
- Sound Detection**
- Headphones OUT1, 26
  - Microphone, Physiological Sounds SS17L, 28
  - Microphone, Speech Frequency SS62L, 30
  - Stethoscope, Electronic Trans. SS30L, 28
  - Tubeophone OUT101, 31
- SS Series (Simple Sensor) Transducers, 26-40**
- Stimulation**
- Clip Leads, BSLCBL7/11/12, 38
  - Human-safe Electrode HSTM01, 28
  - Needle Electrodes, for animal, ELSTM2, 38
  - Nerve Chambers NERVE1/2, 38
  - Stimulators BSLSTM and SS58L, 25
  - Stimulus Presentation, SuperLab, STP35W, 32
- Strip Electrodes, 37**
- Stroboscope TSD122, 32**
- SuperLab System, STP35W, 32**
- Switches**
- Digital Switches SS53L-SS55L, 30
  - Hand Switch (analog) SS10L, 26
- Temperature Sensors**
- Fast-response Thermistor SS6L, 26
  - Waterproof or Liquid Immersion, SS7L-SS8L, 28
- Tissue Bath Stations & Accessories, 33**
- Transducers, SS Series, 26-40**
- Transducer Accessory "TA" Packs, 42**
- Toxicology, 22**
- Ultimate System, 9**
- Upgrade**
- Customer Loyalty Program, contact BIOPAC
  - Lesson 16 & 17 Hardware, BSLUPGI6&17, 40
- USB Interface CBLUSB, 40**
- Variable Assessment Trans. SS43L, 30**
- Y-adaptor CBL204, 39**



# Biopac Student Lab users

A partial list of satisfied Biopac Student Lab users follows. If you would like a list tailored to identify users in your area or your application of interest, contact BIOPAC at [info@biopac.com](mailto:info@biopac.com)

## United States

Alma College  
Augusta State University  
Austin Peay State University  
Bakersfield College  
Barry University  
Baylor College of Medicine  
Baylor University  
Bellarmine University  
Bethel College  
Birmingham Southern College  
Boston University  
Bowling Green State University  
Brevard Community College  
Brigham Young University  
Bristol Heart Institute  
Bucknell University  
Cal State University:  
    Bakersfield; Fullerton;  
    Los Angeles; Northridge;  
    Sacramento; Stanislaus  
California Lutheran Univ.  
California Polytechnic:  
    Pomona; San Luis Obispo  
Case Western Reserve Univ.  
Catawba College  
Claremont McKenna College  
Clemson University  
Cleveland State University  
College of Holy Cross  
College of the Siskiyous  
Columbus State University

Cornell University  
Cosumnes River College  
Cuesta College  
Dartmouth College  
Diablo Valley College  
Dixie College  
Duke University  
Eastern Oregon University  
Elizabethtown College  
Emory University  
Everett Community College  
Evergreen State College  
Florida International University  
Florida State University  
Georgia College & State Univ.  
Georgia Institute of Technology  
Georgia State University  
Golden West College  
Gordon College  
Harvard University  
Hofstra University  
Imperial Valley College  
Indiana State University  
Indiana Univ. - Purdue Univ.  
Iowa State University  
Ithaca College  
Ivy Tech State College  
Jacksonville State University  
Jefferson Community College  
Johns Hopkins Medical School  
Kent State University  
LA Mission College  
Lake Superior College

Lehigh University  
Lenoir-Rhyne College  
Life Chiropractic College West  
Linfield College  
Loma Linda University  
Louisiana State University  
Louisiana Tech University  
Loyola College  
Millaps College  
Mississippi Valley State Univ.  
Monterey Peninsula College  
Mount Holyoke College  
Mount Vernon Nazarene Coll.  
Nassau Community College  
National Institutes of Health  
New York Medical College  
Ohio State  
Ohio University  
Oregon Institute of Technology  
Parker College of Chiropractic  
Pasadena City College  
Pierce College, California  
Pierce College, Washington  
Pima Community College  
Purdue University  
Quinnipiac College  
Rice University  
Rochester Institute of  
    Technology  
Rollins College  
Rutgers University  
Sacred Heart University  
San Diego State University

San Francisco State University  
San Jose State University  
Santa Barbara City College  
Santa Monica College  
Santa Rosa Junior College  
Seattle Pacific University  
Shasta College  
Sierra College  
Skidmore College  
Solano Community College  
South Mountain Community  
    College  
Stanford University  
State University of NY  
Stony Brook University  
Taft College  
Texas A&M  
Tulane Univ. School of Medicine  
Tuskegee University  
University of Arizona  
University of California:  
    Berkeley; Riverside; San  
    Francisco; Santa Barbara  
University of Colorado  
University of Connecticut  
University of Delaware  
University of Florida  
University of Georgia  
University of Hawaii, Hilo  
University of Illinois  
University of Kentucky  
University of Maryland  
University of Massachusetts

University of Memphis  
University of Michigan  
University of Minnesota  
University of Mississippi  
University of Missouri  
University of Montana  
University of Nebraska  
University of New England  
University of New Orleans  
University of North Carolina,  
    Chapel Hill  
University of North Texas  
University of Oklahoma  
University of Pennsylvania  
University of Pittsburgh  
University of San Francisco  
University of South Alabama  
University of South Carolina:  
    Aiken; Columbia; Spartanburg  
University of Texas  
University of Utah  
University of Vermont  
University of Virginia  
University of Wisconsin  
University of Wyoming  
Victor Valley College  
Wake Forest University  
Washington University  
Whittier College  
Yale University  
Yavapai College  
Yuba College

## International

Abant İzzet Baysal Üniversitesi  
Acadia University  
Adnan Menderes Üniversitesi  
All India Institute of Medical  
    Sciences  
American University College of  
    Science & Technology, Beirut  
Ankara University  
Aston University  
Bar-Ilan University  
Baskent University  
Bernier Fachhochschule  
Brunel University  
Bryansk State Petrovsky  
    University  
Camosun College  
Cardinal Tien Nursing School  
CEGEP de Sainte-Foy  
CEGEP de Trois-Rivières  
Chang Gung Nursing  
    Technology College  
Charles Sturt University  
Charles University  
China Medical College  
Christian-Albrechts-Universität  
    zu Kiel  
CHU Reims  
Chung Shan Medical University  
CNAM  
CNRS  
Collège Ahuntsic  
Collège de Sherbrooke  
College of the Fraser Valley  
Concordia University  
D.Y. Patil College of Pharmacy  
Daegu University  
Diele University  
Dokuz Eylül University  
Eotvos Lorand University  
ENSSAT

Ethiopian Health and Nutrition  
    Research Institute  
FH Dortmund  
FH Joanneum, Graz  
FH München  
Faculté Polytechnique de Mons  
Facultés Universitaires Notre  
    Dame de la Paix  
Fu Jen Catholic University  
Gaziantep University  
Hacettepe University  
Hanlim University  
Hiroshima Shudo University  
Högskolan i Jönköping  
INSERM  
Inter American University of  
    Puerto Rico  
International Islamic University  
Istanbul University  
Johann Wolfgang Goethe-  
    Universität Frankfurt am Main  
Kakitaya University  
Khon Kaen University  
King Saud University  
King's University College  
Kirkkale University  
Klaipeda University  
Kobe Gakuin University  
Lady Harding Medical College  
Leeds Metropolitan University  
Linköpings Universitet  
Ludwig-Maximilians-Universität  
    München  
Ludwik Rydygier Medical  
    University  
Max Planck Institute  
McGill University  
Mid Sweden University  
Nanyang Technological University  
National Taipei Nursing College  
National Taiwan University  
National Technical University of  
    Athens

National Tsing Hua University  
National University of  
    Singapore  
National Yang Ming University  
Netaji Subhas Institute of  
    Technology  
NgeeAnn Polytechnic  
Niigata Sangyo University  
Nizams Institute of Medical  
    Sciences  
Northumbria University  
Obirin University  
Osmangazi University  
Pingtung Technology University  
PSG Institute of Medical  
    Sciences & Research  
Queensland University of  
    Technology  
Rajguru College of Applied  
    Sciences  
Rijksuniversiteit Groningen  
Royal College of Surgeons in  
    Ireland  
Sedaya International College  
Semmelweis Medical University  
Showa University  
Singapore Polytechnic  
Taipei Medical University  
Taipei Municipal Physical  
    Education College  
Tamagawa University  
Technological Educational  
    Institute of Heraklion  
Trinity College Dublin  
Ubolratchathane University  
Universidad Alberto Hurtado  
Universidad Autónoma de  
    Centro América  
Universidad Autónoma de  
    Coahuila  
Universidad Autónoma de  
    Madrid

Universidad de Cádiz  
Universidad de Costa Rica  
Universidad de Extremadura  
Universidad de León  
Universidad de Navarra  
Universidad de Oviedo  
Universidad de Puerto Rico,  
    Cayey  
Universidad de San Carlos de  
    Guatemala  
Universidad de Zaragoza  
Universidad del País Vasco  
Universidad Nacional  
    Autónoma de México  
Universidad Nacional de Mar  
    del Plata  
Universidade Agostinho Neto  
Universidade da Coruña  
Universidade de Lisboa  
Universidade do Porto  
Universidade Federal do  
    Espírito Santo  
Universitas Andalas  
Universitas Katolik Widya  
    Mandala Surabaya  
Universitat Autònoma de  
    Barcelona  
Universitat de Barcelona  
Universitat de les Illes Balears  
Universitat de València  
Universität Freiburg  
Universität Graz  
Universität Linz  
Universität Mainz  
Universität Tübingen  
Université Bordeaux 2  
Université de Bretagne  
    Occidentale  
Université de Genève  
Université de Metz  
Université de Moncton  
Université de Savoie

Université des Antilles et de  
    la Guyane  
Université du Maine  
Université du Québec à  
    Chicoutimi  
Université du Québec à  
    Montréal  
Université Faculté de Médecine  
    Angers  
Université Faculté de Médecine  
    Brest  
Université Faculté de Médecine  
    Nîmes  
Université Faculté de Médecine  
    Vandoeuvre  
Université Paris 8  
Universiti Kebangsaan Malaysia  
Universiti Malaya  
Universiti Putra Malaysia  
Universiti Sains Malaysia  
Universiti Teknologi Malaysia  
Universiti Teknologi Mara  
    University College  
    Northampton  
University of Bergen  
University of Cambridge  
University of Crete  
University of Delhi  
University of Derby  
University of Glamorgan  
University of Glasgow  
University of Gloucestershire  
University of Nottingham  
University of Oxford  
University of Surrey  
    Roehampton  
University of the West  
    of England  
University of Toronto  
University of Vienna  
Waseda University  
Yamagata University



**BIOPAC**  
Systems, Inc.

42 Aero Camino, Goleta, CA 93117 USA  
Tel. 805-685-0066 • Fax: 805-685-0067

## Worldwide distribution for a world-class product!

You may contact an authorized international representative  
directly or through [www.biopac.com](http://www.biopac.com)



### ASIA-PACIFIC

#### Australia and New Zealand

SDR Clinical Technology  
213 Eastern Valley Way  
Middle Cove NSW 2068 Australia  
Tel: +61-2-9958-2688  
Fax: +61-2-9958-2655  
E-mail: [sdrc@sdrc.com.au](mailto:sdrc@sdrc.com.au)  
Web: <http://www.sdrc.com.au>

#### China

Upwards Teksystems Ltd.  
Room 1910, Prime Tower,  
No. 22 Chaowai Ave.  
Chaoyang District  
Beijing 100020 China  
Tel: +86-10-65885160  
Fax: +86-10-65885162  
E-mail: [info@upwardstek.com](mailto:info@upwardstek.com)  
Web: <http://www.upwardstek.com>

#### India

GENTECH  
13/10, II Floor, W.E.A.  
Saraswati Marg, Karol Bagh  
New Delhi 110 005 India  
Tel: +91-11-25814297  
Fax: +91-11-25814301  
E-mail: [viplesht@hotmail.com](mailto:viplesht@hotmail.com)

#### Japan

Monte System Corporation  
3-17, Kandasakuma-cho,  
Chiyoda-ku  
Tokyo 101-0025 Japan  
Tel: +81-3-5823-0191  
Fax: +81-3-5823-0190  
E-mail: [info@monte.co.jp](mailto:info@monte.co.jp)  
Web: <http://www.monte.co.jp>

#### Korea

EWHA Commercial Co.  
587-7 Bongduk 1-Dong, Nam-ku  
Taegu Korea  
Tel: +82-53-474-9455  
Fax: +82-53-474-9457  
E-mail: [ewha86@netsgo.com](mailto:ewha86@netsgo.com)  
Web: <http://www.physio.co.kr>

#### Malaysia and Singapore

Dynamedix Sdn Bhd  
Suite 10-13, Level 10, UOA-II  
No. 21, Jalan Pinang  
50450 Kuala Lumpur  
Malaysia  
Tel: +60-3-2164-7118  
Tel: +65-9696-4602  
(Singapore Support Line)  
Fax: +60-3-2164-9171  
E-mail: [info@dyna-medix.com](mailto:info@dyna-medix.com)

#### Taiwan

Upwards Biosystems Ltd.  
2F-6, No. 100, Sec. 2, Hopping E.  
Road  
Taipei Taiwan  
Tel: 886-2-27335111, -14  
Fax: 886-2-27332552  
E-mail: [ubl@ms9.hinet.net](mailto:ubl@ms9.hinet.net)

#### Thailand

General Enterprises Marketing L.P.  
694 Soi Ratchadaniwet 24  
Pracharatbumpheh Road,  
Huaikhwang  
Bangkok 10320 Thailand  
Tel: +662-274-2456  
Fax: +662-274-2443  
E-mail: [gem@inet.co.th](mailto:gem@inet.co.th)

### EUROPE

#### Bulgaria

ELPAK Medical & Special  
Equipment  
12 Ivan Bogorov Str.  
Varna 9002 Bulgaria  
Tel: +359-52-602360  
Fax: +359-52-699060  
E-mail: [elpak@elpak.bg](mailto:elpak@elpak.bg)  
Web: <http://www.elpak.com>

#### France

CEROM  
155, rue du Fauborg Saint-Denis  
75010 Paris France  
Tel: +33-01.39.85.77.19  
Fax: +33-01.39.85.77.19  
E-mail: [cerom@tiscali.fr](mailto:cerom@tiscali.fr)

#### Germany

med-NATIC  
Harthausen Strasse 21  
D-81545 Munich Germany  
Tel: +49-89-6 42 29 58  
Fax: +49-89-6 42 29 78  
E-mail: [med-Natic@t-online.de](mailto:med-Natic@t-online.de)  
Web: <http://www.med-natic.de>

#### Greece

St.Tsimpos Ltd.  
Medical & Scientific Instruments  
94 Constantinoupolleos Str.  
546 42 Thessaloniki Greece  
Tel: +30-2310-888630  
Fax: +30-2310-888631  
E-mail: [tsimpos@tsimpos.gr](mailto:tsimpos@tsimpos.gr)  
Web: <http://www.tsimpos.gr>

#### Italy

NORIS srl  
Piazza Alpini d'Italia, 11  
Pontetaro (Parma) 43015 Italy  
Tel: +39-0521.615157  
Fax: +39-0521.616071  
E-mail: [noris@noris-mdn.com](mailto:noris@noris-mdn.com)  
Web: <http://www.noris-mdn.com>

#### Portugal

Sporteste, Lda.  
Rua S. João de Brito, 9-A  
2795-180 Linda-a-Velha Portugal  
Tel: +351 21 420 97 10  
Fax: +351 21 420 97 11  
E-mail: [joao.rasoilo@sporteste.pt](mailto:joao.rasoilo@sporteste.pt)

#### Romania

Russi Scientific Instruments  
3, Tudor Vianu Str.  
712363 Bucharest 1 Romania  
Tel: +40.21.231 0538  
Fax: +40.21.230 1634  
E-mail: [info@russi.ro](mailto:info@russi.ro)

#### Russia

Transonic Ltd. Moscow  
Profsojuznaja Str. 96-5-37  
117485 Moscow Russia  
Tel: +7-095 247 0210  
Fax: +7-095 330 6157  
E-mail: [info@transonic.ru](mailto:info@transonic.ru)  
Web: <http://www.transonic.ru>

#### Spain

Bionic Iberica, S.A.  
Mèxic 8, 1-1°  
08320 El Masnou (Barcelona)  
Spain  
Tel: +34-93-540-1909  
Fax: +34-93-555-6921  
E-mail: [info@bionic.es](mailto:info@bionic.es)  
Web: <http://www.bionic.es>

#### Sweden

JoR AB  
Rubanksgatan 3  
Knivsta, AR 74171 Sweden  
Tel: +46-18 34 28 20  
Fax: +46-18 38 09 94  
E-mail: [jor@jorse](mailto:jor@jorse)  
Web: <http://www.jorse>

#### Turkey

COMMAT Ltd.  
Cetin Emec Bulv. 74. Sok. No:4/9  
Ovecler Ankara Turkey  
Tel: +90-312 417 40 28  
Fax: +90-312 472 74 18  
E-mail: [info@commat.com.tr](mailto:info@commat.com.tr)  
Web: <http://www.commat.com.tr>

#### United Kingdom and Ireland

Linton Instrumentation  
Unit 11, Forge Business Centre  
Upper Rose Lane, Palgrave  
Diss, Norfolk, England, IP22 1AP  
United Kingdom  
Tel: +44-1379 651344  
Fax: +44-1379 650970  
E-mail: [mail@lintoninst.co.uk](mailto:mail@lintoninst.co.uk)  
Web: <http://www.lintoninst.co.uk>

### MIDDLE EAST

#### Israel

Bargal Analytical Instruments &  
Software Ltd.  
12 Wallenberg Raoul Street  
Tel Aviv 69719 Israel  
Tel: +972-3-649-0823  
Fax: +972-3-649-0826  
E-mail: [sales@bargal.co.il](mailto:sales@bargal.co.il)  
Web: <http://www.bargal.co.il>

#### Saudi Arabia

Raez Environmental Systems &  
Advanced Technologies - RESAT  
P.O. Box 261242  
Riyadh 11342, Kingdom of  
Saudi Arabia  
Tel: +966-1-478-2295  
Fax: +966-1-474-0794  
E-mail: [raez@resat.org](mailto:raez@resat.org)  
Web: <http://www.resat.org>

### NORTH AND SOUTH AMERICA

#### Brazil

Sellex  
Rua Arandu, 205/1105  
04562-031 São Paulo SP Brasil  
Tel: (011) 5506-4646  
Fax: (011) 5505-7433  
E-mail: [washington@sellex.com](mailto:washington@sellex.com)  
Web: <http://www.sellex.com>


#### Canada

Harvard Apparatus Canada  
6010 Vanden Abeele  
St. Laurent, Quebec  
Canada H4S 1R9  
Tel: (514) 335-0792  
Toll-Free in Canada: 1-800-361-1905  
Fax: (514) 335-3482  
E-mail: [sales@harvardapparatus.ca](mailto:sales@harvardapparatus.ca)

#### Mexico

Biocologia Quimica, S.A. de C.V.  
Luis Gonzalez Obregon No. 16  
Cronistas, CD. Satelite  
53100 Naucalpan, Mexico  
Contact: Juan Resendiz  
Tel: +52-55 5572-3009  
Fax: +52-55 5393-3019  
E-mail: [bioquim@prodigy.net.mx](mailto:bioquim@prodigy.net.mx)  
Web: <http://www.teoloyucan.com>

- Full translation (software and Lab Manual) available.
- Software translation in process; contact your rep for an update.

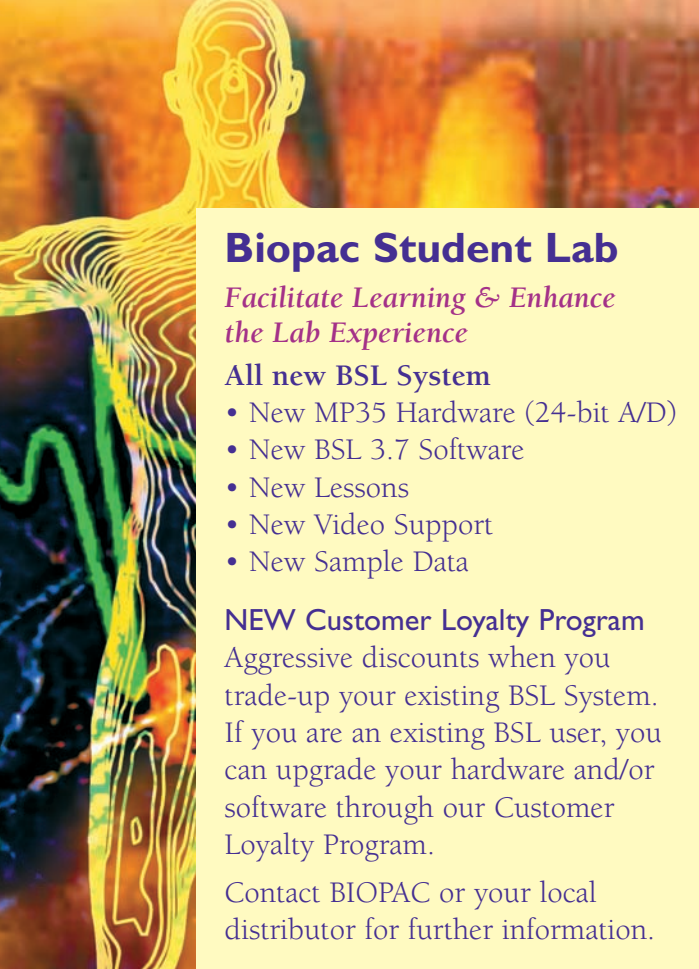

## Biopac Student Lab

*Facilitate Learning & Enhance the Lab Experience*

### All new BSL System

- New MP35 Hardware (24-bit A/D)
- New BSL 3.7 Software
- New Lessons
- New Video Support
- New Sample Data

### NEW Customer Loyalty Program

Aggressive discounts when you trade-up your existing BSL System. If you are an existing BSL user, you can upgrade your hardware and/or software through our Customer Loyalty Program.

Contact BIOPAC or your local distributor for further information.

## The ultimate teaching solution for...

- |                        |                               |
|------------------------|-------------------------------|
| Physiology             | Exercise Phys. & Biomechanics |
| Biology                | Psychophys. & Neurophysiology |
| Biomedical Engineering | Pharmacology & Toxicology     |
| Nursing                | Veterinary Programs           |

## When you get the BSL...they'll get science!

Incorporate clinical techniques & industry standard equipment to teach your students real-world skills.

- |           |                |                    |
|-----------|----------------|--------------------|
| ECG       | Biofeedback    | Nerve Conduction   |
| EEG       | Bioimpedance   | Pulmonary Function |
| EMG       | Cardiac Output | Reaction Time      |
| EOG       | Blood Pressure | Respiration        |
| EDA (GSR) | Heart Sounds   | Temperature        |

**Multi-level learning solutions let you control the material and method of each experiment.**



- Inquiry-based student lessons
- Instructor modified labs
- Advanced analysis tools
- PRO Lessons
- Custom Lessons
- No programming required

*Use for introductory lessons, graduate programs, or advanced research.*

### New hardware includes:

- |  |                                |
|--|--------------------------------|
| Low Voltage Stimulator                           | Airflow Transducers—           |
| Cardiac Output Sensor                            | wide flow range                |
| Tissue Bath Stations                             | Pneumogram Transducer          |
| Nerve Chambers                                   | Signal Processing (Breadboard) |
| CO <sub>2</sub> & O <sub>2</sub> Analysis Module | Variable Response Indicator    |
| Finger Twitch Transducer                         | Speech Freq. Microphone        |
| Differential Pressure Trans.                     | Dissolved O <sub>2</sub> Probe |
|  | Isolated BNC Input Adapter     |

**Plus more! See inside for details or visit [www.biopac.com](http://www.biopac.com)**

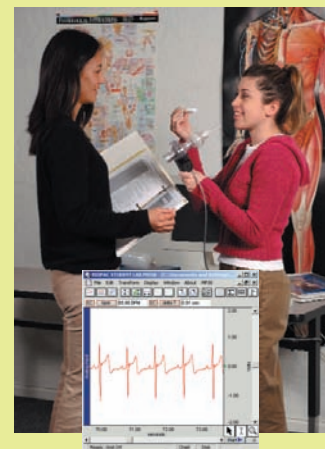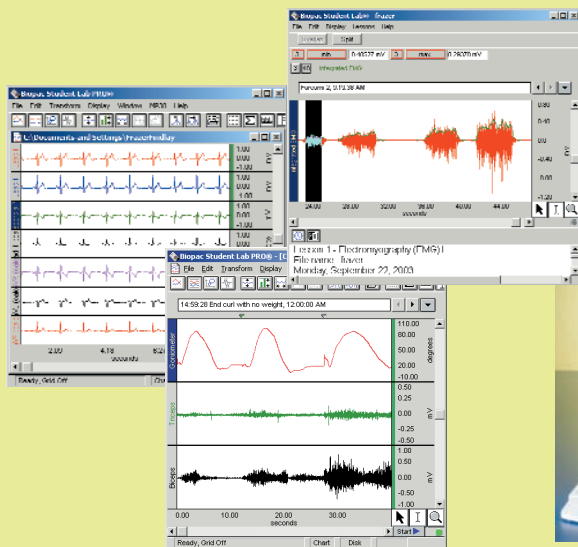

## Complete Solutions for Life Science Education



**BIOPAC**  
Systems, Inc.

42 Aero Camino  
Goleta, CA 93117

See what you've been missing!



www.biopac.com